CONGEN

SureFood[®] QUANTARD Allergen 40

Art. No. S3301
100 rxn

User Manual

February 2020

 Inhalt

1 Allgemeines3

1.1 Beschreibung.....3

1.2 DNA-Präparation4

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1 General Information

1.1 Description

Der SureFood® QUANTARD Allergen 40 is validated for quantitative determination of 1 mg to 400 mg allergenic substance / kg food sample.

The laboratory reference material SureFood® QUANTARD Allergen 40 is based on **corn flour** material. It contains all potentially allergenic food ingredients according to regulation (EU) No 1169/2011 (except for **molluscs, lactose and sulphur dioxide**) in a concentration of **40 mg/kg**.

Content of allergenic substances:

Allergen	Material
Gluten containing cereals	Wheat flour type 40S (<i>Triticum</i> spp.)
Crustacean	Black Tiger shrimps (<i>Penaeus monodon</i>)
Egg	Egg powder , hen (<i>Gallus gallus</i>)
Fish	Humpback salmon (<i>Oncorhynchus gorbuscha</i>)
Peanut	Peanut (<i>Arachis hypogaea</i>), untreated, peeled
Soy	Soybean (<i>Glycine max</i>)
Milk	Milkpowder, cow (<i>Bos taurus</i>)
Edible Nut	Hazelnut (<i>Corylus avellana</i>), peeled Almond (<i>Prunus dulcis</i>), untreated Walnut (<i>Juglans regia</i>), untreated Cashew (<i>Anacardium occidentale</i>), untreated Pecan (<i>Carya illinoensis</i>), untreated Brazil nut (<i>Bertholletia excelsa</i>), peeled Pistachio (<i>Pistacia vera</i>), peeled Macadamia (<i>Macadamia ternifolia</i>), roasted
Celery	Celery seeds (<i>Apium graveolens</i>)
Mustard	Mustard seeds, yellow (<i>Sinapis alba</i>)
Sesame	Sesam seeds (<i>Sesamum indicum</i>)
Lupin	Lupin seeds (<i>Lupinus hartwegii</i>), peeled

The laboratory reference material SureFood® QUANTARD Allergen 40 is validated in combination with the DNA extraction method SureFood® PREP Advanced, protocol 1, and the SureFood® ALLERGEN real-time PCR detection methods.

Further information about the quantification of allergenic substances is given in the manual of the SureFood® ALLERGEN real-time PCR products.

In case of using alternative methods the sample weight and extraction method in combination with the detection system have to be validated.

1.2 DNA-preparation

For DNA-preparation the use of SureFood® PREP Advanced (Art. No. S1053), protocol 1 is recommended.

1.3 Kit components and storage

Reagent	Amount
QUANTARD Allergen 40	1 x 2,0 g

Store all reagents dry at –20°C and protected from light.

Note: To avoid moisture absorption thaws the unopened laboratory reference material at room temperature for at least one hour before use.

1.4 Additionally required equipment and materials

- DNA-Extraction kit (e.g. SureFood® PREP Advanced, Art. No.: S1053)
- Real-time PCR methods (SureFood® ALLERGEN series)

2 Further Information

2.1 Technical Support

For further questions please contact your distributor or send an e-mail to sales@r-biopharm.de.

2.2 Distribution and Ordering

R-Biopharm AG
An der neuen Bergstrasse 17,
64297 Darmstadt, Germany
Phone: +49 (0) 61 51 - 81 02-0
Fax: +49 (0) 61 51 - 81 02-20
E-Mail: orders@r-biopharm.de
www.r-biopharm.com

r-biopharm®