# **RIDA<sup>®</sup> 0.9% NaCl (sterile)**

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# RIDA<sup>®</sup> 0.9% NaCl (sterile)

## 1. Intended use

Sterilized, physiological NaCl solution (0.9%) for activation of all Compact Dry plates prior to application for air sampling or membrane filters (membrane filter samples from liquids).

## 2. General

Beside the suitability of Compact Dry plates for analysis of food samples they are especially suited for incubation of membrane filter samples as well as application as air sampling plates. The plates contain a nutrient pad which has to be moisturized with 1 ml of sterile liquid prior to application.

## 3. Test principle

RIDA<sup>®</sup> 0.9% NaCl (sterile) ampules contain 1 ml of sterilized, physiological saline (0.9%) which can be applied to Compact Dry as moisturizing liquid.

## 4. Reagents provided

Each kit contains sufficient ampules for 150 determinations:

30 x 5 ampules

(containing 1 ml of sterilized, physiological NaCl solution (0.9%))

## 5. Materials required but not provided

- incubator (fixed to 35 °C/95 °F or adjustable in a range between 30 and 40 °C /86 and 104°F)
- Compact Dry plates
- In case of membrane filter samples: sterile membrane filters for microbiological purposes (pore size 0.45 µm, Ø 47 mm); filtration device

## 6. Warnings and precautions for users

RIDA<sup>®</sup> 0.9% NaCl (sterile) is for in-vitro-use only. Do not use it past the expiration date on the kit label.

## 7. Storage instructions

RIDA<sup>®</sup> 0.9% NaCl (sterile) is stable at room temperature until expiration date.

## 8. Indication of instability or deterioration of reagents

Stained or turbid ampules should be discarded.

## 9. Test implementation

1. Open the RIDA<sup>®</sup> 0.9% NaCl ampule (twist off closure)
2. Open lid of the Compact Dry plate and apply the content of the ampule to the nutrient pad
3. Re-cover the Compact Dry plate after application of NaCl solution
4. In case of air sampling:
  - Decontaminate the surface of the sampling area using 70% ethanol solution
  - Place the plate on the sterile surface
  - Open lid and place it on the sterile surface with inner side heading downwards
  - Expose Compact Dry plate to the air for 30 min
  - Close plate, turn it around (upside-down) and incubate following the instructions for use of the respective parameter (s. Compact Dry manuals)
5. In case of membrane filtration:
  - Put membrane filter on top of the filter grit of the filtration device and attach the (decontaminated) filter funnel
  - Pour sample liquid into the funnel and switch on pump of the filtration device
  - After sucking through all liquid, rinse the funnel with 20 ml of sterile water or 0.9 saline and wait until all rinsing liquid has been sucked through, also
  - Remove membrane filter (Caution: vacuum – ventilate collection flask!) and place filter on the moisturized Compact Dry plate with impinged side up
  - Close plate, turn it around (upside-down) and incubate following the instructions for use of the respective parameter (s. Compact Dry manuals)

Manuals containing instructions for use and evaluation of Compact Dry plates are enclosed in each test kit. An information brochure about Compact Dry is available at R-Biopharm.

The data corresponds to our present state of technology and provides information on our products and their uses. R-Biopharm makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. Defective products will be replaced. There is no warranty of merchantability of this product, or of the fitness of the product for any purpose. R-Biopharm shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.

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are registered trademarks of R-Biopharm AG