Promedia ST-25

Art. Nr.: Z0302

Hersteller: ELMEX Ltd., Tokio, Japan

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Promedia ST-25

Art. No.: Z0302

Manufacturer: ELMEX, Ltd., Tokyo, Japan

1. Intended use

Promedia ST-25 is a sampling system consisting of a sterile swab in a tube with 10 ml sterile PBS buffer. The quantitative detection of microorganisms in environmental samples will be enabled in combination with Compact Dry medium plates.

2. General

Promedia ST-25 combines a sampling tool, dilution medium and transportation tool in one device. The swab test allows environmental sampling under sterile conditions and is intended to be used in combination with Compact Dry medium plates for hygiene monitoring.

3. Test principle

The surface sample is taken with the swab stick which is fixed in the cap of the tube. After sampling the PBS buffer serves as dilution medium. The sample is blended by shaking the tube and 1 ml sample aliquots are poured through the drop cap of the tube onto one of the dry medium systems. Inoculated medium plates are incubated specified conditions and examined for typical colonies.

4. Reagents provided

Each package unit contains 10 Promedia ST-25.

5. Materials required but not provided

- Incubator (35 °C / 95 °F minimum): CULTURA[®] Mini-Incubator, Art. No. ZC7140651
- Compact Dry plates (Art. No. HS7311 – HS9802) – in principle all available types of plates are suitable for usage in combination with Promedia ST-25
- Optional: Frames for definition of 100 cm² (Frame 100 cm², Art. No. ZFR1600000)

6. Warnings and precautions for users

The Promedia ST-25 is for in-vitro-use only. After expiration date (s. package label) there is no guarantee for functionality anymore.

7. Storage instructions

The Promedia ST-25 is stable at room temperature until expiration date on the package label.

8. Indication of instability or deterioration of reagents

Tubes with stained or turbid buffer should be discarded.

9. Test implementation

1. Write sample ID on the label of the buffer tube.
2. Open the screw cap of the buffer tube and remove the swab.
3. Wipe the sampling surface (approx. 100 cm², usage of Frame 100 cm² recommended) using the swab and place the swab back into the buffer. Close the screw cap thoroughly.
4. Blend the sample by shaking the tube.
5. Dispense 1 ml of sample buffer onto the dry medium plate using the dropper opening in the tube cap (1 ml graduation on the tube; 1 ml $\hat{=}$ approx. 20 drops). Further 1 ml aliquots can be plated onto different medium plates (e.g. Compact Dry CF for Coliforms, Compact Dry X-SA for *S. aureus*, etc.)
6. Incubate the medium plates under specified conditions (s. individual Compact Dry test instructions).
7. Specify and count the grown colonies according to the corresponding information in the test instructions of the respective Compact Dry product.

Detailed instructions for the use of Compact Dry products are included in the respective test kits.

An overview about all available Compact Dry products, package sizes and product codes is given in the Compact Dry brochure as well as in the R-Biopharm product catalogue.

The data corresponds to our present state of technology and provides information on our products and their uses. R-Biopharm makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. Defective products will be replaced. There is no warranty of merchantability of this product, or of the fitness of the product for any purpose. R-Biopharm shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.

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R-Biopharm AG