# **RIDASCREEN<sup>®</sup> Risk Material 10/5**

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# RIDASCREEN® Risk Material 10/5

## Brief information

RIDASCREEN® Risk Material 10/5 (Art. No.: R6703) is a sandwich enzyme immunoassay for the quantitative analysis of risk material (CNS) in and on raw meat, meat products and on contaminated surfaces.

All reagents required for the enzyme immunoassay - including standards - are contained in the test kit.

The test kit is sufficient for 96 determinations (including standards).

A microtiter plate spectrophotometer is required for quantification.

Technicians need no specialized training to perform the RIDASCREEN® Risk Material 10/5 test, however, free support is offered from the distributor on request, if necessary.

Sample preparation: place the sampling swab into the sample dilution buffer

Time requirement: sample preparation  
(for 10 meat samples)..... approx. 10 min  
test implementation (incubation time)..... 15 min

Limit of detection:  $\leq 0.1$  % for CNS tissue  
in raw meat and meat products

Specificity: The RIDASCREEN® Risk material 10/5 kit allows detection of CNS tissue from cattle, veal, sheep, goat, horse, poultry and pig.  
A specification of the kind of tissue (brain, spinal cord or eye) is not possible.

Note: For the determination of risk material (CNS) in processed (heated) meat and sausages we refer to the RIDASCREEN® Risk Material test , Art. No.: R6701.

## 1. Intended use

RIDASCREEN® Risk Material 10/5 is a sandwich enzyme immunoassay for the quantitative analysis of risk material (CNS) in / on raw meat, meat products and on contaminated surfaces.

## 2. General

BSE (bovine spongiform encephalopathy) has been diagnosed first in Great Britain in 1986. The molecular mechanisms of infection are still unresolved, but it is clear, that the Prion protein (prP<sup>Sc</sup>) plays a major role in the transmission and development of the brain disease. Feeding of insufficient heated animal meal is likely to be the main cause of the spread of disease, which has caused more than 180000 dead cows up to the end of 2000. Application of extracts from infected brain induced infection in different animal species. Similar diseases are now described in cows, sheep, goat, cat and many other mammals. With a high presumption, the new variant of the human Creutzfeldt-Jakob disease (vCJD) is caused by infected beef and beef products. The mortality of this disease is 100 %, with 111 deaths in Great Britain up to May 2002. As a protective measurement, animal meal / protein as animal feed is no longer allowed (2000/766/EG) and potential risk material has been banned from food production by the EU (2000/418/EG) / (2001/999/EG).

## 3. Test principle

The basis of the test is the antigen-antibody reaction. The detection of risk material ensures by the determination of Glial Fibrillary Acidic Protein (GFAP), a cellular marker, which can be found very high concentrated in CNS tissue. The wells of the microtiter strips are coated with specific antibodies against GFAP. If the sample is contaminated with risk material, the contained glial fibrillary acidic proteins will bind to the specific capture antibodies. The bound GFAP is detected by an antibody peroxidase-conjugate directed against GFAP (enzyme conjugate). Any unbound enzyme conjugate is then removed in a washing step. After substrate/chromogen is added to the wells, bound enzyme conjugate converts the chromogen into a blue product. The addition of the stop solution leads to a color change from blue to yellow. The measurement is made photometrically at 450 nm. The absorbance is proportional to the risk material concentration in the sample.

## 4. Reagents provided

Each kit contains sufficient materials for 96 measurements (including standard analyses). Each test kit contains:

- 1 x Microtiter plate (12 strips with 8 removable wells each)  
coated with antibodies against GFAP
- 4 x Standard solutions (1.3 ml each)  
GFAP in aqueous solution  
corresponding to 0 % (zero standard), 0.1 %, 0.2 %, 0.4 %  
risk material (CNS) in samples  
ready to use
- 1 x Conjugate (6 ml).....red cap  
peroxidase conjugated antibody against GFAP  
ready to use
- 1 x Substrate/chromogen, 10 ml ..... brown cap  
stained red, contains tetramethylbenzidine
- 1 x Stop reagent (14 ml).....yellow cap  
contains 1 N sulfuric acid
- 1 x Sample Dilution Buffer (10 ml)..... white cap  
as a 10x concentrate  
contains 0.5 % SDS
- 1 x Washing buffer (Salt)  
for preparation of a 10 mM Phosphate Buffer, pH 7.4  
contains 0.05 % Tween 20
- 2 x Pasteur Pipette  
1 ml, graduated  
for placing sample dilution buffer

## 5. Materials required but not provided

### 5.1. Equipment:

- microtiter plate spectrophotometer (450 nm)
- variable 20 - 200 µl- and 200 - 1000 µl-micropipettes

- swabs
  - test tubes
- } available as part of a Sampling Tool kit  
(Art. No.: Z0010) from your local distributor

## 6. Warnings and precautions for the users

The stop solution contains 1 N sulfuric acid (R36/38, S2-26).

The sample dilution buffer contains SDS. Avoid contact of the reagent with eyes, skin and clothing.

## 7. Storage instructions

Store the kit at 2 - 8 °C (35 - 46 °F). Do not freeze any test kit components.

Return any unused microwells to their original foil bag, reseal them together with the desiccant provided and further store at 2 - 8 °C (35 - 46 °F).

The substrate/chromogen solution is light sensitive, therefore, avoid exposure to direct light.

No quality guarantee is accepted after the expiration date on the kit label.

Do not interchange individual reagents between kits of different lot numbers.

## 8. Indication of instability or deterioration of reagents

- any bluish coloration of the reddish substrate/chromogen solution prior to test implementation
- a value of less than 0.4 absorbance units ( $A_{450\text{ nm}} < 0.4$ ) for standard 4 (0.4 %)

## 9. Preparation of Samples

### 9.1. Raw meat and meat products

- if samples are homogenous, insert the dry swab several times into the sample; if samples are solid, swab the surface thoroughly while rotating the swab
- squeeze the swab thoroughly in a test tube containing 1 ml diluted sample dilution buffer (see 10.1.)  
(alternatively 50 mg of sample can be transferred into diluted sample dilution buffer (see 10.1.))
- use 50 µl per well in the assay

## 9.2. Surface swabbing

- pre-moisten the swab by placing it into diluted sample dilution buffer (see 10.1.)
- swab a representative area (10 x 10 cm, resp. 4" x 4")
- squeeze the swab thoroughly in a test tube containing 1 ml diluted sample dilution buffer (see 10.1.)
- use 50 µl per well in the assay

### Storage of samples:

Diluted in sample dilution buffer the samples can be stored at 2 - 8 °C (36 - 46 °F) for approx. 2 days. Bring samples to room temperature (20 - 25 °C / 68 - 77 °F) and mix shortly before use.

## 10. Test implementation

### 10.1. Preliminary comments

Bring all reagents to room temperature (20 - 25 °C / 68 - 77 °F) before use.

The **sample dilution buffer** is provided as a 10fold concentrate. Therefore the buffer concentrate has to be diluted 1:10 (1+9) with dist. water before use. Dissolve eventually occurring crystals in a water bath at 37 °C (98 °F) before dilution. We recommend dilution of the buffer concentrate in only one step (i.e. 10 ml concentrate + 90 ml dist. water).

The diluted buffer can be used up to the expiry date printed on the buffer label if stored at 2 - 8 °C (36 - 46 °F).

For one determination place 1 ml of the diluted buffer into each test tube by using the graduated pipettes contained in the kit.

A PBS-Tween buffer is needed as **washing buffer**, please use the washing buffer salt (pouch) contained in the kit (see 4.). The contents is dissolved in one liter of distilled water. The ready to use washing buffer expires after approx. 4 - 6 weeks at 2 - 8 °C (36 - 46 °F).

Alternatively: Dissolve the contents of the envelope in 100 ml of distilled water to obtain a 10fold concentrated washing buffer. Use 1 part of this concentrate and dissolve with 9 parts of distilled water to obtain the ready to use washing buffer.

The 10fold concentrate expires after approx. 8 - 12 weeks, store at room temperature (20 - 25 °C / 68 - 77 °F).

## 10.2. Test procedure

Carefully follow the recommended washing procedure. Do not allow microwells to dry between working steps.

1. Insert a sufficient number of wells into the microwell holder for all standards and samples to be run. Record standard and sample positions. Do not use more than 4 strips with 8 wells (28 samples + 4 standards) at a time.
2. Pipet 50 µl of standard solution or prepared sample to separate wells.
3. Add 50 µl of enzyme conjugate (red cap) to the bottom of each well, mix gently by shaking the plate manually and incubate for 10 min (+/- 1 min) at room temperature (20 - 25 °C / 68 - 77 °F)
4. Pour the liquid out of the wells and tap the microwell holder upside down vigorously (three times in a row) against absorbent paper to ensure complete removal of liquid from the wells. Fill all the wells with 250 µl washing buffer (see 10.1.) and pour out the liquid again. Repeat two more times.
5. Add 100 µl of substrate/chromogen (brown cap) to each well. Mix gently by shaking the plate manually and incubate for 5 min (+/- 1 min) at room temperature (20 - 25 °C / 68 - 77 °F) in the dark.
6. Add 100 µl of stop reagent (yellow cap) to each well. Mix gently by shaking the plate manually and measure the absorbance at 450 nm. Read within 15 min after addition of stop reagent.

## 11. Results

The color obtained for the standards serves as the reference value.

A special software, the RIDA<sup>®</sup>SOFT Win (Art. No. Z9999), is available to evaluate the RIDASCREEN<sup>®</sup> enzyme immunoassays. The software allows automated data retrieval from the reader and result calculation by linear regression.

The course of the standard curve is shown in the Quality Assurance Certificate enclosed in the test kit.

### Graphic evaluation:

The absorbance values of the standards are entered in a system of coordinates on graph paper against the risk material contents [%]. The values of the standards are connected with a line and the risk material concentration in % depending on the absorbance of each sample can be read from the calibration curve.

**Remark:**

The limit of detection of the RIDASCREEN® Risk Material 10/5 kit is  $\leq 0.1$  % for CNS tissue (spinal cord) in raw meat and meat products. The limit of quantification is 0,2 % for CNS tissue (spinal cord) in raw meat and meat products. When results are interpreted it has to be considered, that the contents of marker proteins in different risk materials (brain, spinal cord, eye) may vary. For example, brain may give positive results in higher concentrations than spinal cord.

R-Biopharm makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. If any materials are defective, R-Biopharm will provide a replacement product. There is no warranty of merchantability of this product, or of the fitness of the product for any purpose. R-Biopharm shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.

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are registered trademarks of R-Biopharm AG