Compact Dry LS medium for *Listeria spec* / Medium zum Nachweis von Listerien / milieu pour listeria / medio para listeria / medio per listeria / meio para listeria

100 plates/Platten/plaques/placas/lastre/placas

Art-No. HS8811

40 plates/Platten/plaques/placas/lastre/placas

Art-No. HS8812

English

Deutsch

Français

Compact Dry LS is a ready to use, chromogenic plate for detection of Listeria spec.

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Sample pretreatment

Viable count in water or liquid foodstuff

Drop 1 ml of specimen (dilute if necessary) on the middle of the Compact Dry plate.

Viable count in solid foodstuff

Add buffer solution to the sample and homogenize by stomacher®. Drop 1 ml of specimen (dilute if necessary) on the middle of the dry sheet of the Compact Dry plate.

Viable count in swab test specimen

Use a steril wet swab to wipe the surface which should be analyzed and put it back into the device with wiping solution. Drop 1 ml of wiping solution (dilute if necessary) on the middle of the Compact Dry plate. Use the Compact Dry Swab (Art. No. ZCS1002953, content 40 pcs) as single application (1 ml) or Promedia ST-25 (Art. No. Z0302, content 10 pcs) as multiple application (10 ml) for this procedure.

Test instructions

1. Open the cap and drop 1 ml of specimen on the middle of the Compact Dry plate.
2. Specimen diffuses automatically and evenly into the sheet and transforms the dried sheet into a gel within seconds.
3. Put the cap again on the plate and write the information needed on the memorandum section.
4. Turn over the capped plate and put in the incubator.
5. After incubation count the number of colored colonies underneath the plate. White paper placed under the plate helps to count the colonies.

Incubation time 24 ± 2 hours
plus a further 24 ± 2 hours
Incubation temperature 35 or 37 ± 1 °C

Interpretation of the results

Listeria species forms light blue / blue colonies of 1~2 mm in diameter by chromogens contained in a medium.

Storage and shelf life

Keep at room temperature (+ 1 to + 30 °C).
Total shelf life 12 months after manufacturing.

Notes

- It is known that *Listeria ivanovii* tends to grow slowly and *Listeria seeligeri* tends to be inhibited in this plate.
- A combination of selective agents in the medium inhibits the growth of other bacteria.
- High concentrations on plates (> 300 cfu) will cause the entire growth area to become blue / green. In this case dilute the sample.
- After use please follow the current disposal regulations.
- The growth area is 20 cm². The back of the plate has a grid carved of 1 cm x 1 cm to make the colony counting easier. In case of any difficulties to count colonies due to large number of colonies grown, total viable count can be obtained by multiplying 20 by an average number of colonies per grid counted from several grids.
- Compact Dry plates are produced at an ISO 9001 certified site.

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