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VitaFast® Pantothenic Acid Spiking standard

Brief information

Ca-pantothenate standard for the use of spiking food, feed and pharmaceutical products.

1. Reagents provided

3 x Ca-pantothenate spiking standard (solid)

Content: **see Quality Assurance Certificate**

Remark: after the expiry date no guarantee of quality.

2. Required reagents, not provided

Redist. or deionized water for the preparation of the spiking standard solution.

3. Warning and precautions for the user

No precautions necessary.

4. Storage instructions

Store the reagents at 2 - 8 °C (35.6 - 46.4 °F).

5. Preparation of the spiking standard solution

- open the Ca-pantothenate spiking solution bottle, place the lid down with the opening facing upwards
- add **x ml (x = see Quality Assurance Certificate „Spiking standard“ or label „Spiking standard“)** redist. or deionized water;
the content corresponds to 120 µg pantothenic acid per 1 ml (or 12 µg per 100 µl);
this means: 110.4 µg active pantothenic acid per 1 ml (or 11.04 µg per 100 µl)
- close the bottle with the lid and dissolve the standard by shaking = spiking solution (**expiry 1 day in the dark**)
- add the needed volume to the sample (see example below) and prepare the sample according to the product information VitaFast® Pantothensäure / Pantothenic Acid

The sample should be spiked with 50 % or 100 % of the vitamin content in the sample.

Example for a spiking procedure:

Estimated active pantothenic acid content in the sample: 6 mg / 100 g; take 1 g sample, this corresponds with a pantothenic acid content of 60 µg / 1 g sample (the calculation refers to 1 g weighed sample on 40 ml).

100 % spiking:

Spiking with 100 % (of the content of the sample):

Weigh 1 g sample + add 544 µl spiking solution (corresponds 60 µg absolute, respectively 6 mg / 100 g)

Calculation of the estimated total pantothenic acid content:

Estimated content of the sample + content of the 100 % spiking

6 mg / 100 g (estimated content) + 6 mg / 100 g (spike) = 12 mg / 100 g active pantothenic acid (total content)

Prepare the spiked sample according to **6. Sample preparation** of the product information insert VitaFast® Pantothensäure / Pantothenic Acid (Art. No. P1005).