---

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# QuickGEN PCR Kit

## P1 Screening

### 1. Examples of successful detected microorganisms

#### Bacteria

Lactobacillus spp.

Lactococcus lactis

Megasphaera spp.

Micrococcus kristinae

Pectinatus spp.

Pediococcus spp.

#### Yeast

Saccharomyces spp.

Dekkera spp.

Zygosaccharomyces spp.

Rhodotorula spp.

Torulaspora spp.

Kluyveromyces spp.

Debaromyces spp.

Candida spp.

Hanseniaspora spp.

Pichia spp.

## 2. Intended use

Qualitative detection of bacterial and yeast contaminations in filtrated beer and beer mixes. For products with high amount of yeast use the Screening Kit without yeast detection to find bacterial contaminations more sensitive (Q033).

## 3. Test principle

The TaqMan® real-time PCR is based on hot-start-PCR and sequence-specific dual labelled probes (FAM/DQ; HEX/DQ; ROX/DQ) which, when accurately hybridised, emit a measurable fluorescent signal of a defined wavelength in the extension phase. The increase of signal is continuously measured in a real-time PCR detection instrument. To avoid false negative PCR-results an Inhibition Control (CY5/DQ) is amplified together in one reaction vessel with the specific sequence. The system contains dUTP. Optional: the use of Uracil-N-Glycosylase will eliminate any contamination with Uracil containing amplicons from former PCRs (the enzyme is not part of this kit). **The tubes contain lyticase.**

## 4. Kit contents

The kit contains sufficient reagents for 48 PCR reactions:

|   |                |
|---|----------------|
| 1 x Premix  | white cap      |
| 48 x Dye Strips (freeze-dried, incl. IC-DNA and lyticase) | tube strips    |
| 6 x Cap Strips  |                |
| 1 x ddH <sub>2</sub> O                                    | colourless cap |
| 1 x Control-DNA (freeze-dried)                            | yellow cap     |

## 5. Storage conditions

**The Control-DNA is freeze-dried and has to be solved in ddH<sub>2</sub>O prior to use (see 7.1).**

Do **not** freeze the lyophilized Dye Strips and lyophilized Control-DNA.

The PCR reagents should be stored at 2– 8 °C (35 – 46 °F). Keep Premix for storage at - 20 °C (- 4 °F) after opening. Avoid loss of sensitivity by repeating freezing and thawing more than 3 times. For irregular use aliquot the Premix.

The Dye Strips contain the fluorescent labelled probes and should be handled light protected. All reagents are stable for 12 months, if they are stored correctly.

## **6. Materials required, but not provided**

### **6.1. Instruments**

Real-time PCR instrument for low profile PCR tubes

Centrifuge for 1.5 – 2.0 mL reaction tubes

Centrifuge for strips

Pipettes

“Vortex”

### **6.2. Reagents and plastic ware**

sterile ddH<sub>2</sub>O

sterile filter tips

**optional:** Uracil N-Glycosylase (0.01 U/μL added to the PCR reaction)

Colour Compensation Set for LC480 (GEN-IAL<sup>®</sup>, REF Q800)

**LC480: Use a Colour compensation prior to first use of the kit**

## 7. PCR

### 7.1. PCR-Setup

*When using the kit for the first time, the freeze-dried Control-DNA has to be shortly centrifuged and carefully resolved:*

- add 55  $\mu\text{L}$  sterile ddH<sub>2</sub>O to the freeze-dried Control-DNA
- after 15 minutes mix well

**Before every use thoroughly mix all PCR-components and centrifuge briefly.**

Remove the film from the needed tube strips and pipette the PCR-components. After pipetting close the tube strips with the provided cap strips.

PCR-reaction Setup:

| PCR-components | volume ( $\mu\text{L}$ ) |
|----------------|--------------------------|
| Premix         | 15.0                     |
| Sample-DNA     | 5.0*                     |
| Total volume   | <b>20.0</b>              |

\* if using the Simplex Easy DNA- or QuickGEN Yeast Sample Preparation Kit:  
add 2.5  $\mu\text{L}$  DNA and 2.5  $\mu\text{L}$  sterile ddH<sub>2</sub>O

1. Pipette 15  $\mu\text{L}$  of the Premix into each PCR-tube, making sure that, prior to the first filling, the tip has been moistened.
2. Add 5  $\mu\text{L}$  sample DNA, add 5  $\mu\text{L}$  of the Control-DNA for the positive control reaction, add 5  $\mu\text{L}$  of the extraction control and 5  $\mu\text{L}$  of ddH<sub>2</sub>O for the negative control\* reaction. Use a fresh tip with each DNA filling.
3. Close the tubes immediately and centrifuge them shortly.
4. Place the tubes in the PCR-machine and start run.

**Very important: \* Please fill up the negative control with 5  $\mu\text{L}$  ddH<sub>2</sub>O to avoid unspecific amplification.**

**Work swiftly to avoid warming up and keep away from light.**

## 7.2 PCR-Program

### 7.2.1 PCR-Program LC480

1. Click the button **tool** on the right side in the window **LightCycler 480 Software release 1.5.0. SP1**
2. Click the button **Detection formats** at the left side of the menu bar
3. Click **New** in the window **Detection formats** and name the experiment
4. Open the window **Filter Combination Selection** and choose the following filter combinations: 465-510 / 533-580 / 533-610 / 618-660.
5. Open the window **Selected Filter Combination** list and add the following amounts

| Excitation Filter | Emission Filter | Name        | Melt factor | Quant factor | Max. Integ. Time |
|-------------------|-----------------|-------------|-------------|--------------|------------------|
| 465               | 510             | 465-510 Fam | 1           | 10           | 2                |
| 533               | 580             | 533-580 Hex | 1           | 10           | 2                |
| 533               | 610             | 533-610 610 | 1           | 10           | 2                |
| 618               | 660             | 618-660 660 | 1           | 10           | 2                |

6. **Close** the window
7. Click the button **New Experiment** on the right side of the menu bar
8. From the pull-down menu **Detection formats** choose the defined experiment, click the button **Customize** and check the detection formats. All of them have to be activated
9. Click the button **ok**
10. Define the following program:

1. Program Name: **Heat**

Cycles **1** Analysis Mode **None**

| Target (°C) | Acquisition Mode | Hold (hh:mm:ss) | Ramp Rate (°C/s) | Acquisitions (per °C) | Sec Target (°C) | Step size (°C) | Step Delay (cycles) |
|-------------|------------------|-----------------|------------------|-----------------------|-----------------|----------------|---------------------|
| 37          | None             | 00:15:00        | 4.40             |                       | 0               | 0              | 0                   |
| 95          | None             | 00:15:00        | 4.40             |                       | 0               | 0              | 0                   |

2. Program Name: **Ampli**

Cycles **30\*-35** Analysis Mode **Quantification**

| Target (°C) | Acquisition Mode | Hold (hh:mm:ss) | Ramp Rate (°C/s) | Acquisitions (per °C) | Sec Target (°C) | Step size (°C) | Step Delay (cycles) |
|-------------|------------------|-----------------|------------------|-----------------------|-----------------|----------------|---------------------|
| 95          | None             | 00:00:10        | 4.40             |                       | 0               | 0              | 0                   |
| 60          | Single           | 00:00:20        | 2.20             |                       | 0               | 0              | 0                   |

**\* for preenriched samples program 30 cycles**

3. Program name : **Cool**

Cycles **1** Analysis Mode **None**

| Target (°C) | Acquisition Mode | Hold (hh:mm:ss) | Ramp Rate (°C/s) | Acquisitions (per °C) | Sec Target (°C) | Step size (°C) | Step Delay (cycles) |
|-------------|------------------|-----------------|------------------|-----------------------|-----------------|----------------|---------------------|
| 40          | None             | 00:00:20        | 2.20             |                       | 0               | 0              | 0                   |

**Optional:** Saving the program as **run template**. The template button allows to select and apply a template to the currently open object and to save the currently open object as a template. Click the clamp beside the button **Apply Template** . Click **Save as template** and save the file in **templates**.

11. Click the button **Subset editor** on the left side
12. Click the button **+** and **New Subset 1** appears
13. Mark the wells in the **New Subset 1 Settings** window
14. Click the button **Apply**
15. Click the button **Sample editor** on the left side
16. **Very important:** Activate In the window **Step 1 Select Workflow: Abs.Quant**
17. Choose the subset **New Subset** in the window **Step 2 Select Samples**
18. Define your probes in the **Sample table**
19. Switch to the button **Experiment** and start the run with the button **start run**

## 8. Evaluation

The evaluation has to be made according to the data analysis program recommended by the real-time instrument manufacturer.

LC480: For analysis activate Colour Compensation

**Lactobacillus/ Pediococcus-DNA:** FAM-channel (LC480: 465-510)

**Megasphaera/ Pectinatus-DNA:** ROX-channel (LC480: 533-610)

**Yeast-DNA:** HEX-channel (LC480: 533-588)

**Inhibition Control-DNA:** CY5-channel (LC480: 610-660)

A sample is ***Lactobacillus/ Pediococcus* positive**, if there is a detectable fluorescence increase in the **FAM-channel (LC480: 465-510)** and the negative controls show no amplification. The positive controls should have a positive fluorescence signal. The Inhibition Control in the CY5-channel (LC480: 610-660) may be positive or negative (depending on the amount of DNA or inhibitors in the sample reaction). In the negative controls it has to be positive.

A sample is ***Megasphaera/ Pectinatus* positive**, if there is a detectable fluorescence increase in the **ROX-channel (LC480: 533-610)** and the negative controls show no amplification. The positive controls should have a positive fluorescence signal. The Inhibition Control in the CY5-channel (LC480: 610-660) may be positive or negative (depending on the amount of DNA or inhibitors in the sample reaction). In the negative controls it has to be positive.


A sample is **yeast positive**, if there is a detectable fluorescence increase in the **HEX-channel (LC480: 533-588)** and the negative controls show no amplification. The positive controls should have a positive fluorescence signal. The Inhibition Control in the CY5-channel (LC480: 610-660) may be positive or negative (depending on the amount of DNA or inhibitors in the sample reaction). In the negative controls it has to be positive.

A sample is **negative**, if there is no detectable fluorescence increase in the different channels and the positive controls have a positive fluorescence signal. The negative controls show no amplification. The Inhibition Control in the CY5-channel (LC480: 610-660) has to be positive in the sample and in the negative controls, a false negative result due to inhibitory effects is then excluded.

# analysis flowchart

## LC480: analysis after activating Colour Compensation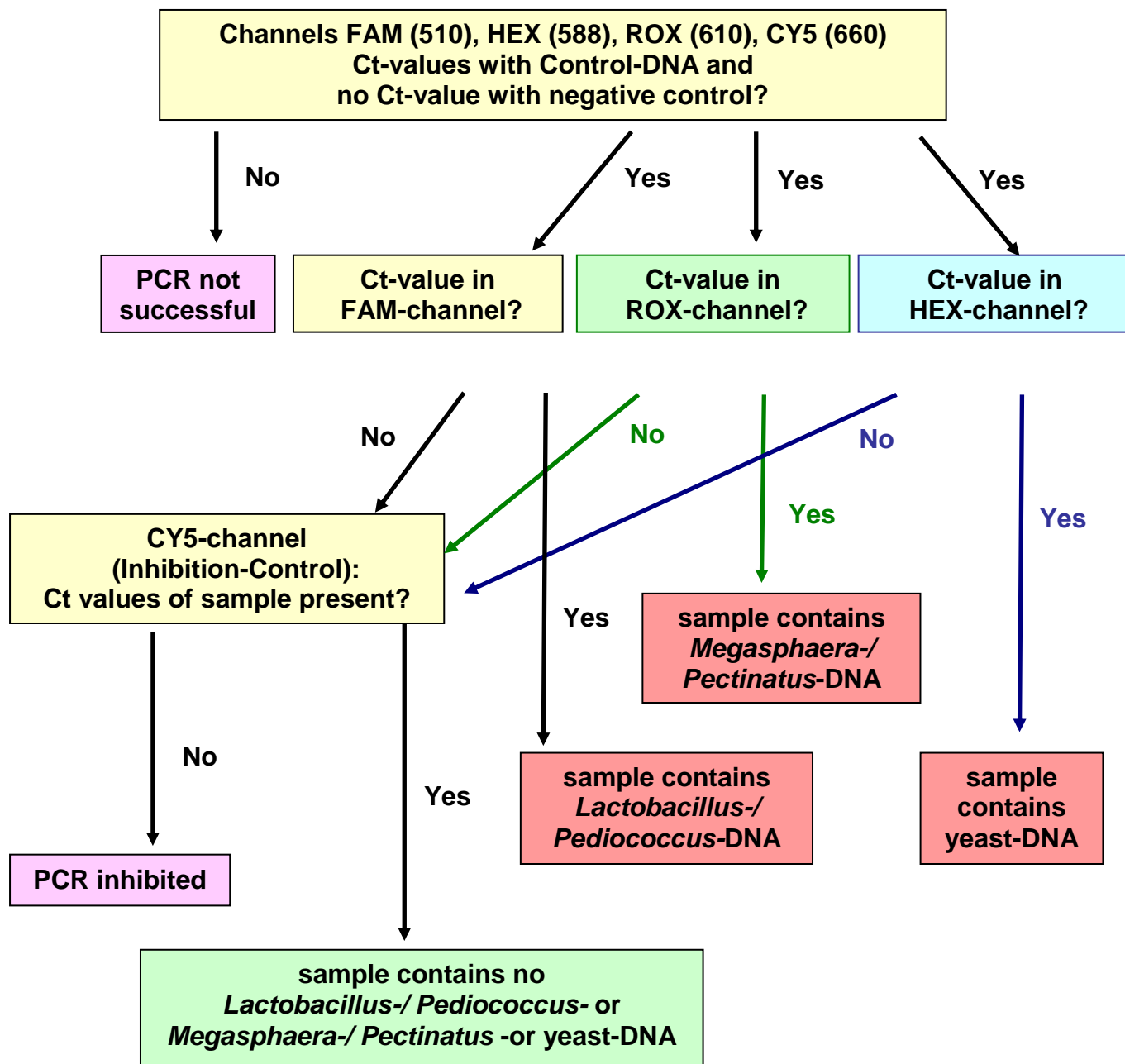

**Note:**

The polymerase-chain reaction (PCR) is protected by patents and requires a licence from Hoffmann-LaRoche Inc..

The provided product does not authorise the purchaser for the commercial use of this method.

GEN-IAL makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. If any materials are defective, GEN-IAL will provide a replacement product. There is no warranty of merchantability of this product, or of the fitness of the product for any purpose. GEN-IAL shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.