
**CONGEN**

# **SureFast<sup>®</sup> Yersinia 3plex**

Art. No. F5132  
100 rxn

## **User Manual**

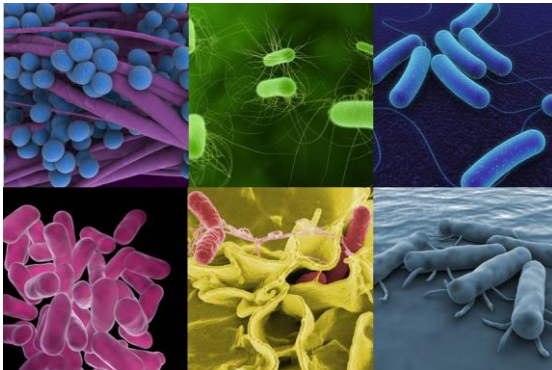**February 2020**

 **Inhalt**

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 **Content**

|       |   |    |
|-------|---|----|
| 1     | General Information .....                           | 8  |
| 1.1   | Description .....                                   | 8  |
| 1.2   | Limit of Detection .....                            | 8  |
| 1.3   | DNA-preparation .....                               | 9  |
| 1.4   | Kit components and storage .....                    | 9  |
| 1.5   | Additionally required equipment and materials ..... | 9  |
| 1.6   | Setup .....   | 9  |
| 1.7   | Detection channel Set-up .....                      | 10 |
| 2     | Qualitative Analysis .....                          | 10 |
| 2.1   | Protocol .....                                      | 11 |
| 2.1.1 | Preparation of the master-mix .....                 | 11 |
| 2.1.2 | Preparation of the real-time PCR-mix .....          | 11 |
| 2.2   | Interpretation of results .....                     | 12 |
| 3     | Further Information .....                           | 12 |
| 3.1   | Product Information .....                           | 12 |
| 3.2   | Technical Support .....                             | 12 |

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## **1 General Information**

### **1.1 Description**

The SureFast® Yersinia 3plex is a real-time PCR for the direct, qualitative detection and differentiation of specific *ail* gene DNA sequences of *Yersinia pseudotuberculosis* and *Yersinia enterocolitica*.

Each reaction contains an internal amplification control (IAC). If the DNA contains PCR inhibiting substances, the signal of the amplification control will be affected or the amplification will be suppressed. Examples for PCR inhibiting substances are alcohols (e.g. ethanol, isopropanol), surfactants (e.g. CTAB, SDS, Triton X100) and salts (e.g. sodium chloride). In addition spices, herbs, algae, cocoa and further sample matrices might have PCR inhibiting effects.

The real-time PCR assay can be performed with commonly used real-time PCR instruments, equipped for detection of four fluorescence emissions at the channels FAM, VIC/HEX and Cy5 at the same time. The technical validation of instruments was performed on Agilent Mx3005P, Agilent AriaDx, Bio-Rad CFX96, Roche LightCycler® 480 II, Applied Biosystems 7500, Qiagen Rotor-Gene Q and R-Biopharm RIDA®CYCLER.

### **1.2 Limit of Detection**

The SureFast® Yersinia 3plex real-time PCR has a limit of detection of  $\leq 5$  DNA copies.

The assay limit of detection depends on sample matrix, processing grade, DNA preparation and DNA content.

The SureFast® PCR systems are very sensitive and therefore even a small amount of target DNA is sufficient for a successful analysis. The concentration of total DNA in the sample does not allow a conclusion on the quantity and quality of the target DNA.

### **1.3 Cross Reactivity**

The detection system of *Yersinia pseudotuberculosis* shows a cross reactivity to *Yersinia pestis*.

**1.4 DNA-preparation**

For DNA-preparation the use of SureFast® PREP Bacteria is recommended. To assess the process of bacterial growth, it is suggested to compare the samples at the beginning and at the end of the culturing (bacterial growth at Cp difference > 3). For cultivation the method of the ISO 18867 is recommended.

**1.5 Kit components and storage**

| Kit Code | Reagent          | Amount      | Lid Color  |
|----------|------------------|-------------|------------|
| 1        | Reaction Mix     | 2 x 1100 µl | Yellow     |
| 2        | Taq Polymerase   | 1 x 80 µl   | Dark Red   |
| 3        | Positive Control | 1 x 200 µl  | Light Blue |

**Store all reagents at –20°C and protected from light. The Taq Polymerase can be stored at +2 to +8°C for multiple uses on the same day.**

**Note: The Taq Polymerase may be in a frozen or unfrozen state. This does not affect the quality of the Taq Polymerase or the performance of the real-time PCR.**

**1.6 Additionally required equipment and materials**

- DNA-Extraction kit (e.g. SureFast® PREP Bacteria Art. No. F1021 )
- real- time PCR instrument with four detection channels (510 nm, 580 nm, 610 nm and 660 nm)
- real-time PCR consumable (plates, tubes, capillaries, foils, caps)
- pipettes with filter tips
- powder-free disposable gloves
- Vortex mixer
- microcentrifuge with a rotor for the reaction tubes

**1.7 Setup**

|  | <b>Blockcycler &amp;<br/>R-Biopharm RIDA®CYCLER</b> | <b>Rotorcycler &amp;<br/>LightCycler® 480 II</b> |
|--|---|--|
| Initial Denaturation (HOLD)            | 1 min, 95°C   | 1 min, 95°C                                      |
| Cycles                                 | 45  | 45   |
| Denaturation                           | 15 sec, 95°C  | 10 sec, 95°C                                     |
| Annealing/Extension (CYCLE)            | 30 sec, 60°C  | 15 sec, 60°C                                     |
| Temperature Transition Rate/ Ramp Rate | Maximum   | Maximum  |

**1.8 Detection channel Set-up**

| Real-time PCR device               | Detection                    | Detection channel | Quencher | Note  |
|------------------------------------|------------------------------|-------------------|----------|---|
| <b>Agilent Mx3005P</b>             | <i>Y. pseudotuberculosis</i> | FAM               | +        |   |
|                                    | IAC                          | HEX               | +        |   |
|                                    | <i>Y. enterocolitica</i>     | Cy5               | +        |   |
| <b>Applied Biosystems 7500</b>     | <i>Y. pseudotuberculosis</i> | FAM               | None     | Check the passive reference option ROX is none. |
|                                    | IAC                          | VIC               | None     |   |
|                                    | <i>Y. enterocolitica</i>     | Cy5               | None     |   |
| <b>Bio-Rad CFX96</b>               | <i>Y. pseudotuberculosis</i> | FAM               | +        |   |
|                                    | IAC                          | VIC / HEX         | +        |   |
|                                    | <i>Y. enterocolitica</i>     | Cy5               | +        |   |
| <b>R-Biopharm RIDA®CYCLER</b>      | <i>Y. pseudotuberculosis</i> | green             | +        |   |
|                                    | IAC                          | yellow            | +        |   |
|                                    | <i>Y. enterocolitica</i>     | red               | +        |   |
| <b>Qiagen Rotor-Gene Q</b>         | <i>Y. pseudotuberculosis</i> | green             | +        |   |
|                                    | IAC                          | yellow            | +        |   |
|                                    | <i>Y. enterocolitica</i>     | red               | +        |   |
| <b>Roche LightCycler® 480 II</b>   | <i>Y. pseudotuberculosis</i> | 465-510           | +        |   |
|                                    | IAC                          | 533-580           | +        |   |
|                                    | <i>Y. enterocolitica</i>     | 618-660           | +        |   |
| <b>Roche cobas® z 480 Analyzer</b> | <i>Y. pseudotuberculosis</i> | 465-510           | +        |   |
|                                    | IAC                          | 540-580           | +        |   |
|                                    | <i>Y. enterocolitica</i>     | 610-670           | +        |   |

## 2 Qualitative Analysis

### 2.1 Protocol

#### 2.1.1 Preparation of the master-mix

Calculate the total number of reactions needed (samples and control reactions) for the specific PCR assay as well as for the inhibition control.

Recommended control reactions for the specific PCR assay: negative control, extraction control, positive control. For the analysis of enrichments additional controls are recommended: zero control (sample before enrichment) and medium control. The reaction mix contains an internal amplification control (IAC) per reaction.

#### **Reactions needed for the qualitative *Y. pseudotuberculosis* and *Y. enterocolitica* detection:**

3 reactions for controls (1x no-template control, 1x extraction control, 1x positive control)

For each sample: at least 1 reaction for each sample DNA

#### **Reactions needed for the qualitative *Y. pseudotuberculosis* and *Y. enterocolitica* detection in enrichments:**

5 reactions for controls (1x no-template control, 1x extraction control, 1x positive control, 1x zero control, 1x medium control)

For each sample: at least 1 reaction for each sample DNA

It is also recommended to prepare the master-mix with 10 % additional volume in order to compensate reagent loss. Allow the reagents to thaw, mix and centrifuge before opening and use.

#### **Example for the calculation and preparation of 10 reactions:**

| Components of the master-mix | Amount per reaction | 10 reactions (with 10% excess) |
|------------------------------|---------------------|--------------------------------|
| Reaction Mix                 | 19.3 µl             | 212.3 µl                       |
| Taq Polymerase               | 0.7 µl              | 7.7 µl                         |
| <b>Total volume</b>          | <b>20 µl</b>        | <b>220 µl</b>                  |

**Mix each master-mix well and centrifuge shortly before use.**

#### 2.1.2 Preparation of the real-time PCR-mix

- Pipette 20 µl of the master-mix into appropriate tubes/wells.
- Close the negative control (the negative control is ready for PCR without any addition).
- Pipette 5 µl of sample DNA into the designated tubes/wells and close them.
- Pipette 5 µl of Positive Control into the designated tubes/wells and close them.
- Centrifuge all tubes/plates or capillaries shortly at low speed.
- Place tubes/plates into the real-time PCR instrument and start the run according to the setup.

## 2.2 Interpretation of results

The evaluation has to be made according to the usual analysis program recommended by the real-time PCR instrument manufacturer.

The control reactions have to show the correct results.

*Y. pseudotuberculosis* DNA is detected in the FAM-channel and *Y. enterocolitica* DNA is detected in the Cy5-channel (see table). In the VIC/HEX-channel the amplification control is detected.

A sample is stated **positive** for the respective parameter, if the sample DNA shows amplification in the respective channel.

A sample is stated **negative** for the respective parameter, if the sample DNA shows no amplification in the respective channel and if the internal control (VIC/HEX-channel) of the sample is **positive** with a shift in Cp-value  $\leq 2$  compared to the negative control.

If the sample DNA in the VIC/HEX-Channel shows **no amplification** or a shift in Cp-value  $> 2$  compared to the negative control, it contains PCR inhibiting substances. A significant decrease in the fluorescence signal can also show the presence of PCR inhibiting substances. Under these circumstances DNA isolation and purification of the sample need to be improved. Alternatively the DNA can be diluted (recommendation 1:2 in PCR-water) and analysed again for inhibition. Please note that the dilution factor also affects the detection limit of the specific *Y. pseudotuberculosis* or *Y. enterocolitica* PCR assay.

| Result in the respective channel            |   |                        | Interpretation                            |
|---|---|------------------------|---|
| FAM channel<br><i>Y. pseudotuberculosis</i> | Cy5 channel<br><i>Y. enterocolitica</i> | VIC/HEX channel<br>IAC |   |
| <b>positive</b>                             | negative                                | <b>positive</b>        | <i>Y. pseudotuberculosis</i> DNA detected |
| negative                                    | <b>positive</b>                         | <b>positive</b>        | <i>Y. enterocolitica</i> DNA detected     |
| negative                                    | negative                                | negative               | invalid                                   |

## 3 Further Information

### 3.1 Product Information

- Detailed information about setup of several real-time PCR devices  
(Download: [www.congen.de/en/company/downloads](http://www.congen.de/en/company/downloads))
- Validation Report upon request

### 3.2 Technical Support

For further questions please send an e-mail to [info@congen.de](mailto:info@congen.de).