CONGEN

SureFood[®] GMO ID DAS-40278-9 Corn

Art. No. S2140
100 rxn

User Manual

June 2020

 **Inhalt**

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Content

1	General Information	8
1.1	Description	8
1.2	Limit of Detection	8
1.3	DNA-preparation	9
1.4	Kit components and storage	9
1.5	Additionally required equipment and materials	9
1.6	Setup.....	10
1.7	Detection channel Set-up	10
2	Qualitative Analysis	11
2.1	Protocol	11
2.1.1	Preparation of the master-mix	11
2.1.2	Preparation of the real-time PCR-mix	11
2.2	Interpretation of results	12
3	Further Information	12
3.1	Product Information.....	12
3.2	Technical Support	12
3.3	Distribution and Ordering	12

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1 General Information

1.1 Description

The SureFood® GMO ID DAS-40278-9 Corn is a real-time PCR for the direct, qualitative detection of a specific genetically modified corn DNA sequence. Therefore the kit contains an event specific real-time PCR system for the DAS-40278-9 corn (OECD unique identifier DAS-40278-9).

The specific detection is according to the validated method of the European Commission.

The real-time PCR assay can be performed with commonly used real-time PCR instruments. The technical validation of instruments was performed on LightCycler® 480 II, Qiagen Rotor-Gene Q, Bio-Rad CFX96, R-Biopharm RIDA®CYCLER, LTF MyGo Pro, Agilent AriaDx and Agilent Mx3005P.

1.2 Limit of Detection

The SureFood® GMO ID DAS-40278-9 Corn real-time PCR has a limit of detection of ≤ 5 DNA copies.

This is equivalent to approx. 0.01 % for unprocessed corn grain.

The assay limit of detection depends on sample matrix, processing grade, DNA preparation and DNA content.

The SureFood® PCR systems are very sensitive and therefore even a small amount of target DNA is sufficient for a successful analysis. The concentration of total DNA in the sample does not allow a conclusion on the quantity and quality of the target DNA.

1.3 DNA-preparation

For DNA-preparation of raw material the use of SureFood® PREP Basic (Art. No. S1052) and for highly processed food and feed the use of SureFood® PREP Advanced (Art. No. S1053) is recommended. SureFood® PREP Add On (Art. No. S1055) is intended to be used for the extraction of DNA from raw materials as well as processed food and feed with sample weight of 2 g. It is used in conjunction with the SureFood® PREP Basic.

1.4 Kit components and storage

Kit Code	Reagent	Amount	Lid Color
1	Reaction Mix	2 x 1100 µl	Yellow
2	Taq Polymerase	1 x 80 µl	Dark Red
3	Positive Control	1 x 200 µl	Light Blue

Store all reagents at –20°C and protected from light. The Taq Polymerase can be stored at +2 to +8°C for multiple uses on the same day.

Note: The Taq Polymerase may be in a frozen or unfrozen state. This does not affect the quality of the Taq Polymerase or the performance of the real-time PCR.

1.5 Additionally required equipment and materials

- DNA-Extraction kit
(e.g. SureFood® PREP Basic Art. No. S1052 / SureFood® PREP Advanced Art. No. S1053 / SureFood® PREP Add On Art. No. S1055)
- real-time PCR instrument
- real-time PCR consumable (plates, tubes, foils, caps)
- pipettes with filter tips
- powder-free disposable gloves
- Vortex mixer
- micro centrifuge with a rotor for the reaction tubes

1.6 Setup

	Blockcycler & R-Biopharm RIDA®CYCLER	Rotorcycler & LTF MyGo Pro
Initial Denaturation (HOLD)	5 min, 95°C	1 min, 95°C
Cycles	45	45
Denaturation	15 sec, 95°C	10 sec, 95°C
Annealing/Extension (CYCLE)	30 sec, 60°C	15 sec, 60°C
Temperature Transition Rate/ Ramp Rate	Maximum	Maximum

1.7 Detection channel Set-up

Real-time PCR device	Detection	Detection channel	Quencher	Note
Agilent Mx3005P	DAS-40278-9 corn	FAM	+	
Agilent AriaDx	DAS-40278-9 corn	FAM	+	
Bio-Rad CFX96	DAS-40278-9 corn	FAM	+	
R-Biopharm RIDA®CYCLER	DAS-40278-9 corn	green	+	
LTF MyGo Pro	DAS-40278-9 corn	FAM	+	
Qiagen Rotor- Gene Q	DAS-40278-9 corn	green	+	Note: Please use only 0.1 ml reaction tube. The gain settings must be set to 5 (factory default) for all channels.
Roche LightCycler® 480 II	DAS-40278-9 corn	465-510	+	

2 Qualitative Analysis

2.1 Protocol

2.1.1 Preparation of the master-mix

Calculate the total number of reactions needed (samples and control reactions) for the specific PCR assay as well as for the inhibition control.

Recommended control reactions for the specific PCR assay: negative control, extraction control, positive control and inhibition control.

For the preparation of the inhibition control the use of the SureFood® GMO SCREEN 4plex 35S/NOS/FMV+IAC kit (Art. No. S2126), the SureFood® GMO SCREEN 35S/NOS/FMV kit (Art. No. S2026) or the SureFood® GMO Plant 4plex Corn/Soya/Canola+IAC (Art. Nr. S2158), respectively, is recommended.

Reactions needed for the qualitative DAS-40278-9 corn detection:

3 reactions for controls (1x no-template control, 1x extraction control, 1x positive control)

For each sample: at least 1 reaction for each sample DNA

It is also recommended to prepare the master-mix with 10 % additional volume in order to compensate reagent loss. Allow the reagents to thaw, mix and centrifuge before opening and use.

Example for the calculation and preparation of 10 reactions:

Components of the master-mix	Amount per reaction	10 reactions (with 10% excess)
Reaction Mix	19.3 µl	212.3 µl
Taq Polymerase	0.7 µl	7.7 µl
Total volume	20 µl	220 µl

Mix each master-mix well and centrifuge shortly before use.

2.1.2 Preparation of the real-time PCR-mix

- Pipette 20 µl of the master-mix into appropriate tubes/wells.
- Close the negative control (the negative control is ready for PCR without any addition).
- Pipette 5 µl of sample DNA into the designated tubes/wells and close them.
- Pipette 5 µl of Positive Control into the designated tubes/wells and close them.
- Centrifuge all tubes/plates shortly at low speed.
- Place tubes/plates into the real-time PCR instrument and start the run according to the setup.

2.2 Interpretation of results

The evaluation has to be made according to the usual analysis program recommended by the real-time PCR instrument manufacturer.

The control reactions have to show the correct results.

A sample is stated **positive**, if the sample DNA shows amplification.

A sample is stated **negative**, if the sample DNA shows no amplification and if the external inhibition control of the sample is **positive** with a shift in Cp-Value ≤ 2 compared to the negative control.

If the sample DNA in the external inhibition control shows **no amplification** or a shift in Cp-value > 2 compared to the negative control, it contains PCR inhibiting substances. A significant decrease in the fluorescence signal can also show the presence of PCR inhibiting substances. Under these circumstances DNA isolation and purification of the sample need to be improved. Alternatively the DNA can be diluted (recommendation 1:2 in PCR-water) and analysed again for inhibition. Please note that the dilution factor also affects the detection limit of the specific DAS-40278-9 corn PCR assay.

3 Further Information

3.1 Product Information

- Detailed information about setup of several real-time PCR devices
(Download: www.congen.de/en/company/downloads)
- Validation Report upon request

3.2 Technical Support

For further questions please contact your distributor or send an e-mail to sales@r-biopharm.de.

3.3 Distribution and Ordering

R-Biopharm AG
An der neuen Bergstrasse 17,
64297 Darmstadt, Germany
Phone: +49 (0) 61 51 - 81 02-0
Fax: +49 (0) 61 51 - 81 02-20
E-Mail: orders@r-biopharm.de
www.r-biopharm.com