bioavid Absorptionspuffer / Absorbent buffer

Art. Nr. BS810-15

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Art. Nr. BS810-15

INTENDED USE

The absorbent buffer (Art. No. BS810-15) is an extraction buffer that can be used for the processing of food samples in the bioavid rapid allergen tests. It is particularly recommended for samples containing polyphenols (e.g. coffee, cocoa, red wine) or samples containing high amounts of oil. It is also suitable for neutralizing samples with a slightly acidic or alkaline pH (the pH value of the extract should be in the neutral range (pH 6 - 8) for testing).

It can be assumed that the absorbent buffer is also suitable for extracting other foods; this must be checked by the user before use.

1. STORAGE

Store the test kit at 2 - 25 °C (35 - 77 °F).

2. REAGENTS PROVIDED

Each kit contains sufficient materials for 15 measurements.

Content	Description	Amount
Sample tube	Containing 9 mL absorbent puffer each	15 x

3. WARNINGS AND PRECAUTIONS FOR THE USERS

Bring the buffer to room temperature (20 - 25 °C / 68 - 77 °F) before use.

The product / test is only suitable within the scope of its intended use.

Do not use the test kit after the expiration date (see test kit label).

All reagents and materials must be recovered or disposed after use at customer's own responsibility according to the protection of human health and the environment. Please observe the applicable national regulations concerning waste disposal (e.g. Waste Management Act, Regulations on Dangerous Chemicals, etc.).

4. PREPARATION OF SAMPLES

Shake the test tube before use to suspend the insoluble components.

1. Homogenize (grind thoroughly to powder and mix well or mix well a solution respectively) a sufficient amount (e.g. 50 g or 50 mL) to ensure taking a representative test portion of sample.
2. Transfer 1 g (or in case of liquid samples 1 mL) of homogenized sample into a sample tube.
3. Mix for 10 - 15 minutes (overhead shaker, rotator or similar).
4. Centrifuge at 2,000 - 3,000 x g for 3 - 5 minutes or filter.
5. After centrifugation or filtration, a sediment of approx. 0.5 - 1 mL volume is visible.
6. 0.1 mL of the supernatant is used in the test for testing for allergens (for further procedure, see product-specific instruction for use chapter "Test procedure").

Note: As a use to neutralize liquid samples with a slightly acidic or alkaline pH, no incubation (step 3) is required. Immediately after addition and mixing of 1 mL sample, centrifugation or filtration can be performed.

5. LIMITATION

Due to the large amount of different foods, it is not possible for the manufacturer of a test kit to validate all foods for the test. The suitability of the respective foodstuff for use with the absorbent buffer must be verified by the user.

Version overview

Version number	Chapter and title
2022-12-15	Release version
2024-02-01	Revision
2024-08-21	Current version Changes made: <ul style="list-style-type: none">– Adaptation Intended use: Neutralization only for slightly acidic and alkaline samples– Change of “Disclaimer”

DISCLAIMER

1. In conformance with the German Civil Code (“BGB”) R-Biopharm AG provides a limited warranty (“Gewährleistung”) against defects in design and manufacture present at delivery that make the Product unsuitable or unsafe for the contractually specified Product use and location through properly qualified personnel. Such limited warranty warrants Product title and against such material Product defects for 12 months, or if the Product is provided with a shelf life, the length of the declared shelf life, calculated from the date of risk transfer pursuant to delivery terms, and further provided customer provides timely and proper notice of the defect.
ALL OTHER WARRANTIES OR GUARANTIES, EXPRESS OR IMPLIED, OF ANY KIND ARE EXCLUDED, WHETHER IMPLIED BY CUSTOM, PRACTICE, THE COURSE OF DEALING BETWEEN THE PARTIES OR OTHERWISE. The responsibility for any express or implied warranties or extensions of R-Biopharm’s own warranty made by dealers, distributors, or resellers is solely with such dealer, distributor, or reseller, and R-Biopharm AG expressly disclaims liability for such warranties.
2. R-Biopharm Products are designed for sole use by properly trained and qualified personnel applying appropriate industry standards and practices. Defects are covered only to the extent that such defects are demonstrably the result of defective material, design, parts, or faulty workmanship, which affect safety or result in substandard performance below specifications. R-Biopharm AG is not liable for any improper use nor the consequences of
 - a. the failure to read, understand and follow use and safety instructions, or use proper preparation and storage;
 - b. the failure to utilize trained and unqualified personnel, suitable samples or sampling techniques;
 - c. chemical, electromagnetic, mechanical, or electrolytic influences outside R-Biopharm AG provided standard perimeters through product specific data sheets, written product descriptions, or as otherwise expressly agreed upon in writing; or
 - d. any combination thereof.
3. R-Biopharm AG is also not liable for any changes or modifications, not carried out by R-Biopharm AG, nor consequences of use, applications, or processing which are unsafe or otherwise inconsistent with intended Product purpose as limited by written Product descriptions and specifications of R-Biopharm AG.
4. R-Biopharm AG’s liability for ordinary breach of contract is limited to repair, replacement, other substitute performance, or refund. The choice of remedies is within R-Biopharm AG’s sole discretion. R-Biopharm AG is not contractually liable for any incidental, or consequential damages, including but not limited to purchaser’s expenses, losses, or damages from loss of good will, frustrated business purposes, sales expectations, investment loss, actual or anticipated lost profits, End User indemnity, or any other business expenditures.

5. The foregoing limited warranty is solely intended to fulfill the warranty requirements (“Gewährleistung”) implied by German Civil Code. It is not intended to extend the scope or period of such Gewährleistung or provide additional warranties. Nor is this Disclaimer otherwise intended to limit or extent mandatory liability for damages in tort resulting from injury to life, body, or health, for intentional or grossly negligent acts, fraud, or due to strict liability under applicable product liability laws. A reversal of the burden of proof or rules of contract interpretation is not intended.