CompactDry EC

medium for *E. coli* and coliforms / Medium für *E. coli* und Coliforme / milieu pour *E. coli* et coliformes

100 plates / Platten / plaques
40 plates / Platten / plaques


Art. No. HS8781
Art. No. HS8782

CompactDry™

Simple and Easy Dry Media for Microbial Count and Detection

COMPACTDRY EC

Instructions

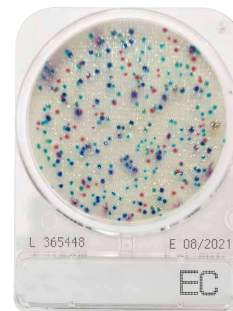Storage up
to 30 °C

Sample Preparation

Viable count in water or liquid foodstuff

1 mL of sample (diluted if necessary) is deposited on the CompactDry EC plate. Pre-treat the samples according to ISO 6887 or NMKL 91.

Viable count in solid food

Add buffer solution (e.g. 0.9 % saline solution, PBS, buffered peptone water, etc.) to the sample and homogenize using a Stomacher® (as per manufacturer instructions). 1 mL of sample (diluted if necessary) is dispensed in the centre on the CompactDry EC plate.

Viable count in swab test sample

Use the swab to wipe the surface, put the swab back into the tube and screw tightly. Shake the tube to mix. 1 mL of sample (diluted if necessary) is dispensed in the centre on the CompactDry EC plate.

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CompactDry Protocol

1 Open the cap. **2** Dispense 1 mL of sample in the middle of the CompactDry plate OR add filter as instructed. **3** Sample diffuses passively and evenly across the dehydrated media sheet transforming it into a gel within seconds. **4** Replace the cap, label and turn the plate over. **5** Incubate for the required time and temperature. **6** Following incubation, count the number of coloured colonies.

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Interpretation

E.coli forms blue colonies. Red and blue colonies together make up the total coliform group count.

General information

E.Coli 0157 forms pink/red/purple colonies.

The growth area is 20 cm². The back of the plate has a 1 cm² grid to ease colony counting. If there are a large number of colonies, the total viable count can be obtained by multiplying the average number of colonies in several grids by 20.

High concentrations (>250 CFU) will cause the entire growth area to become white/pink. In this case dilute the sample.

After use please follow local disposal regulations.

Storage and shelf life

Store at room temperature (+1 to +30 °C). Shelf life 24 months after manufacturing.

Inkubationstemperatur:
37 ±1 °C (Microval/NordVal)
35 ±2 °C (AOAC)
Inkubationszeit:
24 ±2 Stunden

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