COMPACTDRY ETB

Instructions

Storage up
to 30 °C

Sample Preparation

Viable count in water or liquid foodstuff

1 mL of sample (diluted if necessary) is deposited on the CompactDry ETB plate. Pre-treat the samples according to ISO 6887 or NMKL 91.

Viable count in solid food

Add buffer solution (e.g. 0.9 % saline solution, PBS, buffered peptone water, etc.) to the sample and homogenize using a Stomacher® (as per manufacturer instructions). 1 mL of sample (diluted if necessary) is dispensed in the centre on the CompactDry ETB plate.

Viable count in swab test specimen

Use the swab to wipe the surface, put the swab back into the tube and screw tightly. Shake the tube to mix. 1 mL of sample (diluted if necessary) is dispensed in the centre on the CompactDry ETB plate.

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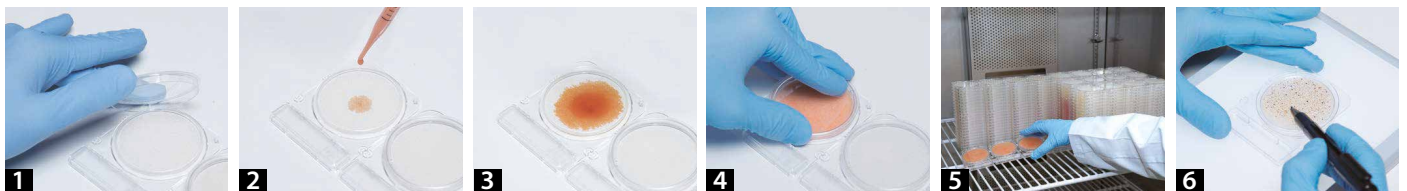
CompactDry Protocol

1 Open the cap. **2** Dispense 1 mL of sample in the middle of the CompactDry plate OR add filter as instructed. **3** Sample diffuses passively and evenly across the dehydrated media sheet transforming it into a gel within seconds. **4** Replace the cap, label and turn the plate over. **5** Incubate for the required time and temperature. **6** Following incubation, count the number of coloured colonies.

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Incubation time: 24 ±2 hours

Interpretation

Enterobacteriaceae are red/purple coloured. Red/purple coloured colonies plus other coloured colonies together form the total gram negative bacteria count.

General information

Some colonies might not be clearly red/pink coloured.

The growth area is 20 cm². The back of the plate has a 1 cm² grid to ease colony counting. If there are a large number of colonies, the total viable count can be obtained by multiplying the average number of colonies in several grids by 20.

High concentrations (>250 CFU) will cause the entire growth area to become red/pink. In this case dilute the sample.

After use please follow local disposal regulations.

Storage and shelf life

Store at room temperature (+1 to +30 °C). Shelf life 16 months after manufacturing.

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