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PolyBIND®

1. Characteristics

Surface activated particles of polyethylene

2. Intended use

PolyBIND® particles make a fast and easy isolation of microorganism from large sample amounts and viscous or solid loaded liquids possible. The statistical probability to determine trace contamination is radical increased with this product.

3. Form of delivery

Polymer powder with defined particle size (100-200 µm)

4. Preventive measures

Basically careful handling with chemicals. Do not inhale or swallow. Avoid skin or eye contact.

5. Content of package

The kit contains PolyBIND® powder for 50 analyses:
50 x 100 mg in glass bottles

6. Matrices

- beer, beer-based mixed drinks
- red wine
- clear and cloudy juices without pulp
- barrel soda
- mineral water

7. Principle

- PolyBIND® particles bind the given microorganisms in a liquid under acidic conditions (**pH-value <4.0**)
- separation of polymer particles with corresponding hardware:
 - A. with Granulab-19 Apparatus (s. www.rschuett.de)
 - B. with the inline proceed Granuinline (s. www.rschuett.de)
- in dependence of the cell count the PolyBIND® bonded microorganisms have to pre-concentrated in corresponding culture media or have to be directly analyzed
- microbiological diagnostic: visual control of turbidity, microscope or real-time PCR

8. Disposal

The pre-enriched culture medium with PolyBIND® powder has to be autoclaved after use.

Attention: PolyBIND® powder must not get into wastewater. therefore it has to be separated with a filter and disposed with residual waste.

Please mail to info@gen-ial.de for further questions about applications

GEN-IAL makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. If any materials are defective, GEN-IAL will provide a replacement product. There is no warranty of merchantability of this product, or of the fitness of the product for any purpose. GEN-IAL shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.