---

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## Enterobacteriaceae

### 1. Intended use

Qualitative PCR-detection of enterobacteriaceae as indicators for hygiene and postprocessing contaminations. The system contains an integrated inhibition control.

### 2. Test principle

The TaqMan<sup>®</sup> real-time PCR is based on hot-start-PCR and sequence-specific dual labelled probes (FAM/DQ) which, when accurately hybridised, emit a measurable fluorescent signal of a defined wavelength in the extension phase. The increase of signal is continuously measured in a real-time PCR detection instrument. To avoid false negative PCR-results an Inhibition Control (HEX/DQ) is amplified together in one reaction vessel with the specific sequence. The system contains dUTP. Optional: the use of Uracil-N-Glycosylase will eliminate any contamination with Uracil containing amplicons from former PCRs (the enzyme is not part of this kit).

### 3. Kit contents

The kit contains sufficient reagents for 50 PCR reactions:

1 x Premix	white cap
1 x Dye Mix (freeze-dried, incl. IC-DNA)	dark vial, red cap
1 x ddH <sub>2</sub> O	colourless cap
1 x Control-DNA (freeze-dried)	yellow cap

#### 4. Storage conditions

The Dye Mix and the Control-DNA are freeze-dried, they have to be solved in ddH<sub>2</sub>O prior to use (see 6.1).

Do **not** freeze the lyophilized Dye Mix and lyophilized Control-DNA.

The PCR reagents should be stored at 2 – 8 °C (35 – 46 °F).

Keep Premix for storage at - 20 °C (- 4 °F) **after arrival**. Avoid loss of sensitivity by repeating freezing and thawing more than 3 times. For irregular use aliquot the Premix.

The Dye Mix contains the fluorescent labelled probes and should be handled light protected. All reagents are stable for 12 months, if they are stored correctly.

#### 5. Materials required, but not provided

##### 5.1. Instruments

Real-time PCR instrument with channels FAM and HEX

Centrifuge for 1.5 – 2.0 mL reaction tubes

Centrifuge for multiwell plates or strips

Pipettes

“Vortex”

##### 5.2. Reagents and plastic ware

sterile ddH<sub>2</sub>O

sterile reaction vessels 1.5 – 2.0 mL

sterile optical tubes, plates, capillaries

sterile filter tips

**optional:** Uracil N-Glycosylase (0.01 U/μL added to the PCR reaction mix)

## 6. PCR

### 6.1. PCR-setup

*When using the kit for the first time, the freeze-dried kit components have to be shortly centrifuged and carefully resolved:*

- add 80  $\mu\text{L}$  sterile ddH<sub>2</sub>O to freeze dried Dye Mix
- add 55  $\mu\text{L}$  sterile ddH<sub>2</sub>O to the freeze-dried Control-DNA
- after 15 minutes mix well

**Before every use thoroughly mix all PCR-components and centrifuge briefly.**

PCR-reaction per sample:

PCR-Components	amount ( $\mu\text{L}$ )
Premix	16.0
Dye Mix	1.5
Sample-DNA	up to 2.5
Total volume	<b>20.0</b>

1. Prepare a mastermix by mixing Premix and Dye Mix
2. Multiply said volumes with the number of PCR preparations including controls (positive control, negative control, extraction control), taking into account pipette reserves of approximately 5-10 %.
3. Divide 17.5  $\mu\text{L}$  of the PCR-mastermix among the individual reaction vessels, making sure that, prior to the first filling, the tip of the pipette has been moistened.
4. Add 2.5  $\mu\text{L}$  sample DNA, add 2.5  $\mu\text{L}$  of the Control-DNA for the PCR positive control, add 2.5  $\mu\text{L}$  of the extraction control and 2.5  $\mu\text{L}$  of ddH<sub>2</sub>O for the negative control\* reaction. Use a fresh tip with each DNA filling.
5. Close the tubes immediately and centrifuge them shortly.
6. Place the tubes in the PCR-machine and start run.

**Very important: \* Please fill up the negative control with 2.5  $\mu\text{L}$  ddH<sub>2</sub>O to avoid unspecific amplification.**

**Work swiftly to avoid warming up and keep away from light.**

## 6.2 PCR-Program

### 6.2.1 PCR-Program LC480

1. Click the button **tool** on the right side in the window **LightCycler 480 Software release 1.5.0. SP1**
2. Click the button **Detection formats** at the left side of the menu bar
3. Click **New** in the window **Detection formats** and name the experiment
4. Open the window **Filter Combination Selection** and choose the following filter combinations: 465-510 / 533-580
5. Open the window **Selected Filter Combination List** and add the following amounts

Excitation Filter	Emission Filter	Name	Melt factor	Quant factor	Max. Integ. Time
465	510	465-510 FAM	1	10	2
533	580	533-580 HEX	1	10	2

6. **Close** the window
7. Click the button **New Experiment** on the right side of the menu bar
8. From the pull-down menu **Detection formats** choose the defined experiment, click the button **Customize** and check the detection formats. All of them have to be activated.
9. Click the button **ok**
10. Define the following program:

1. Program Name: **Heat**

Cycles **1** Analysis Mode **None**

Target (°C)	Acquisition Mode	Hold (hh:mm:ss)	Ramp Rate (°C/s)	Acquisitions (per °C)	Sec Target (°C)	Step size (°C)	Step Delay (cycles)
95	None	00:15:00	4.40		0	0	0

2. Program Name: **Ampli**

Cycles **40** Analysis Mode **Quantification**

Target (°C)	Acquisition Mode	Hold (hh:mm:ss)	Ramp Rate (°C/s)	Acquisitions (per °C)	Sec Target (°C)	Step size (°C)	Step Delay (cycles)
95	None	00:00:10	4.40		0	0	0
60	Single	00:00:20	2.20		0	0	0

3. Program Name: **Cool**

Cycles **1** Analysis Mode **None**

Target (°C)	Acquisition Mode	Hold (hh:mm:ss)	Ramp Rate (°C/s)	Acquisitions (per °C)	Sec Target (°C)	Step size (°C)	Step Delay (cycles)
40	None	00:00:20	2.20		0	0	0

**Optional:** Saving the program as **run template**. The template button allows to select and apply a template to the currently open object and to save the currently open object as a template. Click the clamp beside the button **Apply Template**. Click **Save as template** and save the file in **templates**.

11. Click the button **Subset editor** on the left side

12. Click the button **+** and **New Subset 1** appears

13. Mark the wells in the **New Subset 1 Settings** window

14. Click the button **Apply**

15. Click the button **Sample editor** on the left side

16. **Very important:** Activate In the window **Step 1 Select Workflow: Abs.Quant**

17. Choose the subset **New Subset** in the window **Step 2 Select Samples**

18. Define your probes in the **Sample table**

19. Switch to the button **Experiment** and start the run with the button **start run**

## 6.2.2 PCR-Program for other real-time machines

The Premix contains no ROX. This must be considered according to the settings of the real-time machine. **ABI 7500:** Under "Assign Targets and Samples" in `Select the dye to use as the passive reference` choose „none“.

**Mx3500P:** Under Instruments in `Filter Set Gain Settings` change the Filter set gain multiplier to CY5 x1, ROX x1, HEX x4 und FAM x8.

For the use of UNG the thermal cycler program has to be changed according to manufacturers instructions.

Step	Time	Temp.	
Initial denaturation of DNA	15 min	95 °C	
Cycling Denaturation	10 sec	95 °C	Cycle 40 x
Cycling Annealing/ Elongation	20 - 30 sec*	60 °C	

\* please program 20 sec, if the real-time-PCR machine allows it

## For MyGo Pro User

### Software 3.5 / 3.6 MyGo Pro and ESR

Step	Time	Temp.	
Initial denaturation of DNA	15 min	95 °C	
Cycling Denaturation	10 sec	95 °C	Cycle 40 x
Cycling Annealing/ Elongation	30 sec	60 °C	

**Advanced Settings**

Integration time (s)

0.9

## 7. Evaluation

The evaluation has to be done according to the data analysis program recommended by the real-time instrument manufacturer.

For LC480: For analysis activate Colour Compensation

**Enterobacteriaceae-DNA:** FAM-channel (LC480: 465-510 nm)

**Inhibition Control-DNA:** HEX-channel (LC480: 533-580 nm)


A sample is **Enterobacteriaceae positive**, if there is a detectable fluorescence increase in the **FAM-channel (465-510 nm)** and the negative controls show no amplification. The positive controls should have a positive fluorescence signal. The Inhibition Control in the HEX-channel (533-580 nm) may be positive or negative (depending on the amount of DNA or inhibitors in the sample reaction). For negative controls it has to be positive.

A sample is negative, if there is no detectable fluorescence increase in the different channels and the positive controls have a positive fluorescence signal. The negative controls show no amplification. The Inhibition Control in the HEX-channel (533-580 nm) has to be positive in the sample and in the negative controls, a false negative result due to inhibitory effects is then excluded.

# analysis flowchart

## LC480: analysis after activating Colour Compensation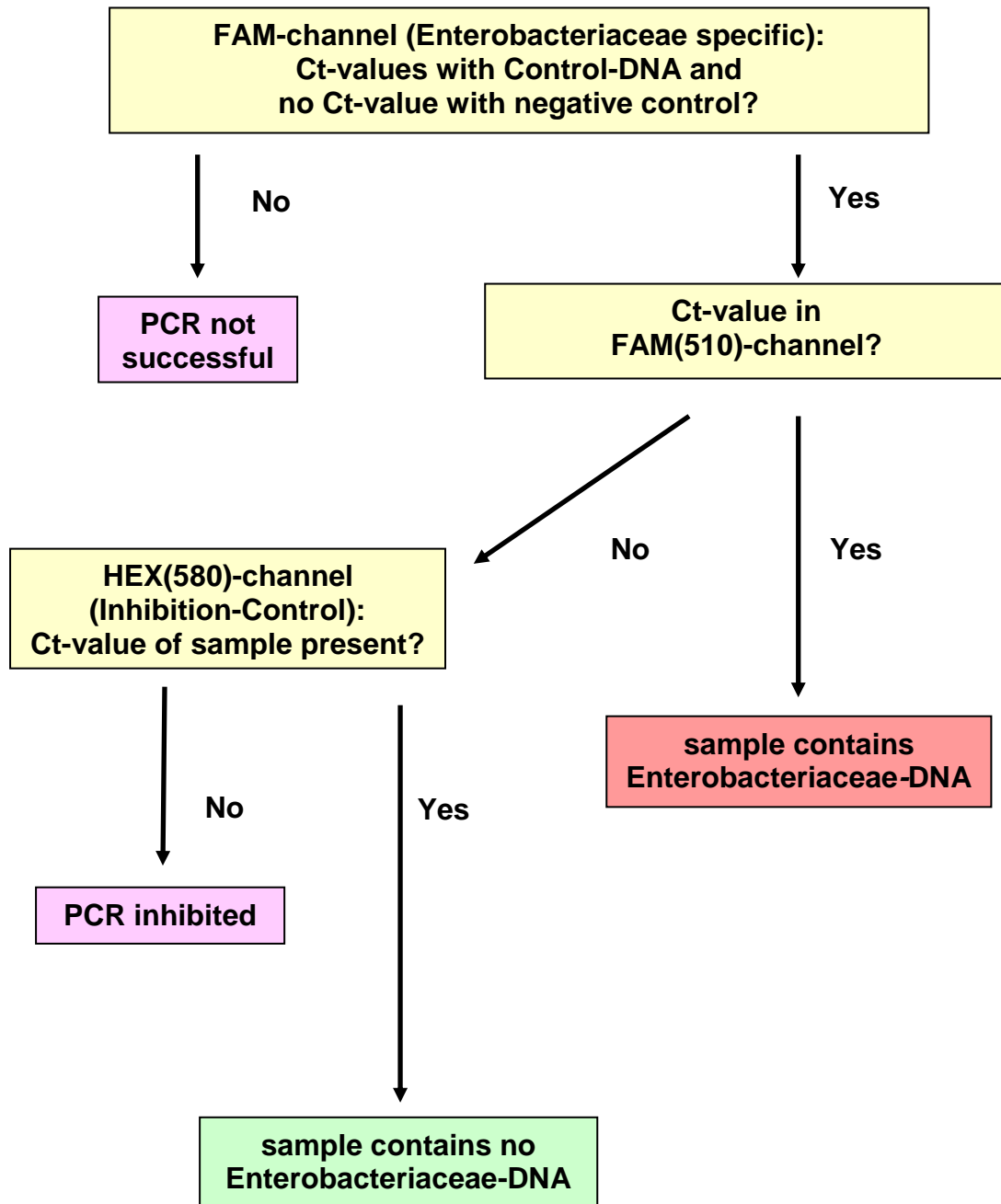

**Note:**

The polymerase-chain reaction (PCR) is protected by patents and requires a licence from Hoffmann-LaRoche Inc..

The provided product does not authorise the purchaser for the commercial use of this method.

GEN-IAL makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. If any materials are defective, GEN-IAL will provide a replacement product. There is no warranty of merchantability of this product, or of the fitness of the product for any purpose. GEN-IAL shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.