---

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## Screening and differentiation

### 1. Intended use

Screening and differentiation of beer spoilage bacteria in beer and beer mixing drinks. The following bacteria are differentiated:

Well	FAM	JOE	ROX
1	Negative control		Inhibition control
2	Positive control		
3	<i>L. rossiae</i>	<i>L. backii</i>	Inhibition control
4	<i>L. brevis/parabrevis/brevisimilis</i>	<i>L. lindneri</i>	<i>L. casei/paracasei</i>
5	Pediococcus spp.**	Lactobacillus spp.*	Pectinatus./ Megasphaera

\* **Lactobacillus spp.:** *L. acetotolerans*, *L. collinoides/ paracollinoides*, *L. coryniformis*, *L. plantarum/ paraplantarum*, *L. perolens/ harbinensis*, *L. buchneri/ parabuchneri*

\*\* **Pediococcus spp.** *P. acidilactici*, *P. parvulus*, *P. inopinatus*, *P. pentosaceus*, *P. damnosus*, *P. clausenii*

### 2. Test principle

The TaqMan® real-time PCR is based on hot-start-PCR and sequence-specific dual labelled probes (FAM/DQ; JOE/DQ; ROX/DQ) which, when accurately hybridised, emit a measurable fluorescent signal of a defined wavelength in the extension phase. The increase of signal is continuously measured in a real-time PCR detection instrument. To avoid false negative PCR-results an Inhibition control (ROX/DQ) is amplified together in one tube with the specific sequence. The system contains dUTP. Optional: the use of Uracil-N-Glycosylase will eliminate any contamination with Uracil containing amplicons from former PCRs (the enzyme is not part of this kit).

### 3. Kit contents

The kit contains sufficient reagents for 96 PCR reactions (24 samples):

1 x Premix	white cap
12 x Dye Strips (freeze-dried)	tube strips
1 x ddH <sub>2</sub> O (for PCR negative controls)	colourless cap

#### 4. Storage conditions

Do **not** freeze the lyophilized Dye Strips.

The PCR reagents should be stored at 2 – 8 °C (35 – 46 °F).

Keep Premix for storage at - 20 °C (- 4 °F) **after arrival**. Avoid loss of sensitivity by repeating freezing and thawing more than 3 times. For irregular use aliquot the Premix.

The Dye Strips contain the fluorescent labelled probes and should be handled light protected.

All reagents are stable for 12 months, if they are stored correctly.

#### 5. Materials required, but not provided

##### 5.1. Instruments

Real-time PCR machine for low profile tubes clear

Centrifuge for tube strips

Pipettes

“Vortex”

##### 5.2. Reagents and plastic ware

sterile ddH<sub>2</sub>O

sterile filter tips

**optional:** Uracil N-Glycosylase (0.01 U/μL added to the PCR reaction mix)

## 6. PCR

### 6.1. PCR-Setup

Before every use thoroughly mix Premix and centrifuge briefly.

PCR-reaction setup:

PCR-Components	Volume (µL)
Premix	15.0
Sample-DNA	5.0*
Total volume	<b>20.0</b>

\* if using the Simplex Easy DNA- or QuickGEN Yeast Sample Preparation Kit:  
add 2.5 µL DNA and 2.5 µL sterile ddH<sub>2</sub>O

Pipetting scheme:

Tube 1*	Premix 15.0 µL	<b>NTC, ddH<sub>2</sub>O 5.0 µL,</b> <b><u>Attention:</u> pipette no sample-DNA</b>
Tube 2*	Premix 15.0 µL	5,0 µL ddH <sub>2</sub> O, PTC already included in the tube
Tube 3*	Premix 15.0 µL	sample-DNA 5.0 µL (sample 1)
Tube 4*	Premix 15.0 µL	sample-DNA 5.0 µL (sample 1)
Tube 5*	Premix 15.0 µL	sample-DNA 5.0 µL (sample 1)
Tube 6*	Premix 15.0 µL	sample-DNA 5.0 µL (sample 2)
Tube 7*	Premix 15.0 µL	sample-DNA 5.0 µL (sample 2)
Tube 8*	Premix 15.0 µL	sample-DNA 5.0 µL (sample 2)

\* please notice the orientation of the tubes: they have to be opened to the right side

## 6.2 PCR-Program

**Note:** For the use of UNG the thermal cycler program has to be changed according to manufacturers` instructions.

### Software 3.5 / 3.6 ESR

Step	Time	Temp.	
Initial denaturation	15 min	95 °C	
Cycling Denaturation	15 sec	95 °C	Cycle 35 x
Cycling Annealing/ Elongation	30 sec	60 °C	

#### Advanced Settings

Integration time (s)

0.7

## 7. Evaluation

The evaluation has to be made according to the data analysis program recommended by the real-time instrument manufacturer.

The negative control (NTC, tube 1) has to be negative in the FAM- and JOE-channel and positive in the ROX-channel (Inhibition Control). Is the Inhibition Control in the ROX-channel negative and/or the results in the FAM/JOE-channel positive, PCR has to be repeated.

The Inhibition Control (tube 3 and 6) has to be positive in the ROX-channel. Is the Inhibition Control in the ROX-channel negative, inhibitors or high amount of DNA are in the sample reaction.

The Positive Control (tube 2) has to be positive in the FAM-channel. If it is negative, PCR has to be repeated.

## PCR-Analysis:

please notice the orientation of the tubes: they have to be opened to the right side

	Channel	Species	Signal
<b>Tube 1</b>	FAM	NTC	negative
	JOE	NTC	negative
	ROX	Inhibition Control	positive
<b>Tube 2</b>	FAM	Positive Control	positive
	JOE	-	
	ROX	-	
<b>Tube 3</b>	FAM	<i>Lactobacillus rossiae</i>	
	JOE	<i>Lactobacillus backii</i>	
	ROX	Inhibitionskontrolle	positive
<b>Tube 4</b>	FAM	<i>Lactobacillus brevis/ parabrevis/ brevisimilis</i>	
	JOE	<i>Lactobacillus lindneri</i>	
	ROX	<i>Lactobacillus casei/ paracasei</i>	
<b>Tube 5</b>	FAM	Pediococcus spp.	
	JOE	Lactobacillus spp.	
	ROX	Pectinatus spp./ Megasphaera spp.	
<b>Tube 6</b>	FAM	<i>Lactobacillus rossiae</i>	
	JOE	<i>Lactobacillus backii</i>	
	ROX	Inhibition Control	positive
<b>Tube 7</b>	FAM	<i>Lactobacillus brevis/ parabrevis/ brevisimilis</i>	
	JOE	<i>Lactobacillus lindneri</i>	
	ROX	<i>Lactobacillus casei/ paracasei</i>	
<b>Tube 8</b>	FAM	Pediococcus spp.	
	JOE	Lactobacillus spp.	
	ROX	Pectinatus spp./ Megasphaera spp.	

**Example for interpretation:** If the samples show positive JOE-signals in tubes 4 and/or 7 *Lactobacillus lindneri* is identified in the samples. If the sample contains mixtures, you have positive signals in different channels.

Note:

The polymerase-chain reaction (PCR) is protected by patents and requires a licence from Hoffmann-LaRoche Inc.. The provided product does not authorise the purchaser for the commercial use of this method. GEN-IAL makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. If any materials are defective, GEN-IAL will provide a replacement product. There is no warranty of merchantability of this product, or of the fitness of the product for any purpose. GEN-IAL shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.