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Pipettierschema:

Tube 1*	Premix 15,0 µL	ddH₂O 5,0 µL, NTC <u>Achtung:</u> keine Proben-DNA pipettieren
Tube 2*	Premix 15,0 µL	ddH₂O 5,0 µL, PTC bereits im tube enthalten <u>Achtung:</u> keine Proben-DNA pipettieren
Tube 3*	Premix 15,0 µL	Proben-DNA 5,0 µL (Probe 1)
Tube 4*	Premix 15,0 µL	Proben-DNA 5,0 µL (Probe 1)
Tube 5*	Premix 15,0 µL	Proben-DNA 5,0 µL (Probe 1)
Tube 6*	Premix 15,0 µL	Proben-DNA 5,0 µL (Probe 2)
Tube 7*	Premix 15,0 µL	Proben-DNA 5,0 µL (Probe 2)
Tube 8*	Premix 15,0 µL	Proben-DNA 5,0 µL (Probe 2)

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Advanced Settings MyGo Pro	Integration time (s)	0.7
Advanced Settings MyGo Pro ESR	Integration time (s)	1.0

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Screening and differentiation

1. Intended use

Screening and differentiation of beer spoilage bacteria in beer and beer mixing drinks. The following bacteria are differentiated:

Well	FAM	HEX	ROX
1	Negative control		Inhibition control
2	Positive control		
3	<i>L. rossiae</i>	<i>L. backii</i>	Inhibition control
4	<i>L. brevis/parabrevis/brevisimilis</i>	<i>L. lindneri</i>	<i>L. casei/paracasei</i>
5	Pediococcus spp.**	Lactobacillus spp.*	Pectinatus./ Megasphaera

* **Lactobacillus spp.:** *L. acetotolerans*, *L. collinoides/ paracollinoides*, *L. coryniformis*, *L. plantarum/ paraplantarum*, *L. perolens/ harbinensis*, *L. buchneri/ parabuchneri*

** **Pediococcus spp.** *P. acidilactici*, *P. parvulus*, *P. inopinatus*, *P. pentosaceus*, *P. damnosus*, *P. claussenii*

2. Test principle

The TaqMan® real-time PCR is based on hot-start-PCR and sequence-specific dual labelled probes (FAM/DQ; HEX/DQ; ROX/DQ) which, when accurately hybridised, emit a measurable fluorescent signal of a defined wavelength in the extension phase. The increase of signal is continuously measured in a real-time PCR detection instrument. To avoid false negative PCR-results an Inhibition control (ROX/DQ) is amplified together in one tube with the specific sequence.

3. Kit contents

The kit contains sufficient reagents for 96 PCR reactions (24 samples):

1 x Premix	white cap
12 x Dye Strips (freeze-dried)	tube strips
1 x ddH ₂ O (for PCR negative controls)	colourless cap

4. Storage conditions

Do **not** freeze the lyophilized Dye Strips.

The PCR reagents should be stored at 2 – 8 °C (35 – 46 °F).

Keep Premix for storage at - 20 °C (- 4 °F) **after arrival**. Avoid loss of sensitivity by repeating freezing and thawing more than 3 times. For irregular use aliquot the Premix.

The Dye Strips contain the fluorescent labelled probes and should be handled light protected.

All reagents are stable for 12 months, if they are stored correctly.

5. Materials required, but not provided

5.1. Instruments

Real-time PCR machine for low profile tubes clear

Centrifuge for tube strips

Pipettes

“Vortex”

5.2. Reagents and plastic ware

sterile ddH₂O

sterile filter tips

6. PCR

6.1. PCR-Setup

Before every use thoroughly mix Premix and centrifuge briefly.

PCR-reaction setup:

PCR-Components	Volume (µL)
Premix	15.0
Sample-DNA	5.0*
Total volume	20.0

* if using the Simplex Easy DNA- or QuickGEN Yeast Sample Preparation Kit:
add 2.5 µL DNA and 2.5 µL sterile ddH₂O

Pipetting scheme:

Tube 1*	Premix 15.0 µL	ddH₂O 5.0 µL, NTC <u>Attention:</u> pipette no sample-DNA
Tube 2*	Premix 15.0 µL	ddH₂O 5.0 µL, PTC already included in the tube <u>Attention:</u> pipette no sample-DNA
Tube 3*	Premix 15.0 µL	sample-DNA 5.0 µL (sample 1)
Tube 4*	Premix 15.0 µL	sample-DNA 5.0 µL (sample 1)
Tube 5*	Premix 15.0 µL	sample-DNA 5.0 µL (sample 1)
Tube 6*	Premix 15.0 µL	sample-DNA 5.0 µL (sample 2)
Tube 7*	Premix 15.0 µL	sample-DNA 5.0 µL (sample 2)
Tube 8*	Premix 15.0 µL	sample-DNA 5.0 µL (sample 2)

* please notice the orientation of the tubes: they are marked with A - H

6.2 PCR-Program

Software MyGo Pro/ ESR 3.5 / 3.6

Step	Time	Temp.	
Initial denaturation	15 min	95 °C	
Cycling Denaturation	15 sec	95 °C	Cycle 35 x
Cycling Annealing/ Elongation	30 sec	60 °C	

Advanced Settings MyGo Pro Integration time (s) 0.7

Advanced Settings MyGo Pro ESR Integration time (s) 1.0

7. Evaluation

The evaluation has to be made according to the data analysis program recommended by the real-time instrument manufacturer.

The negative control (NTC, tube 1) has to be negative in the FAM- and HEX-channel and positive in the ROX-channel (Inhibition Control). Is the Inhibition Control in the ROX-channel negative and/or the results in the FAM/HEX-channel positive, PCR has to be repeated.

The Inhibition Control (tube 3 and 6) has to be positive in the ROX-channel. Is the Inhibition Control in the ROX-channel negative, inhibitors or high amount of DNA are in the sample reaction.

The Positive Control (tube 2) has to be positive in the FAM-channel. If it is negative, PCR has to be repeated.

PCR-Analysis:

please notice the orientation of the tubes: they are marked with A - H

	Channel	Species	Signal
Tube 1	FAM	NTC	negative
	HEX	NTC	negative
	ROX	Inhibition Control	positive
Tube 2	FAM	Positive Control	positive
	HEX	-	
	ROX	-	
Tube 3	FAM	<i>Lactobacillus rossiae</i>	
	HEX	<i>Lactobacillus backii</i>	
	ROX	Inhibitionskontrolle	positive
Tube 4	FAM	<i>Lactobacillus brevis/ parabrevis/ brevisimilis</i>	
	HEX	<i>Lactobacillus lindneri</i>	
	ROX	<i>Lactobacillus casei/ paracasei</i>	
Tube 5	FAM	Pediococcus spp.	
	HEX	Lactobacillus spp.	
	ROX	Pectinatus spp./ Megasphaera spp.	
Tube 6	FAM	<i>Lactobacillus rossiae</i>	
	HEX	<i>Lactobacillus backii</i>	
	ROX	Inhibition Control	positive
Tube 7	FAM	<i>Lactobacillus brevis/ parabrevis/ brevisimilis</i>	
	HEX	<i>Lactobacillus lindneri</i>	
	ROX	<i>Lactobacillus casei/ paracasei</i>	
Tube 8	FAM	Pediococcus spp.	
	HEX	Lactobacillus spp.	
	ROX	Pectinatus spp./ Megasphaera spp.	

Example for interpretation: If the samples show positive HEX-signals in tubes 4 and/or 7 *Lactobacillus lindneri* is identified in the samples. If the sample contains mixtures, you have positive signals in different channels.

Note:

The polymerase-chain reaction (PCR) is protected by patents and requires a licence from Hoffmann-LaRoche Inc.. The provided product does not authorise the purchaser for the commercial use of this method. GEN-IAL makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. If any materials are defective, GEN-IAL will provide a replacement product. There is no warranty of merchantability of this product, or of the fitness of the product for any purpose. GEN-IAL shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.