r-biopharm®

RIDASCREEN® Nitrofuran AHD

REF R3713

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RIDASCREEN® Nitrofurantoin (AHD)

Brief information

RIDASCREEN® Nitrofurantoin (AHD) (Art. No. 3713) is a competitive enzyme immunoassay for the quantitative analysis of the nitrofurantoin metabolite AHD in shrimp and fish (see chapter 1. Intended use).

All reagents required for the enzyme immunoassay, including standards, are contained in the test kit. The test kit is sufficient for a maximum of 96 determinations (including standards). A microtiter plate spectrophotometer is required for quantification.

Sample preparation:	homogenization, derivatization, extraction, centrifugation, evaporation and defatting
Time requirement:	Sample preparation for 10 samples Sample preparation part 1.....approx. 1.5 h Incubation.....3 h Sample preparation part 2approx. 1.5 h Test implementation (incubation time) .. 1 h 15 min
Detection limit: (corresponding to the standard substance)	Shrimp approx. 200 ng/kg (ppt) Fish..... approx. 76 ng/kg (ppt)
Detection capability (cc β): (corresponding to the standard substance)	Shrimp 250 ng/kg
Recovery rate: (corresponding to the standard substance)	Shrimp 94 % Fish..... 82 %
Specificity:	Nitrophenyl- (NP) AHD 100 % NP-AMAZ, NP-AOZ, NP-SEM < 1 % Nitrofurantoin approx. 32 % Furaltadone, furazolidone, nitrofurazone..... < 2 % AHD, AMAZ, AOZ, SEM..... < 1 %

The specificity of the RIDASCREEN® Nitrofurantoin (AHD) test was determined by analyzing the cross reactivities to corresponding substances in buffer system. In samples, the specificity may deviate from those determined in the buffer system due to matrix effects. Prior to the analysis of cross-reactive substances, the user has to determine the Limit of Detection and the Recovery for the substance in the respective sample matrix. The test cannot discriminate between analytes and cross-reactive substances.

In order to increase the quality of assessment when performing ELISA procedures, we refer additionally to our Good ELISA Practice brochure. It lists minimum standards and conditions that are required when using test kits of R-Biopharm AG to perform ELISA analysis. The brochure can be retrieved, printed and downloaded from the website

<https://food.r-biopharm.com/media/technical-guides/>.

Related products and accessories for nitrofurantoin determination

RIDASCREEN® Nitrofurantoin (AOZ) (Art. No. R3703)

RIDASCREEN® Nitrofurantoin (AMTZ) (Art. No. R3722)

RIDASCREEN® Nitrofurantoin (SEM) (Art. No. R3724)

RIDA® Nitrofurantoin (AMTZ) Spiking Solution (Art. No. R3799)

RIDA® Nitrofurantoin (AOZ) Spiking Solution (Art. No. R3798)

RIDA® Nitrofurantoin (AHD) Spiking Solution (Art. No. R3796)

RIDA® Nitrofurantoin (SEM) Spiking Solution (Art. No. R3797)

1. Intended use

RIDASCREEN® Nitrofurantoin (AHD) (Art. No. 3713) is a competitive enzyme immunoassay for the quantitative analysis of the nitrofurantoin metabolite AHD in shrimp and fish.

2. General information

Nitrofurans are synthetic broad-spectrum antibiotics, which are frequently used in animal production due to their excellent antibacterial and pharmacokinetic properties. They have also been used as growth promoters during the production of Shrimp, poultry and pigs. Long term animal experiments have shown that the parent compounds and their metabolites have carcinogenic and mutagenic characteristics. This led to the prohibition of nitrofurans for the treatment of animals used for food production. In 1993, the EU banned the

nitrofurans furaltadone, nitrofurantoin and nitrofurazone for use in animals used as sources of food, and in 1995 the use of furazolidone was also prohibited.

The analysis of nitrofurans are based on the detection of the tissue bound metabolites of nitrofurans. The parent compounds are difficult to detect accurately since they are metabolized very rapidly after treatment. The tissue bound nitrofuran metabolites however are present for a long time after administration and they are used to detect nitrofuran abuse. Prior to analysis, the metabolites have to be derivatized by incubation with 2-Nitrobenzaldehyde (2NBA) into Nitrophenyl- (NP) AHD, NP-AMOZ, NP-AOZ and NP-SEM.

parent compound	metabolite		after derivatization
Nitrofurantoin	1-Aminohydantoin	AHD	NP-AHD
Furaltadon	3-Amino-5-morpholino methyl-2-oxazolidinon	AMOZ	NP-AMOZ
Furazolidon	3-Amino-2-oxazolidinon	AOZ	NP-AOZ
Nitrofurazon	Semicarbazide	SEM	NP-SEM

3. Test principle

The basis of the test is the antigen-antibody reaction. The microtiter wells are coated with capture antibodies directed against anti-AHD antibodies. AHD standards or sample solutions, AHD conjugate and anti-AHD antibodies are added. Free AHD and AHD conjugate compete for the anti-AHD antibody binding sites (competitive enzyme immunoassay). At the same time, the anti-AHD antibodies are also bound by the immobilized capture antibodies. Any unbound conjugate is then removed in a washing step. Substrate/chromogen is added to the wells and incubated. Bound conjugate converts the chromogen into a blue product. The addition of the stop solution leads to a color change from blue to yellow. The measurement is made photometrically at 450 nm. The absorption is inversely proportional to the AHD concentration in the sample.

4. Reagents provided

Each kit contains sufficient materials for a maximum of 96 measurements (including standard analyses). Each test kit contains:

Component	Cap Colour	Format	Volume
Microtiter plate	-	Ready to use	96 wells
2-Nitrobenzandehyd	-	solid	100 mg
Standard 1	White	Ready to use	0 ng/L 1.3 mL
Standard 2	White	Ready to use	100 ng/L 1.3 mL
Standard 3	White	Ready to use	300 ng/L 1.3 mL
Standard 4	White	Ready to use	900 ng/L 1.3 mL
Standard 5	White	Ready to use	2700 ng/L 1.3 mL
Standard 6	White	Ready to use	8100 ng/L 1.3 mL
Wash buffer salt Tween	-	Salt for dissolving	
Conjugate	Red	Concentrate	11x 0.7 mL
Antibody	Black	Concentrate	11x 0.7 mL
Substrate/Chromogen Substrat/Chromogen Red Chromogen Pro	Brown	Ready to use	10 mL
Stop solution	Yellow	Ready to use	14 mL

5. Reagents required but not provided

5.1 Equipment

- Gloves
- Scale
- Centrifuge + centrifugal vials with cap (e.g. 50 mL centrifuge tubes from Greiner Art. No. 227261)
- Shaker (Vortex)
- Mixer (Stomacher, Ultraturrax)
- Rotary evaporator
- Magnetic stirrer
- Pasteur pipettes
- Graduated pipettes
- Variable 20 - 200 µL and 200 - 1000 µL micropipettes
- If necessary: Multistepper or 8-channel pipette for 50 and 100 µL
- Microtiter plate spectrophotometer (450 nm)
- Optional: RIDASOFT® Win.NET (Art. No. Z9996FF)

5.2 Reagents

- Distilled water (dist. water) or deionized water
- 1 M HCl
- 1 M NaOH
- 0.1 M K₂HPO₄
- Ethyl acetate
- n-Hexane
- Dimethylsulfoxide (DMSO)

6. Warnings and precautions for the users

This test should only be carried out by trained laboratory personnel. The instruction for use must be strictly followed.

This kit may contain hazardous substances. For hazard notes on the contained substances please refer to the appropriate material safety data sheets (SDS) for this product, available online at www.r-biopharm.com.

Do not reuse wells of the microtiter strips (coated microtiter plate, see chapter 10.2.). Use separate pipette tips for each standard and each sample extract to avoid cross contamination.

All reagents and materials must be recovered or disposed after use at customers own responsibility according to the protection of human health and the environment. Please observe the applicable national regulations concerning waste disposal (e.g. Waste Management Act, Regulations on Dangerous Chemicals, etc.).

7. Storage instructions

Store the kit at 2 - 8 °C (35 - 46 °F). Do not freeze any test kit components.

To avoid moisture inside the wells, open the foil bag for withdrawal of microwells only after having reached room temperature (20 - 25 °C / 68 - 77 °F).

Return any unused microwells to their original foil bag, reseal them together with the desiccant provided and further store at 2 - 8 °C (35 - 46 °F).

The reddish substrate/chromogen is light sensitive. Therefore, avoid exposure to direct light.

Do not use the test kit after the expiration date (see test kit label).

Do not interchange individual reagents between kits of different lot numbers.

8. Indication of instability or deterioration of reagents

- Bluish coloration of the reddish substrate/chromogen prior to test implementation
- Value of less than 0.6 absorbance units ($A_{450\text{ nm}} < 0.6$) for standard 1

9. Sample preparation

The samples should be stored in a cool place, protected against light.

- Homogenize the sample completely, use stomacher or mixer
- Prepare 10 mM 2-nitrobenzaldehyde (see chapter 4.) in dimethyl sulfoxide (DMSO)
(Note: this solution has to be prepared directly before use, i.e. dissolve 7.6 mg 2-Nitrobenzaldehyde in 5 mL DMSO.)

9.1. Shrimp

- Mix 1 g of the homogenized sample with 4 mL demineralized water, 0.5 mL 1 M HCl and 100 μ L 10 mM 2-nitrobenzaldehyde (in DMSO) by shaking vigorously
- Derivatize as described in chapter 9.3

9.2. Fish

- Mix 1 g of the homogenized sample with 3.9 mL demineralized water, 0.5 mL 1 M HCl and 200 μ L 10 mM 2-Nitrobenzaldehyde (in DMSO) by shaking vigorously
- Derivatize as described in chapter 9.3

9.3. Derivatization

- Incubate at 50 °C (122 °F) for 3 hours
- Add 5 mL 0.1 M K_2HPO_4 , 0.4 mL 1 M NaOH and 5 mL ethyl acetate, shake vigorously for 30 s
- Optional: sample incubation up to 50 °C (122 °F) can be used for better phase separation
- Centrifuge: 10 min / 3000 g / room temperature (20 - 25 °C / 68 - 77 °F)
- Transfer 2.5 mL of the ethyl acetate layer (upper layer) into a new vial and evaporate
- Dissolve the residue in 1 mL n-hexane and mix properly with 1 mL sample buffer (see chapter 10.1)
- Centrifuge: 10 min / 3000 g / room temperature (20 - 25 °C / 68 - 77 °F)
- Use 50 μ L of the lower, aqueous phase per well in the assay

10. Test procedure

10.1 Test preparation

Bring all reagents to room temperature (20 - 25 °C / 68 - 77 °F) before use.

The **anti-AHD antibody** is provided as a concentrate. Since the diluted antibody solution has a limited stability, only the amount which actually is needed should be reconstituted. Before pipetting, the antibody concentrate should be shaken thoroughly. For reconstitution, the antibody concentrate is diluted 1:11 (1+10) in wash buffer (e. g. 200 µL concentrate + 2 mL wash buffer, ready to use sufficient for 4 microtiter strips).

The **AHD conjugate** is provided as a concentrate. Since the diluted conjugate solution has a limited stability, only the amount which actually is needed should be reconstituted. Before pipetting, the conjugate concentrate should be shaken carefully. For reconstitution, the conjugate concentrate is diluted 1:11 (1+10) in wash buffer (e. g. 200 µL concentrate + 2 mL wash buffer, ready to use sufficient for 4 microtiter strips).

A PBS-Tween buffer is needed as **wash, sample, antibody and conjugate buffer**, please use the wash buffer salt (pouch) contained in the kit (see chapter 4.). Dissolve the total content of the pouch in one liter of demineralized water. The ready to use wash buffer expires after approx. 4 weeks at 2 - 8 °C (36 - 46 °F).

Alternative: Dissolve the contents of the pouch in 100 mL of demineralized water to obtain a 10-fold concentrated wash buffer. Dilute the 10-fold concentrated wash buffer 1:10 (1+9) with demineralized water to obtain the ready to use wash buffer. This solution expires after approx. 8 weeks, stored at room temperature (20 - 25 °C / 68 - 77 °F).

Components should be stored immediately at 2 - 8 °C (35 - 46 °F) when no longer required.

10.2 Test procedure

Carefully follow the recommended washing procedure to obtain unambiguous results. Do not allow microwells to dry between work steps.

It is recommended to pipette the conjugate, the antibody, the substrate/chromogen and the stop solution with a multi-channel or stepper pipette to avoid a time shift over the plate.

Avoid direct sunlight during all incubations. Therefore cover the microtiter plates.

1. Insert a sufficient number of wells into the microwell holder for all standards and samples to be run in duplicate. Record standard and sample positions.
2. Add 50 µL of each standard or prepared sample to separate duplicate wells. Use a new pipette tip for each standard or sample.
3. Add 50 µL of the diluted conjugate to the bottom of each well.
4. Add 50 µL of the diluted antibody to each well, mix gently by shaking the plate manually and incubate for 1 h at room temperature (20 - 25 °C / 68 - 77 °F).
5. Pour the liquid out of the wells and tap the microwell holder upside down vigorously (three times in a row) against absorbent paper to ensure complete removal of liquid from the wells. Fill all the wells with 250 µL wash buffer (see chapter 10.1.) and pour out the liquid again. Repeat two more times.
6. Add 100 µL of substrate/chromogen to each well. Mix gently by shaking the plate manually and incubate for 15 min at room temperature (20 - 25 °C / 68 - 77 °F) in the dark.
7. Add 100 µL of the stop solution to each well. Mix gently by shaking the plate manually and measure the absorbance at 450 nm. Read within 30 minutes after addition of stop solution.

11. Evaluation

Special software, **RIDASOFT® Win.NET (Art. No. Z9996FF)**, is optional available for evaluation of the RIDASCREEN® enzyme immunoassays.

For the evaluation it should be clarified, that all quality criteria are fulfilled for the current test run. The course of the standard curve is shown in the Quality Assurance Certificate (certificate of analysis) enclosed in the test kit.

Remark for the calculation without software:

$$\frac{\text{absorbance standard (or sample)}}{\text{absorbance zero standard}} \times 100 = B/B_0 (\%)$$

The zero standard is thus made equal to 100 % and the absorbance values are quoted in percentages. The values calculated for the standards are entered in a system of coordinates semilogarithmic against the AHD concentration [ng/L].

12. Result interpretation

In order to obtain the AHD concentration in ng/L (ng/kg) actually contained in a sample, the concentration read from the calibration curve must be further multiplied by the corresponding dilution factor. When working in accordance with the regulation stated, the **dilution factor is 2**.

13. Limits of the method

Test results may vary depending on the sample matrix, the actual test procedure and the laboratory environment.

Detection and quantification limits depend on the respective sample matrix, the degree of processing and the extraction method.

14. Recommendation

In order to ensure a high analytical performance we recommend to analyze each sample material in duplicates. Each laboratory may decide to perform the test in single determinations after a qualified risk management analysis. This has no influence on the function of the test kit. However, it should be noted that this increases the risk of overlooking errors in the performance of the test (e.g. pipetting errors). Moreover, a higher result variation will occur when pipetting in single determinations.

In order to ensure a high analytical performance we recommend:

- Pre-flush pipette tips with standard or sample extract prior to pipetting.
- To contact sales@r-biopharm.de if automates (e.g. ThunderBolt® / Bolt™) are used.

15. Further application notes

Application notes are available on request.

For further product information and applications, please contact your local distributor or R-Biopharm at this address: sales@r-biopharm.de.

Version overview

Version number	Chapter and title
2010-10-20	Release version
2015-10-01	General revision
2022-10-26	Current version General revision Changes made: <ul style="list-style-type: none">– Brief information: detection capability supplemented– Chapter 6 and section “Related products and accessories for nitrofurans determination” updated– Chapters 13, 14 and 15 added

Explanation of symbols

General symbols:

Follow the instructions for use

Batch number

Expiry date (YYYY-MM)

Storage temperature

Article number

Number of test determinations

Manufacturing date (YYYY-MM)

Manufacturer + address

Disclaimer

The user assumes all risk in using R-Biopharm AG's products and services.

R-Biopharm AG will warrant that its products and services meet all quality control standards set by R-Biopharm AG, and R-Biopharm AG will, at its option, replace or repair any components, product or repeat services which prove to be defective in workmanship or material within product specific warranty periods or expiration dates and which our examination shall disclose to our satisfaction to be defective as such.

This warranty is expressly in lieu of all other warranties, expressed or implied, as to quality, description, fitness for any particular purpose, merchantability, productiveness, or any other matter. R-Biopharm AG shall be in no way responsible for the proper use of its products and hereby disclaims all other remedies, warranties, guarantees or liabilities, expressed or implied, arising by law or otherwise, and it shall have no liability for any lost profits or damage, direct, indirect or otherwise, to person or property, in connection with the use of any of its products or services.

This warranty shall not be extended, altered or varied except by a written instrument signed by an authorized representative of R-Biopharm AG.

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are registered trademarks of R-Biopharm AG.