RIDA® QUICK Fumonisin RQS ECO

REF R5606

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RIDA®QUICK Fumonisin RQS ECO

Brief information

RIDA®QUICK Fumonisin RQS ECO (Art. No. R5606) is an immunochromatographic test for the quantitative determination of fumonisin in corn (see chapter 1. Intended use).

The test kit contains 20 test strips, each of which can be used for one analysis. All reagents required for the assay are contained in the test kit. Results are evaluated with

- the RIDA®SMART APP software (Art. No. ZRSAM) and an approved smartphone or
- the RIDA®SMART APP software (Art. No. ZRSAM) installed on an Android smartphone and the RIDA®SMART BOX (Art. No. ZRSA-SB)

Sample preparation: homogenization and extraction

Time requirement: sample preparation (for 10 samples)... approx. 10 min
test implementation (incubation time)..... 5 min

Limit of detection (LoD): > 0.3 mg/kg (ppm)

Measuring range: 0.3 - 3 mg/kg and 3 - 10 mg/kg
Quantitative evaluation with the RIDA®SMART APP

Specificity: The RIDA®QUICK Fumonisin RQS ECO test reacts with fumonisin in corn samples.

Further information is contained in the validation report.

The specificity of the RIDA®QUICK Fumonisin RQS ECO test was determined by analyzing the cross reactivities to corresponding substances in buffer system. In samples, the specificity may deviate from those determined in the buffer system due to matrix effects. Prior to the analysis of cross-reactive substances, the user has to determine the Limit of Detection and the Recovery

for the substance in the respective sample matrix. The test cannot discriminate between analytes and cross-reactive substances.

Related products and accessories for fumonisin determination

RIDASCREEN® Fumonisin (Art. No. R3411)

RIDASCREEN®FAST Fumonisin (Art. No. R5602)

RIDA®QUICK Mycotoxin ECO Extractor (Art. No. R5000)

RIDA®SMART APP software (Art. No. ZRSAM)

RIDA®SMART BOX (Art. No. ZRSA-SB)

Trilogy® certified reference material and standards for fumonisin (ISO 17034)

1. Intended use

RIDA®QUICK Fumonisin RQS ECO (Art. No. R5606) is an immunochromatographic test in strip format for the quantitative determination of fumonisin in corn.

2. General information

Fumonisin are carcinogenic, highly toxic metabolites of *Fusarium verticillioides*, which grows hostspecific on corn. The toxic concentrations of fumonisin differ significantly depending on the animal species.

3. Test principle

The immunochromatographic test in the form of a test strip is based on an antigen-antibody reaction. A specific anti-fumonisin antibody detects fumonisin in the sample. During incubation of the test strip, a continuous band pattern (test line and control line) needs to be formed that is used to determine the concentration of fumonisin. The intensity of the test line depends on the fumonisin concentration of the sample. It increases as the fumonisin concentration increases. A weak test band can also occur with a negative sample (contaminated below the detection limit). It must be possible to detect the control line after the reaction has taken place in order to verify the functioning of the test. The control line weakens as the fumonisin concentration in the sample increases.

The test strip is evaluated by smartphone using the RIDA®SMART APP software (Art. No. ZRSAM) or in combination with RIDA®SMART BOX (Art. No. ZRSA-SB).

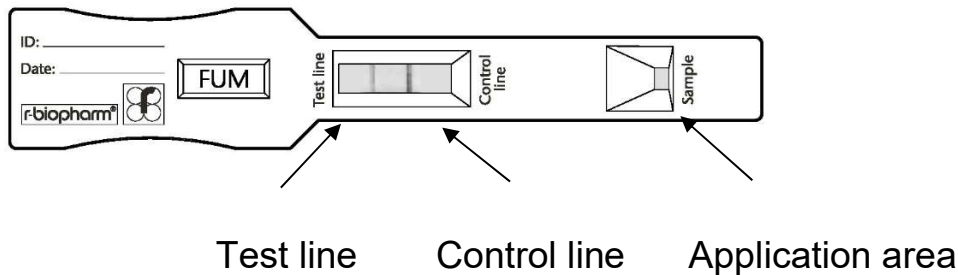
Fig. 1: RIDA®QUICK Fumonisin RQS ECO test strip.

4. Reagents provided

Each kit contains sufficient materials for max. 20 determinations. Each test kit contains:

Component	Cap color	Format	Volume
20 x Test strip	-	Ready to use	One determination per test strip (separately packed)
1 x Extraction buffer	Transparent	Concentrate 5x	100 mL
1 x RIDA®SMART APP cover	-	Ready to use	Evaluation cover for use with RIDA®SMART APP

5. Reagents required but not provided

5.1 Equipment

- Gloves
- Laboratory grinder
- Weighing scale
- Graduated cylinder
- Optional: Shaker
- Optional: Centrifuge (≥ 2000 g) or Filter paper (e.g. Whatman No. 1 or equivalent)
- 100 µl, 300 µl and 700 µl pipette
- Latest version of the RIDA®SMART APP software (Art. No. ZRSAM) activated and installed on a compatible smartphone
- Optional: RIDA®SMART BOX (Art. No. ZRSA-SB)

5.2 Reagents

- Distilled water (dist. Water) or deionized water

6. Warnings and precautions for the users

The product / test is only suitable within the scope of its intended use.

This test should only be carried out by trained laboratory personnel. The instruction for use must be strictly followed.

This kit may contain hazardous substances. For hazard notes on the contained substances please refer to the appropriate material safety data sheets (SDS) for this product, available online at www.r-biopharm.com.

All reagents and materials must be recovered or disposed after use at customers own responsibility according to the protection of human health and the environment. Please observe the applicable national regulations concerning waste disposal (e.g. Waste Management Act, Regulations on Dangerous Chemicals, etc.).

7. Storage instructions

Store the kit at 2 - 8 °C (36 - 46 °F). Do not freeze the test strips or any test kit components.

The test strips are sensitive to moisture. Moist test strips can have a negative impact on the test result. For this reason, it is important to protect the strips against moisture. Do not remove the test strips from the test strip packaging until immediately prior to use in the test.

Do not use the test kit after the expiration date (see test kit label).

Do not interchange individual reagents between kits of different lot numbers.

8. Sample preparation

Bring all components necessary for sample preparation to room temperature (20 - 25 °C / 68 - 77 °F) before use in the test and perform sample preparation at room temperature.

The sample should be stored in a cool place, protected against light.

Dilute the extraction buffer 1:5 (1+4) with distilled water. Store this ready to use buffer at 2 - 8 °C (36 - 46 °F). The expiry is not affected by dilution.

A representative sample (according to accepted sampling techniques) should be ground and thoroughly mixed prior to proceeding with the extraction procedure.

8.1 Measuring range 0.3 - 3 mg/kg

- Weigh 5 g of ground and homogenized sample into a suitable container and add 20 mL ready to use extraction buffer
- Close the tube and shake the sample vigorously for 5 min (manually or with shaker / vortex)
- Let the solution settle (≥ 3 min), filtrate or centrifuge
- Mix 100 μ L of the clear supernatant with 700 μ L ready to use extraction buffer (= dilution I)
- Apply 100 μ L of the dilution I on the application area of the test strip (see Fig. 1)

Note:

If necessary, the sample weight can be increased, e.g. 10 g sample in 40 mL ready to use extraction buffer. In this case the volume of ready to use extraction buffer has to be adapted.

8.2 Measuring range 3 - 10 mg/kg

For the measuring range of 3 - 10 mg/kg, the dilution I from sample preparation 8.1 must be further diluted.

- Mix 100 μ L of dilution I with 300 μ L ready to use extraction buffer (= dilution II)
- Apply 100 μ L of the dilution II on the application area of the test strip (see Fig. 1)

9. Test procedure

Bring all components necessary for sample preparation to room temperature (20 - 25 °C / 68 - 77 °F) before use in the test and perform sample preparation at room temperature.

- Remove a test strip from its foil pouch.
- Apply 100 µL of dilution I or dilution II (see chapter 8) on the application area of the test strip (see Fig. 1).
- Incubate the test strip at room temperature.
- Evaluate the result after 5 min (+ max. 15 s) incubation time with the RIDA®SMART APP.

10. Evaluation

The left line in the reaction area is the test band (test line, see Fig. 1). Its appearance and intensity depends on the fumonisin concentration of the sample. The sample is contaminated with fumonisin, if the control band (control line) is visible and the test band (test line) is also visible.

The right line in the reaction area is a control band (control line) and must be present after each test procedure. If this line is not visible, the test was not performed correctly or the reagents were not in order. In this case, the test should be repeated with a new test strip. If the right band is missing again, please inform your local distributor.


Please note:

- Fragmented bands:

Results will be invalid and test strips are inoperative if (testing or control) bands appear fragmented.


- Band detection:

The result field in the RIDA®SMART APP must always be visually checked. The correct two test- and control-bands (red bands) have to be framed as shown in Fig. 2. Results are invalid if another band is framed (see Fig. 3).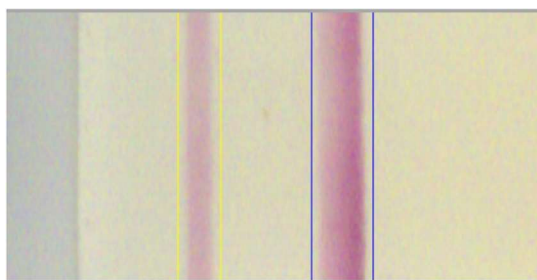

Test line Control line

Fig. 2: Band detection valid

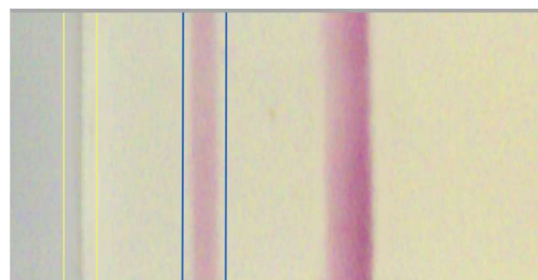Test line Control line

Fig. 3: Band detection invalid

– Reported result:

If the sample is highly contaminated with fumonisin, the control line will be only faintly visible. In this case the test result might be reported as “> 3 ppm or invalid result” (RIDA®SMART APP). In this instance, dilute the sample as described in chapter 8.2 and repeat the test with a new test strip.

To obtain comparable results, the test strips must always be evaluated after 5 min.

10.1 Evaluation with RIDA®SMART APP software in combination with an approved smartphone

First, please read the instructions for use of the RIDA®SMART APP software (Art. No. ZRSAM) attentively. The test strip evaluation has to be performed by using the RIDA®SMART APP software according to these guidelines. Please find a short description (Quick Starter Guide) of how to use the RIDA®SMART APP on our website www.r-biopharm.de. To analyze this test with RIDA®SMART APP software you must update to the latest software version. On the Certificate of analysis you find the latest version by scanning the QR Code.

For the test strip evaluation the RIDA®SMART APP cover is needed which is included in the test kit.

RIDA®SMART APP software applications for evaluation:

Matrix	Measuring range	RIDA®SMART APP application
Corn	0.3 - 3.0 mg/kg	0.3 – 3.0 ppm
Corn	3.0 - 10.0 mg/kg	3.0 – 10.0 ppm

10.2 Evaluation with RIDA®SMART APP software in combination with the RIDA®SMART BOX

First, please read the instructions for use of the RIDA®SMART APP software and of the RIDA®SMART BOX attentively. The test strip evaluation with RIDA®SMART APP software in combination with the RIDA®SMART BOX has to be performed according to these guidelines. Please find a short description (Quick Starter Guide) of how to use the RIDA®SMART APP on our website www.r-biopharm.com. To analyze this test with RIDA®SMART APP software you must update to the latest software version. On the Certificate of analysis you find the latest version by scanning the QR Code.

The RIDA®SMART software applications that need to be selected can be found in chapter 10.1.

11. Recommendation

In order to ensure a high analytical performance we recommend:

- In case of extremely acidic or basic samples, adjust the sample's pH value (pH 6.5 - 7.5) to neutral prior to extraction.
- In order to ensure accurate and correct test procedure, please analyze artificially contaminated samples

12. Sensitivity

This test determines fumonisin contaminations in the measurement range of 0.3 - 10 mg/kg in combination with the RIDA®SMART APP software.

13. Further application notes

Further application notes are available on request.

For further product information and applications, please contact your local distributor or R-Biopharm at this address: sales@r-biopharm.de.

Version overview

Version number	Chapter and title
2018-10-10	Release version
2021-10-27	General revision
2024-02-27	General revision
2025-05-05	Current version General revision Changes made: - 10. Evaluation RIDA®QUICK Scan deleted/RIDA®SMART BOX added - Change of article number RIDA®SMART APP New Disclaimer

Explanation of symbols

General symbols:

-  Follow the instructions for use
-  Batch number
-  Expiry date (YYYY-MM-DD)
-  Storage temperature
-  Article number
-  Number of test determinations
-  Manufacturing date (YYYY-MM-DD)
-  Manufacturer + address

Disclaimer

1. In conformance with the German Civil Code ("BGB") R-Biopharm AG provides a limited warranty ("Gewährleistung") against defects in design and manufacture present at delivery that make the Product unsuitable or unsafe for the contractually specified Product use and location through properly qualified personnel. Such limited warranty warrants Product title and against such material Product defects for 12 months, or if the Product is provided with a shelf life, the length of the declared shelf life, calculated from the date of risk transfer pursuant to delivery terms, and further provided customer provides timely and proper notice of the defect.
ALL OTHER WARRANTIES OR GUARANTIES, EXPRESS OR IMPLIED, OF ANY KIND ARE EXCLUDED, WHETHER IMPLIED BY CUSTOM, PRACTICE, THE COURSE OF DEALING BETWEEN THE PARTIES OR OTHERWISE. The responsibility for any express or implied warranties or extensions of R-Biopharm's own warranty made by dealers, distributors, or resellers is solely with such dealer, distributor, or reseller, and R-Biopharm AG expressly disclaims liability for such warranties.
2. R-Biopharm Products are designed for sole use by properly trained and qualified personnel applying appropriate industry standards and practices. Defects are covered only to the extent that such defects are demonstrably the result of defective material, design, parts, or faulty workmanship, which affect safety or result in substandard performance below specifications. R-Biopharm AG is not liable for any improper use nor the consequences of
 - a. the failure to read, understand and follow use and safety instructions, or use proper preparation and storage;
 - b. the failure to utilize trained and unqualified personnel, suitable samples or sampling techniques;
 - c. chemical, electromagnetic, mechanical, or electrolytic influences outside R-Biopharm AG provided standard perimeters through product specific data sheets, written product descriptions, or as otherwise expressly agreed upon in writing; or
 - d. any combination thereof.
3. R-Biopharm AG is also not liable for any changes or modifications, not carried out by R-Biopharm AG, nor consequences of use, applications, or processing which are unsafe or otherwise inconsistent with intended Product purpose as limited by written Product descriptions and specifications of R-Biopharm AG.
4. R-Biopharm AG's liability for ordinary breach of contract is limited to repair, replacement, other substitute performance, or refund. The choice of remedies is within R-Biopharm AG's sole discretion. R-Biopharm AG is not contractually liable for any incidental, or consequential damages, including but not limited to purchaser's expenses, losses, or damages from loss of good will, frustrated business purposes, sales expectations, investment loss, actual or anticipated lost profits, End User indemnity, or any other business expenditures.
5. The foregoing limited warranty is solely intended to fulfill the warranty requirements ("Gewährleistung") implied by German Civil Code. It is not intended to extend the scope or period of such Gewährleistung or provide additional warranties. Nor is this Disclaimer otherwise intended to limit or extent mandatory liability for damages in tort resulting from injury to life, body, or health, for intentional or grossly negligent acts, fraud, or due to strict liability under applicable product liability laws. A reversal of the burden of proof or rules of contract interpretation is not intended.

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are registered trademarks of R-Biopharm AG.