




# RIDA®SMART BOX connection: Access Point – quick starter guide

**1** Open your settings menu on your phone and go to WiFi settings.

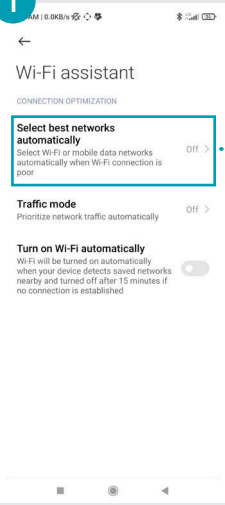


Before you connect the RIDA®SMART BOX to your phone in std. mode (access point), we recommend you to disable the WiFi setting „select best networks automatically“.

**2** Turn on the RIDA®SMART BOX, as soon as the blue light is on continuously you can select the box in the WiFi menu and establish a connection.

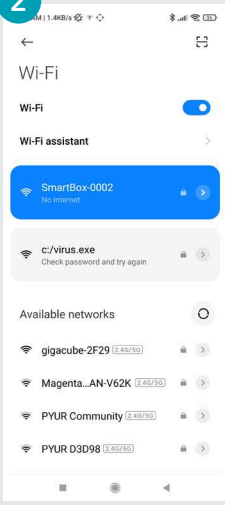



**Important:** only the first connection will ask for a password: **smartBoxSS1D**


**3** Now open the RIDA®SMART APP and go to the settings menu.

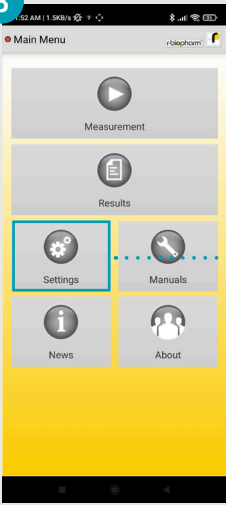

**4** Then enter the general settings menu.

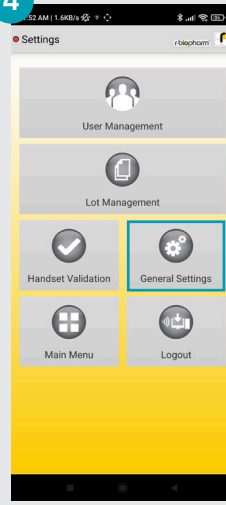


**5** Here you will find the RIDA®SMART BOX settings. Activate the use of the device by checking the box „use smart box“.

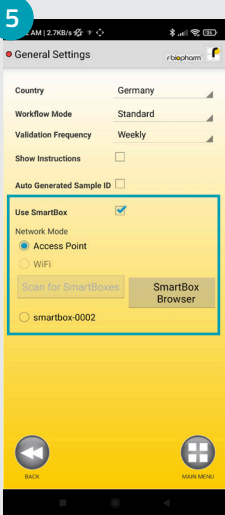

Select the access point mode and click „Scan for SmartBoxes“

The connected box appears, select it.


**6** When the RIDA®SMART BOX is successfully connected, dot in the top left corner turns green and the message “connection established with smartbox” appears.

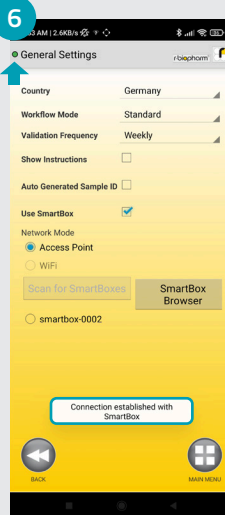


**Important:** the connection establishment may take a few seconds.

Go back to main menu and enter measurement menu.

**7**

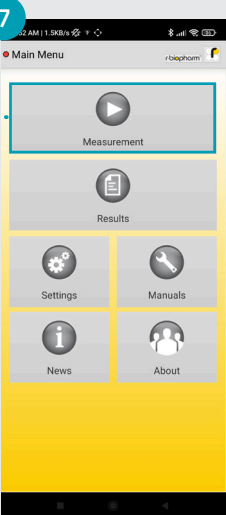

**8** Enter your sample data and choose your application.

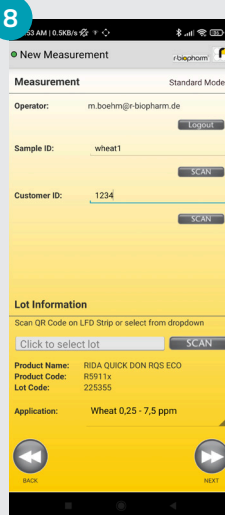

Click next.

The measurement starts and LED strip of the smart box turns purple. After measurement, the result appears as usual in the RIDA®SMART APP.

**9**

