**CONGEN**

# **SureFood® ANIMAL QUANT Turkey**

Art. No. S1015  
2 x 50 rxn

## **User Manual**

**September 2025**

 **Inhalt**

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## **Content**

|       |                                                      |    |
|-------|------------------------------------------------------|----|
| 1     | General Information .....                            | 10 |
| 1.1   | Description .....                                    | 10 |
| 1.2   | Limit of Detection and Limit of Quantification ..... | 10 |
| 1.3   | DNA-preparation .....                                | 11 |
| 1.4   | Kit components and storage .....                     | 11 |
| 1.5   | Additionally required equipment and materials .....  | 11 |
| 1.6   | Precautions for users .....                          | 11 |
| 1.7   | Setup.....                                           | 12 |
| 1.8   | Detection channel Set-up .....                       | 12 |
| 2     | Qualitative Analysis .....                           | 13 |
| 2.1   | Protocol .....                                       | 13 |
| 2.1.1 | Preparation of the master-mix .....                  | 13 |
| 2.1.2 | Preparation of the standard DNA dilutions .....      | 14 |
| 2.1.3 | Preparation of the real-time PCR-mix .....           | 14 |
| 2.2   | Interpretation of results .....                      | 15 |
| 3     | Limitations of the method .....                      | 16 |
| 4     | Further Information .....                            | 16 |
| 4.1   | Product Information .....                            | 16 |
| 4.2   | Technical Support .....                              | 16 |
| 4.3   | Distribution and Ordering .....                      | 16 |

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## **1 General Information**

### **1.1 Description**

The SureFood® ANIMAL QUANT Turkey real-time PCR detects the relative turkey (*Meleagris gallopavo*) DNA amount to the total animal (amniots) DNA content in meat samples. Therefore, the kit contains two PCR systems, one for detection of a turkey-specific gene (**Turkey** detection gene) and one for the detection of an animal gene (**Reference** gene).

The real-time PCR assay can be used with established real-time PCR instruments. The internal technical verification of instruments was performed on Agilent AriaDx, Agilent Mx3005P, Applied Biosystems 7500, Bio-Rad CFX96, Qiagen Rotor-Gene Q, R-Biopharm RIDA®CYCLER and Roche LightCycler® 480 II.

### **1.2 Limit of Detection and Limit of Quantification**

The turkey specific PCR has a limit of detection of  $\leq 5$  DNA-copies.

The assay limit of detection depends on sample matrix, processing grade, DNA-preparation and DNA-content.

The limit of quantitation depends on the concentration of the sample DNA used in the analysis. For example, if 50,000 target-sequence copies of the reference gene are present, the relative quantitation limit for turkey DNA is 0.1%.

The SureFood® PCR systems are very sensitive and therefore even a small amount of target DNA is sufficient for a successful analysis. The concentration of total DNA in the sample does not allow a conclusion on the quantity and quality of the target DNA.

# SureFood® ANIMAL QUANT Turkey

## (2 x 50 rxn)

Art. No. S1015

September 2025

### 1.3 DNA-preparation

For DNA-preparation of meat samples the use of SureFood® PREP Basic (Art. No. S1052) or SureFood® PREP Advanced (Art. No. S1053) is recommended.

### 1.4 Kit components and storage

| Kit Code | Reagent                         | Amount      | Lid Color  |
|----------|---------------------------------|-------------|------------|
| 1        | Reference Reaction Mix          | 2 x 500 µl  | Orange     |
| 2        | Turkey Reaction Mix             | 2 x 500 µl  | Yellow     |
| 3        | Taq Polymerase                  | 1 x 11 µl   | Red        |
| 4        | Dilution Buffer                 | 1 x 1300 µl | White      |
| 5        | Standard DNA                    | 1 x 45 µl   | Dark Blue  |
| 6        | Positive Control (100 % turkey) | 1 x 80 µl   | Light Blue |

**Store all reagents at -28 to -16°C and protected from light. The tube of the Taq Polymerase should not be thawed.**

### 1.5 Additionally required equipment and materials

- DNA-Extraction kit  
(e.g. SureFood® PREP Basic Art. No. S1052 / SureFood® PREP Advanced Art. No. S1053)
- real- time PCR instrument
- real-time PCR consumable (plates, tubes, foils, caps)
- pipettes with filter tips
- powder-free disposable gloves
- Vortex mixer
- micro centrifuge with a rotor for the reaction tubes

### 1.6 Precautions for users

- Extraction, PCR preparation and the PCR run should be separated in different rooms to avoid cross-contaminations.
- This test must only be performed by laboratory personnel trained in molecular biology methods.
- Strictly follow the working instructions.
- When handling samples, wear disposable gloves. After finishing the test, wash your hands.
- Do not smoke, eat or drink in areas where samples or test reagents are being used.
- Do not use the kit after the expiration date.
- All reagents and materials used have to be disposed properly after use. Please refer to the relevant national regulation for disposal.

# SureFood® ANIMAL QUANT Turkey (2 x 50 rxn)

Art. No. S1015

September 2025

## 1.7 Setup

|                                        | Blockcycler &<br>R-Biopharm RIDA®CYCLER | Rotorcycler       |
|----------------------------------------|-----------------------------------------|-------------------|
| Initial Denaturation (HOLD)<br>Cycles  | 5 min, 95°C<br>45                       | 1 min, 95°C<br>45 |
| Denaturation                           | 15 sec, 95°C                            | 10 sec, 95°C      |
| Annealing/Extension (CYCLE)            | 30 sec, 60°C                            | 15 sec, 60°C      |
| Temperature Transition Rate/ Ramp Rate | Maximum                                 | Maximum           |

## 1.8 Detection channel Set-up

| Real-time PCR device             | Detection | Detection channel | Quencher | Note                                                                                                                                                                                                           |
|----------------------------------|-----------|-------------------|----------|----------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| <b>Agilent Mx3005P</b>           | Turkey    | FAM               | +        |                                                                                                                                                                                                                |
|                                  | Reference | FAM               | +        |                                                                                                                                                                                                                |
| <b>Agilent AriaDx /Mx</b>        | Turkey    | FAM               | +        |                                                                                                                                                                                                                |
|                                  | Reference | FAM               | +        |                                                                                                                                                                                                                |
| <b>Applied Biosystems 7500</b>   | Turkey    | FAM               | None     | Check the passive reference option ROX is none.                                                                                                                                                                |
|                                  | Reference | FAM               | None     |                                                                                                                                                                                                                |
| <b>Bio-Rad CFX96 / Dx / Opus</b> | Turkey    | FAM               | +        | Baseline Settings:<br>• Baseline subtracted curve fit<br>• Apply fluorescence drift correction                                                                                                                 |
|                                  | Reference | FAM               | +        |                                                                                                                                                                                                                |
| <b>R-Biopharm RIDA®CYCLER</b>    | Turkey    | green             | +        | <b>Ignore cycles before</b> , if there is a significant deviation in the baseline at the start of the run. Please see page 45 of the cyclor operating instructions, section 12.1.2 Cycling analysis parameter. |
|                                  | Reference | green             | +        |                                                                                                                                                                                                                |
| <b>Qiagen Rotor-Gene Q</b>       | Turkey    | green             | +        | <b>Note:</b> Please use only 0.1 ml reaction tubes. The gain settings must be set to 5 (factory default) for all channels.                                                                                     |
|                                  | Reference | green             | +        |                                                                                                                                                                                                                |
| <b>Roche LightCycler® 480 II</b> | Turkey    | 465-510           | +        | For the quantitative analysis use the type "Fit Points".                                                                                                                                                       |
|                                  | Reference | 465-510           | +        |                                                                                                                                                                                                                |

## **2 Qualitative Analysis**

### **2.1 Protocol**

#### **2.1.1 Preparation of the master-mix**

Calculate the total number of reactions needed (samples, control reactions and standards) for the specific PCR assay.

#### **Required reactions for the evaluation of the reference detection (see 2.2):**

5 reactions for the standard curve

3 reactions for controls\* (1x negative control, 2x Positive Control)

For each sample: at least 1 reaction for each sample DNA

#### **Required reactions for the evaluation of turkey detection (see 2.2):**

5 reactions for the standard curve

3 reactions for controls\* (1x negative control, 2x Positive Control)

For each sample: at least 1 reaction for each sample DNA

It is also recommended to prepare the master-mix with 10 % additional volume in order to compensate reagent loss. Allow the reagents to thaw, mix by vortexing and centrifuge before opening and use.

The tube of the Taq Polymerase should not be mixed by vortexing.

#### **Example for the calculation and preparation of 10 reactions:**

| <b>Components of the master-mix</b> | <b>Amount per reaction</b> | <b>10 reactions (with 10 % excess)</b> |
|-------------------------------------|----------------------------|----------------------------------------|
| Reaction Mix                        | 19.9 µl                    | 218.9 µl                               |
| Taq Polymerase                      | 0.1 µl                     | 1.1 µl                                 |
| <b>Total volume</b>                 | <b>20 µl</b>               | <b>220 µl</b>                          |

**Mix each master-mix well and centrifuge shortly before use.**

#### **\* Description of the controls**

- Negative control: only master-mix
- Positive Control: master-mix and within the kit's provided Positive Control

### 2.1.2 Preparation of the standard DNA dilutions

Dilute the Standard DNA (**Code 5**) in 1:10 steps in Dilution Buffer (**Code 4**) in order to prepare different DNA concentrations for the standard curves of the reference gene (**Reference**) and the detection gene (**Turkey**). Prepare 5 dilutions of the supplied Standard DNA (**Code 5**) with the supplied Dilution Buffer (**Code 4**). Prepare 5 reaction tubes (labeled S1 to S5) and add 45 µl Dilution Buffer (**Code 4**) each.

**The following procedure is recommended:**

| Standard | Dilutions                                 | Copy number per µl | Final copy number per reaction** |
|----------|-------------------------------------------|--------------------|----------------------------------|
| S1       | 45 µl Dilution Buffer + 5 µl Standard DNA | 100,000 copies     | 500,000 copies                   |
| S2       | 45 µl Dilution Buffer + 5 µl DNA of S1    | 10,000 copies      | 50,000 copies                    |
| S3       | 45 µl Dilution Buffer + 5 µl DNA of S2    | 1,000 copies       | 5,000 copies                     |
| S4       | 45 µl Dilution Buffer + 5 µl DNA of S3    | 100 copies         | 500 copies                       |
| S5       | 45 µl Dilution Buffer + 5 µl DNA of S4    | 10 copies          | 50 copies                        |

**\*\*Note:** 5 µl of standard DNA are used for each calibration point. The final copy number per reaction is to be entered in the analysis software of the real-time PCR detection system.

### 2.1.3 Preparation of the real-time PCR-mix

- Pipette 20 µl of the master-mix into appropriate tubes/wells.
- Close the negative control.
- Pipette 5 µl of sample DNA into the designated tubes/wells and close them.
- Pipette 5 µl of Positive Control and the standard dilutions into the designated tubes/wells and close them.
- Centrifuge all tubes/plates shortly at low speed.
- Place tubes/plates into the real-time PCR instrument and start the run according to the setup.

## 2.2 Interpretation of results

The calculation for both reaction systems (**Reference**, **Turkey**) has to be made separately. Mark the standards, the controls and the samples for the specific system (**Turkey**) and make the evaluation according to the usual analysis program recommended by the real-time PCR instrument manufacturer. Repeat the same procedure for the reference gene system (**Reference**). The value for the slope of the standard curve has to be between -3.1 and -3.6 and the correlation coefficient  $R^2 > 0.98$ . In case of different values for the standard curve, it should not be used for calculation.

By using the calculated copy numbers for **Reference** and **Turkey** the relative turkey content of the sample DNA and the Positive Control can be determined in the following way (example):

|                         |               |                                   |              |
|-------------------------|---------------|-----------------------------------|--------------|
| Sample <b>Turkey</b>    | 1,350 copies  | Positive Control <b>Turkey</b>    | 900 copies   |
| Sample <b>Reference</b> | 45,000 copies | Positive Control <b>Reference</b> | 1,100 copies |

Divide the copy number of the specific system by the copy number of the reference gene system and multiply by 100 to obtain the percentage.

**Turkey** DNA content = **Turkey** copy number \* 100 / **Reference** copy number

sample **Turkey** DNA content =  $1,350 * 100 / 45,000$       sample **Turkey** DNA content = 3 %

For the given example the numbers lead to a **Turkey** DNA content of 3.0 %, for the sample and 81.8 % for the Positive Control with the same calculation.

For a final calculation the use of a correction factor K for the correction of run-to-run fluctuations is necessary. The correction factor is the relation of the true percentage value of the Positive Control (100% **Turkey** content) and the measured percentage of the Positive Control.

$K = \text{true value} / \text{measured value}$        $K (\text{example}) = 100\% / 81.8\% = 1.2$

The measured DNA content for the sample is multiplied with K to obtain a corrected DNA content.

sample DNA content = measured sample DNA content \* K      sample (example) =  $3.0\% * 1.2 = 3.6\%$

For that example the **Turkey DNA content is 3.6 %**.

## **3 Limitations of the method**

- The presence of PCR inhibitors may cause invalid results.
- Extremely low levels of target below the limit of detection (LoD) may be detected, but results may not be reproducible.
- Innards, particularly kidney or liver, show a higher DNA content per weight unit when compared to muscle tissue. This could lead to over-estimation of the species content of the animal from which the innards come from. For meat products with innards content > 5 % care must be taken when relative species contents are reported.
- In highly processed samples, the limit of detection may be shifted. Factors such as high pressures, mechanical stresses, chemical treatment, extreme temperatures and/or extreme pH values during manufacturing process – such as in canning production – can damage or degrade nucleic acids. This means that the sensitivity of the test kit may be reduced and not all original components may be detected.

## **4 Further Information**

### **4.1 Product Information**

- Microsoft Excel template of calculation and detailed information about setup of several real-time PCR devices (Download: [www.congen.de/en/downloads](http://www.congen.de/en/downloads))
- Product-related documents (Download: [www.congen.de/en/eifu/](http://www.congen.de/en/eifu/))
- Validation Report upon request

### **4.2 Technical Support**

For further questions please contact your distributor or send an e-mail to [sales@r-biopharm.de](mailto:sales@r-biopharm.de).

### **4.3 Distribution and Ordering**

R-Biopharm AG  
An der neuen Bergstrasse 17,  
64297 Darmstadt, Germany  
Phone: +49 (0) 61 51 - 81 02-0  
Fax: +49 (0) 61 51 - 81 02-20  
E-Mail: [orders@r-biopharm.de](mailto:orders@r-biopharm.de)  
[www.r-biopharm.com](http://www.r-biopharm.com)