**CONGEN**

# SureFood<sup>®</sup> PREP Add-On

Art. No. S1055  
15 extractions

## User Manual

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# 1 General Information

## 1.1 Description

This kit is intended to be used for the extraction of animal and plant DNA (deoxyribonucleic acid) from raw materials as well as processed food and feed with sample weight of 2 g. It is used in conjunction with the SureFood® PREP Basic (Art. No. S1052) or SureFood® PREP Advanced (Art. No. S1053).

## 1.2 Kit components and storage

Kit Code	Reagent /Material	Amount (per box)
<b>K</b>	Proteinase K	2 Tubes
<b>H</b>	PCR grade water	2 x 1,1 ml
<b>L</b>	Lysis Buffer	3 x 60 ml

The Proteinase K should be stored at -28 bis -16°C. All other reagents of the kit should be stored dry and at room temperature (14 – 25 °C).

Before every extraction make sure that all components have room temperature. If there are any precipitates within the provided solutions, solve them by warming carefully (up to 30°C).

## 1.3 Additionally required equipment and materials

- DNA-Extraction kit (e.g. SureFood® PREP Advanced Art. No. S1053 / SureFood® PREP Basic Art. No. S1052)
- suitable equipment for sample comminution and homogenization
- micro balance and spatula for weighing the samples
- suitable reaction tubes free from DNA and DNase (in minimum 15 ml)
- waterproof pen and tags for labeling the reaction tubes
- unpowdered disposable gloves
- pipettes with filter tips
- Vortex mixer
- Thermomixer/ heating block for suitable centrifugation tubes (up to 65°C)
- centrifuge for suitable centrifugation tubes (up to 7,000 rpm)
- micro centrifuge (up to 12,000 rpm)
- disposal bags or waste bin

## 2 Protocol

### 2.1 Principle

1. Preparation of the basic material
2. Lysis of the basic material
3. Pre-filtration and setting of optimal binding conditions

### 2.2 Preparations

Add 1 ml PCR grade water (**Code H**) to the Proteinase K (**Code K**) and mix thoroughly.

### 2.3 Extraction protocol

#### 1. Preparation of the basic material

Transfer 2 g homogenized sample material into a suitable centrifugation tube (not provided with the kit).

#### 2. Lysis of the basic material

Add 10 ml of Lysis Buffer (**Code L**) and 100 µl of Proteinase K (**Code K**) to the sample.

Mix it well on the vortex mixer.

Incubate on a Thermomixer/heating block under continuous shaking for 30 minutes at 65°C.

#### 3. Pre-filtration and setting of optimal binding conditions

Centrifuge the sample lysate for 1 min at 7,000 rpm.

For duplicates place two Spin Filters (**Code F – from SureFood® PREP Basic or SureFood® PREP Advanced**) into two 2.0 ml clear Receiver Tubes (**Code R from SureFood® PREP Basic or SureFood® PREP Advanced**).

Transfer 400 µl of the liquid supernatant directly onto each Spin Filter (**Code F**).

Centrifuge the Spin Filters with the Receiver Tubes for 1 min at 12,000 rpm. After centrifugation discard the Spin Filters.

**Note:** For non-processed animal material the usage of the Spin Filter (**Code F**) is superfluous. Transfer the liquid supernatant directly into two 2.0 ml Receiver Tubes (**Code R**).

For the following extraction steps (see 2.3/3.3 Extraction protocol) use the SureFood® PREP Basic or SureFood® PREP Advanced beginning with point 4 (Binding of the nucleic acids on a Spin Filter).

## 3 Further Information

### 3.1 Product Information

- Material Safety Data Sheet (Download: [www.congen.de/en/eifu/](http://www.congen.de/en/eifu/))

### 3.2 Technical Support

For further questions please contact your distributor or send an e-mail to [sales@r-biopharm.de](mailto:sales@r-biopharm.de).

### 3.3 Distribution and ordering

R-Biopharm AG  
An der neuen Bergstrasse 17,  
64297 Darmstadt, Germany  
Phone: +49 (0) 61 51 - 81 02-0  
Fax: +49 (0) 61 51 - 81 02-20  
E-Mail: [orders@r-biopharm.de](mailto:orders@r-biopharm.de)  
[www.r-biopharm.com](http://www.r-biopharm.com)

## 4 Safety information

### Lysis Buffer

#### Warning

H319-H412  
P264-P273-P280-P305+P351+P338-  
P337+P313-P501

### Proteinase K

#### Danger

H315-H319-H334-H335  
P261-P264-P280-P304+P340-P312-  
P321-P337+313-P342+P311

<b>H315:</b>	Causes skin irritation.
<b>H319:</b>	Causes serious eye irritation.
<b>H334:</b>	May cause allergy or asthma symptoms or breathing difficulties if inhaled.
<b>H335:</b>	May cause respiratory irritation.
<b>H412:</b>	Harmful to aquatic life with long lasting effects.
<b>P261:</b>	Avoid breathing dust/fumes/gas/mist/vapours/spray.
<b>P264:</b>	Wash hands, forearms and face thoroughly after handling.
<b>P273:</b>	Avoid release to the environment.
<b>P280:</b>	Wear protective gloves/protective clothing/face protection.
<b>P304+P340:</b>	IF INHALED: Remove person to fresh air and keep comfortable for breathing.
<b>P305+P351+P338:</b>	IF IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing.
<b>P312:</b>	Call a POISON CENTER/ doctor if you feel unwell
<b>P321:</b>	Specific treatment (see supplemental first aid instruction on this label).
<b>P337+P313:</b>	If eye irritation persists get medical advice/attention.
<b>P342+P311:</b>	If experiencing respiratory symptoms: Call a POISON CENTER / doctor.
<b>P501:</b>	Dispose of contents/container to hazardous or special waste collection point, in accordance with local, regional, national and/or international regulation.

For further information we offer a Material Safety Data Sheet, see at [www.congen.de/en/eifu/](http://www.congen.de/en/eifu/).  
Alternatively please contact your distributor or send an e-mail to [sales@r-biopharm.de](mailto:sales@r-biopharm.de).