CONGEN

SureFood[®] PREP Add-On

Art. No. S1055
15 extractions

User Manual

Efficient DNA preparation from food and feed

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Content

1	General Information	7
1.1	Description	7
1.2	Kit components and storage	7
1.3	Additionally required equipment and materials	7
2	Protocol	8
2.1	Principle	8
2.2	Preparations	8
2.3	Extraction protocol	8
	1. Preparation of the basic material	8
	2. Lysis of the basic material	8
	3. Pre-filtration and setting of optimal binding conditions	8
3	Further Information	9
3.1	Product Information	9
3.2	Technical Support	9
3.3	Distribution and ordering	9
4	Safety information	10

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1 General Information

1.1 Description

This kit is intended to be used for the extraction of animal and plant DNA (deoxyribonucleic acid) from raw materials as well as processed food and feed with sample weight of 2 g. It is used in conjunction with the SureFood® PREP Basic (Art. No. S1052) or SureFood® PREP Advanced (Art. No. S1053).

1.2 Kit components and storage

Kit Code	Reagent /Material	Amount (per box)
K	Proteinase K	2 Tubes
H	PCR grade water	2 x 1,1 ml
L	Lysis Buffer	3 x 60 ml

The Proteinase K should be stored at -20°C. All other reagents of the kit should be stored dry and at room temperature (14 – 25 °C).

Before every extraction make sure that all components have room temperature. If there are any precipitates within the provided solutions, solve them by warming carefully (up to 30°C).

1.3 Additionally required equipment and materials

- DNA-Extraction kit (e.g. SureFood® PREP Advanced Art. No. S1053 / SureFood® PREP Basic Art. No. S1052)
- suitable equipment for sample comminution and homogenization
- micro balance and spatula for weighing the samples
- suitable reaction tubes free from DNA and DNase (in minimum 15 ml)
- waterproof pen and tags for labeling the reaction tubes
- unpowdered disposable gloves
- pipettes with filter tips
- Vortex mixer
- Thermomixer/ heating block for suitable centrifugation tubes (up to 65°C)
- centrifuge for suitable centrifugation tubes (up to 7,000 rpm)
- micro centrifuge (up to 12,000 rpm)
- ethanol (purity ≥ 96 %) for preparation of Pre-Wash Buffer and Wash Buffer
- disposal bags or waste bin

2 Protocol

2.1 Principle

1. Preparation of the basic material
2. Lysis of the basic material

2.2 Preparations

Add 1 ml PCR grade water (**Code H**) to the Proteinase K (**Code K**) and mix thoroughly.

2.3 Extraction protocol

1. Preparation of the basic material

Transfer 2 g homogenized sample material into a suitable centrifugation tube (not provided with the kit).

2. Lysis of the basic material

Add 10 ml of Lysis Buffer (**Code L**) and 100 µl of Proteinase K (**Code K**) to the reaction tube and mix it well on the vortex mixer. Incubate on a heating block under continuous shaking for 30 minutes at 65°C.

3. Pre-filtration and setting of optimal binding conditions

Centrifuge the sample lysate for 1 min at 7,000 rpm.

For duplicates place two green Spin Filters (**Code F – from SureFood® PREP Basic or SureFood® PREP Advanced**) into two clear 2.0 ml Receiver Tubes (**Code R from SureFood® PREP Basic or SureFood® PREP Advanced**). Transfer 400 µl of the liquid supernatant directly onto each Spin Filter.

Centrifuge the Spin Filters with the Receiver Tubes for 1 min at 12,000 rpm. After centrifugation discard the Spin Filters.

Note: For non-processed animal material the usage of the green Spin Filter is superfluous. Transfer the liquid supernatant directly into two clear 2.0 ml Receiver Tubes (**Code R**).

For the following extraction steps use the SureFood® PREP Basic or SureFood® PREP Advanced beginning with point 4 (Binding of the nucleic acids on a Spin Filter).

3 Further Information

3.1 Product Information

- Material Safety Data Sheet

3.2 Technical Support

For further questions please contact your distributor or send an e-mail to sales@r-biopharm.de.

3.3 Distribution and ordering

R-Biopharm AG
An der neuen Bergstrasse 17,
64297 Darmstadt, Germany
Phone: +49 (0) 61 51 - 81 02-0
Fax: +49 (0) 61 51 - 81 02-20
E-Mail: orders@r-biopharm.de
www.r-biopharm.com

4 Safety information

Lysis Buffer

Warning

H319 H412

P280 P273 P305+P351+P338

Proteinase K

Danger

H315 H319 H334 H335

P280 P305+P351+P338 P261 P271 P337+P313

P342+P311 P304+P340 P312 P302+P352

P332+P313 P362

- H315:** Causes skin irritation.
- H319:** Causes serious eye irritation.
- H334:** May cause allergy or asthma symptoms or breathing difficulties if inhaled.
- H335:** May cause respiratory irritation.
- H412:** Harmful to aquatic life with long lasting effects.
- P273:** Avoid release to the environment.
- P261:** Avoid breathing dust/fumes/gas/mist/vapours/spray.
- P264:** Wash hands thoroughly after handling.
- P271:** Use only outdoors or in a well-ventilated area.
- P280:** Wear protective gloves/protective clothing/eye protection/face protection.
- P305+P351+P338:** If in eyes: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing.
- P337+P313:** If eye irritation persists get medical advice/attention.
- P342+P311:** If experiencing respiratory symptoms: Call a poison center or doctor/physician.
- P304+P340:** If inhaled: Remove person to fresh air and keep comfortable for breathing.
- P312:** Call a poison center/ doctor if you feel unwell.
- P302+P352:** If on skin: Wash with plenty of soap and water.
- P332+P313 P362:** If skin irritation occurs: Get medical advice/attention. Take off contaminated clothing.

For further information we offer a Material Safety Data Sheet. Please contact your distributor or send an e-mail to sales@r-biopharm.de.