CONGEN

SureFood® GMO ID 4plex Soya II

Art. No. S2162
100 rxn

User Manual

March 2026

 Inhalt

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Content

1	General Information	10
1.1	Description	10
1.2	Limit of Detection	10
1.3	DNA-preparation	11
1.4	Kit components and storage	11
1.5	Additionally required equipment and materials	11
1.6	Precautions for users	11
1.7	Setup.....	12
1.8	Detection channel Set-up	12
2	Qualitative Analysis	13
2.1	Protocol	13
2.1.1	Preparation of the master-mix	13
2.1.2	Preparation of the real-time PCR-mix	13
2.2	Interpretation of results	14
3	Limitations of the method	16
4	Further Information	16
4.1	Product Information	16
4.2	Technical Support	16
4.3	Distribution and Ordering	16

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1 General Information

1.1 Description

The SureFood® GMO ID 4plex Soya II is a real-time PCR for the direct, qualitative detection and differentiation of following specific genetically modified soya DNA sequences:

- Roundup Ready Soya (RR Soya) (OECD unique identifier MON-Ø4Ø32-6)
- Roundup Ready 2 Yield Soya (RR2Y Soya) (OECD unique identifier MON-89788-1)
- A2704-12 Soya (OECD unique identifier ACS-GMØØ5-3)
- A5547-127 Soya (OECD unique identifier ACS-GMØØ6-4)

This kit can be used for the screening of genetically modified organisms (GMOs) in food, feed and seeds.

The detections are according to the validated method of the European Commission respectively official collection of detection methods of §64 German food law.

The real-time PCR assay can be performed with commonly used real-time PCR instruments, equipped for detection of four fluorescence emissions at the channels FAM, VIC/HEX, ROX and Cy5 at the same time. The internal technical verification of instruments was performed on Agilent Mx3005P, Applied Biosystems 7500, Bio-Rad CFX96, R-Biopharm RIDA®CYCLER, Roche LightCycler® 480 II, Roche cobas® z 480 Analyzer and Qiagen Rotor-Gene Q.

1.2 Limit of Detection

The SureFood® GMO ID 4plex Soya II real-time PCR has a limit of detection of ≤ 5 DNA copies. This is equivalent to approx. 0.01 % for unprocessed soybean.

The assay limit of detection depends on sample matrix, processing grade, DNA preparation and DNA content.

The SureFood® PCR systems are very sensitive and therefore even a small amount of target DNA is sufficient for a successful analysis. The concentration of total DNA in the sample does not allow a conclusion on the quantity and quality of the target DNA.

Note: Inconsistent mixing ratios* may cause a loss of sensitivity in the low concentration channel in mixed samples especially with high amplicon concentrations (Cp value < 27).

* e.g. 99.9 % RR Soya and 0.1 % A5547-127 Soya

1.3 DNA-preparation

For DNA-preparation of raw material the use of SureFood® PREP Basic (Art. No. S1052), SureFast® Mag PREP Food (Art. No. F1060) and for highly processed food and feed the use of SureFood® PREP Advanced (Art. No. S1053) is recommended. SureFood® PREP Add On (Art. No. S1055) is intended to be used for the extraction of DNA from raw materials as well as processed food and feed with sample weight of 2 g. It is used in conjunction with the SureFood® PREP Basic.

1.4 Kit components and storage

Kit Code	Reagent	Amount	Lid Color
1	Reaction Mix	2 x 1050 µl	Yellow
2	Taq Polymerase	1 x 80 µl	Dark Red
3	Positive Control	1 x 200 µl	Light Blue

Store all reagents at -28°C to -16°C and protected from light. The Taq Polymerase can be stored at +2 to +8°C for multiple uses on the same day.

Note: The Taq Polymerase may be in a frozen or unfrozen state. This does not affect the quality of the Taq Polymerase or the performance of the real-time PCR.

1.5 Additionally required equipment and materials

- DNA-Extraction kit
(e.g. SureFood® PREP Basic Art. No. S1052 / SureFood® PREP Advanced Art. No. S1053 / SureFood® PREP Add On Art. No. S1055 / SureFast® Mag PREP Food Art. No. F1060)
- real- time PCR instrument with four detection channels (510 nm, 580 nm, 610 nm and 660 nm)
- real-time PCR consumable (plates, tubes, foils, caps)
- pipettes with filter tips
- powder-free disposable gloves
- Vortex mixer
- micro centrifuge with a rotor for the reaction tubes

1.6 Precautions for users

- Extraction, PCR preparation and the PCR run should be separated in different rooms to avoid cross-contaminations.
- This test must only be performed by laboratory personnel trained in molecular biology methods.
- Strictly follow the working instructions.
- When handling samples, wear disposable gloves. After finishing the test, wash your hands.
- Do not smoke, eat or drink in areas where samples or test reagents are being used.
- Do not use the kit after the expiration date.
- All reagents and materials used have to be disposed properly after use. Please refer to the relevant national regulation for disposal.

SureFood® GMO ID 4plex Soya II (100 rxn)

Art. No. S2162

March 2026

1.7 Setup

	Blockcycler & R-Biopharm RIDA®CYCLER	Rotorcycler
Initial Denaturation (HOLD)	5 min, 95°C	1 min, 95°C
Cycles	45	45
Denaturation	15 sec, 95°C	10 sec, 95°C
Annealing/Extension (CYCLE)	30 sec, 60°C	15 sec, 60°C
Temperature Transition Rate/ Ramp Rate	Maximum	Maximum

1.8 Detection channel Set-up

For additional details see cycler operating instructions for real-time PCR device.

Real-time PCR device	Detection	Detection channel	Quencher	Note
Agilent Mx3005P	RR-Soya	FAM	+	
	RR2Y-Soya	HEX	+	
	A2704-12-Soya	ROX	+	
	A5547-127-Soya	Cy5	+	
Applied Biosystems 7500	RR-Soya	FAM	None	Check the passive reference option ROX is none.
	RR2Y-Soya	VIC	None	
	A2704-12-Soya	ROX	None	
	A5547-127-Soya	Cy5	None	
Bio-Rad CFX96/Dx/Opus	RR-Soya	FAM	+	Baseline Settings: - Baseline subtracted curve fit - Apply fluorescence drift correction
	RR2Y-Soya	VIC/HEX	+	
	A2704-12-Soya	ROX	+	
	A5547-127-Soya	Cy5	+	
R-Biopharm RIDA®CYCLER	RR-Soya	green	+	Ignore cycles before , if there is a significant deviation in the baseline at the start of the run. Please see page 45 of the cycler operating instructions, section 12.1.2 Cycling analysis parameter.
	RR2Y-Soya	yellow	+	
	A2704-12-Soya	orange	+	
	A5547-127-Soya	red	+	
Qiagen Rotor-Gene Q	RR-Soya	green	+	Note: Please use only 0.1 ml reaction tube. The gain settings must be set to 5 (factory default) for all channels.
	RR2Y-Soya	yellow	+	
	A2704-12-Soya	orange	+	
	A5547-127-Soya	red	+	
Roche LightCycler® 480 II	RR-Soya	465-510	+	The SureCC Color Compensation Kit I (Art. No. F4009) is required.
	RR2Y-Soya	533-580	+	
	A2704-12-Soya	533-610	+	
	A5547-127-Soya	618-660	+	
Roche cobas® z 480 Analyzer	RR-Soya	465-510	+	The SureCC Color Compensation Kit I (Art. No. F4009) is required.
	RR2Y-Soya	540-580	+	
	A2704-12-Soya	540-610	+	
	A5547-127-Soya	610-670	+	

2 Qualitative Analysis

2.1 Protocol

2.1.1 Preparation of the master-mix

Calculate the total number of reactions needed (samples and control reactions) for the specific PCR assay as well as for the inhibition control.

The following control reactions are needed for the specific PCR assay: negative control, extraction control, Positive Control and an external inhibition control per sample.

For the preparation of the inhibition control, it is recommended to use the SureFood® GMO Plant PLUS (Art. No. S2049) or the SureFood® GMO SCREEN 4plex 35S/NOS/FMV+IAC (Art. No. S2126).

Reactions needed for the qualitative RR-Soya, RR2Y-Soya, A2704-12-Soya and A5547-127-Soya detection:

3 reactions for controls* (1x negative control, 1x extraction control, 1x Positive Control)

For each sample: at least 1 reaction for each sample DNA

It is also recommended to prepare the master-mix with 10 % additional volume in order to compensate reagent loss. Allow the reagents to thaw, mix and centrifuge before opening and use.

Example for the calculation and preparation of 10 reactions:

Components of the master-mix	Amount per reaction	10 reactions (with 10 % excess)
Reaction Mix	19.3 µl	212.3 µl
Taq Polymerase	0.7 µl	7.7 µl
Total volume	20 µl	220 µl

Mix each master-mix well and centrifuge shortly before use.

2.1.2 Preparation of the real-time PCR-mix

- Pipette 20 µl of the master-mix into appropriate tubes/wells.
- Close the negative control.
- Pipette 5 µl of sample DNA into the designated tubes/wells and close them.
- Pipette 5 µl of Positive Control into the designated tubes/wells and close them.
- Centrifuge all tubes/plates shortly at low speed.
- Place tubes/plates into the real-time PCR instrument and start the run according to the setup.

*** Description of the controls**

- Negative control: only master-mix
- Extraction control: the extraction is performed without the sample – components from used Prep Kit
- Positive Control: master-mix and within the kit's provided Positive Control

2.2 Interpretation of results

The evaluation has to be made according to the usual analysis program recommended by the real-time PCR instrument manufacturer.

The control reactions have to show the correct results.

RR-Soya DNA is detected in the FAM-channel, RR2Y-Soya DNA is detected in the VIC/HEX -channel, A2704-12-Soya DNA is detected in the ROX-channel and A5547-127-Soya DNA is detected in the Cy5-channel (see table).

A sample is stated **positive** for the respective parameter, if the sample DNA shows amplification in the respective channel.

A sample is stated **negative** for the respective parameter, if the sample DNA shows no amplification in the respective channel and if the external inhibition control of the sample is **positive** with a shift in Cp-value ≤ 2 compared to the negative control.

If the sample DNA in the external inhibition control shows **no amplification** or a shift in Cp-value > 2 compared to the negative control, it contains PCR inhibiting substances. A significant decrease in the fluorescence signal can also show the presence of PCR inhibiting substances. Under these circumstances DNA isolation and purification of the sample need to be improved. It is recommended to repeat the extraction process. Alternatively the DNA can be diluted (recommendation 1:2 in PCR-water) and analysed again for inhibition. Please note that the dilution factor also affects the detection limit of the specific PCR assay.

SureFood® GMO ID 4plex Soya II (100 rxn)

Art. No. S2162

March 2026

Result in the respective channel				Result	Interpretation
FAM channel RR-Soya	VIC/HEX channel RR2Y-Soya	ROX channel A2704-12-Soya	Cy5 channel A5547-127-Soya	external inhibition- control	
positive	negative	negative	negative	positive	RR-Soya DNA detected
negativ	positive	negative	negative	positive	RR2Y-Soya DNA detected
negativ	negative	positive	negative	positive	A2704-12-Soya DNA detected
negativ	negative	negative	positive	positive	A5547-127-Soya DNA detected
negativ	negative	negative	negative	positive	Negative, no target DNA detected
negativ	negative	negative	negative	negative	invalid

Note: The results displayed in the table above represent merely an example. Additional combinations are also possible.

The following table shows the specification ranges of the kit controls

	Specification range
Positive Control (FAM – RR-Soya)	$25 \leq Cp \leq 33$
Positive Control (VIC – RR2Y-Soya)	$25 \leq Cp \leq 33$
Positive Control (ROX – A2704-12-Soya)	$25 \leq Cp \leq 33$
Positive Control (CY5 – A5547-127-Soya)	$25 \leq Cp \leq 33$

3 Limitations of the method

- The presence of PCR inhibitors may cause invalid results.
- Extremely low levels of target below the limit of detection (LoD) may be detected, but results may not be reproducible.
- In highly processed samples, the detection limit may be shifted. Factors such as high pressures, mechanical stresses, chemical treatment, extreme temperatures and/or extreme pH values during manufacturing process – such as in canning production – can damage or degrade nucleic acids. This means that the sensitivity of the test kit may be reduced and not all original components may be detected.

4 Further Information

4.1 Product Information

- Detailed information about setup of several real-time PCR devices
(Download: www.congen.de/en/company/downloads)
- Product-related documents (Download: www.congen.de/en/eifu/)
- Validation Report upon request

4.2 Technical Support

For further questions please contact your distributor or send an e-mail to sales@r-biopharm.de.

4.3 Distribution and Ordering

R-Biopharm AG
An der neuen Bergstrasse 17,
64297 Darmstadt, Germany
Phone: +49 (0) 61 51 - 81 02-0
Fax: +49 (0) 61 51 - 81 02-20
E-Mail: orders@r-biopharm.de
www.r-biopharm.com