CONGEN

SureFood® GMO ID 4plex Soya IV

Art. No. S2165
100 rxn

User Manual

March 2025

 **Inhalt**

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Content

1	General Information	9
1.1	Description	9
1.2	Limit of Detection	9
1.3	DNA-preparation	10
1.4	Kit components and storage	10
1.5	Additionally required equipment and materials	10
1.6	Precautions for users	10
1.7	Setup.....	11
1.8	Detection channel Set-up	11
2	Qualitative Analysis	12
2.1	Protocol	12
2.1.1	Preparation of the master-mix	12
2.1.2	Preparation of the real-time PCR-mix	12
2.2	Interpretation of results	13
3	Limitations of the method	14
4	Further Information	14
4.1	Product Information	14
4.2	Technical Support	14
4.3	Distribution and Ordering	14

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1 General Information

1.1 Description

The SureFood® GMO ID 4plex Soya IV is a real-time PCR for the direct, qualitative detection and differentiation of following specific genetically modified soya DNA sequences:

- MON87705 soya (OECD unique identifier MON-87705-6)
- DAS-81419-2 soya (OECD unique identifier DAS-81419-2)
- MON87751 soya (OECD unique identifier MON-87751-7)
- SYHT0H2 soya (OECD unique identifier SYN-000H2-5)

This kit can be used for the identification of genetically modified organisms (GMOs) in food, feed and seeds.

The detections are according to the validated method of the European Commission.

The real-time PCR assay can be performed with commonly used real-time PCR instruments, equipped for detection of four fluorescence emissions at the channels FAM, VIC/HEX, ROX and Cy5 at the same time. The technical verification of instruments was performed on Roche LightCycler® 480 II, Qiagen Rotor-Gene Q, Bio-Rad CFX96 Dx and R-Biopharm RIDA®CYCLER.

1.2 Limit of Detection

The SureFood® GMO ID 4plex Soya IV real-time PCR has a limit of detection of ≤ 5 DNA copies. This is equivalent to approx. 0.01 % for unprocessed soybean.

The assay limit of detection depends on sample matrix, processing grade, DNA preparation and DNA content.

The SureFood® PCR systems are very sensitive and therefore even a small amount of target DNA is sufficient for a successful analysis. The concentration of total DNA in the sample does not allow a conclusion on the quantity and quality of the target DNA.

Note: Inconsistent mixing ratios* may cause a loss of sensitivity in the low concentration channel in mixed samples especially with high amplicon concentrations (Cp value < 20).

* e.g. 99.9 MON87705 soya and 0.1 % DAS-81419-2 soya

1.3 DNA-preparation

For DNA-preparation of raw material the use of SureFood® PREP Basic (Art. No. S1052) and for highly processed food and feed the use of SureFood® PREP Advanced (Art. No. S1053) is recommended.

1.4 Kit components and storage

Kit Code	Reagent	Amount	Lid Color
1	Reaction Mix	2 x 1050 µl	Yellow
2	Taq Polymerase	1 x 80 µl	Dark Red
3	Positive Control	1 x 190 µl	Light Blue

Store all reagents at -28°C to -16°C and protected from light. The Taq Polymerase can be stored at +2 to +8°C for multiple uses on the same day.

Note: The Taq Polymerase may be in a frozen or unfrozen state. This does not affect the quality of the Taq Polymerase or the performance of the real-time PCR.

1.5 Additionally required equipment and materials

- DNA-Extraction kit
(e.g. SureFood® PREP Basic Art. No. S1052 / SureFood® PREP Advanced Art. No. S1053)
- real-time PCR instrument with four detection channels (510 nm, 580 nm, 610 nm and 660 nm)
- real-time PCR consumable (plates, tubes, foils, caps)
- pipettes with filter tips
- powder-free disposable gloves
- Vortex mixer
- micro centrifuge with a rotor for the reaction tubes

1.6 Precautions for users

- Extraction, PCR preparation and the PCR run should be separated in different rooms to avoid cross-contaminations.
- This test must only be performed by laboratory personnel trained in molecular biology methods.
- Strictly follow the working instructions.
- When handling samples, wear disposable gloves. After finishing the test, wash your hands.
- Do not smoke, eat or drink in areas where samples or test reagents are being used.
- Do not use the kit after the expiration date.

1.7 Setup

	Blockcycler & R-Biopharm RIDA®CYCLER	Rotorcycler
Initial Denaturation (HOLD) Cycles	5 min, 95°C 45	1 min, 95°C 45
Denaturation Annealing/Extension (CYCLE)	15 sec, 95°C 30 sec, 60°C	10 sec, 95°C 15 sec, 60°C
Temperature Transition Rate/ Ramp Rate	Maximum	Maximum

1.8 Detection channel Set-up

Real-time PCR device	Detection	Detection channel	Quencher	Note
Bio-Rad CFX96/Dx/Opus	MON87705 soya	FAM	+	
	DAS-81419-2 soya	VIC/HEX	+	
	MON87751 soya	ROX	+	
	SYHT0H2 soya	Cy5	+	
R-Biopharm RIDA®CYCLER	MON87705 soya	green	+	Ignore cycles before, if there is a significant deviation in the baseline at the start of the run. Please see page 45 of the cycler operating instructions, section 12.1.2 Cycling analysis parameter.
	DAS-81419-2 soya	yellow	+	
	MON87751 soya	orange	+	
	SYHT0H2 soya	red	+	
Qiagen Rotor- Gene Q	MON87705 soya	green	+	Note: Please use only 0.1 ml reaction tube. The gain settings must be set to 5 (factory default) for all channels.
	DAS-81419-2 soya	yellow	+	
	MON87751 soya	orange	+	
	SYHT0H2 soya	red	+	
Roche LightCycler® 480 II	MON87705 soya	465-510	+	The SureCC Color Compensation Kit I (Art. No. F4009) is required.
	DAS-81419-2 soya	533-580	+	
	MON87751 soya	533-610	+	
	SYHT0H2 soya	618-660	+	

2 Qualitative Analysis

2.1 Protocol

2.1.1 Preparation of the master-mix

Calculate the total number of reactions needed (samples and control reactions) for the specific PCR assay as well as for the inhibition control.

Recommended control reactions for the specific PCR assay: negative control, extraction control, Positive Control and an inhibition control per sample.

For the preparation of the inhibition control, it is recommended to use the SureFood® GMO Plant PLUS (Art. No. S2049) or the SureFood® GMO SCREEN 4plex 35S/NOS/FMV+IAC (Art. No. S2126).

Reactions needed for the qualitative MON87705 soya, DAS-81419-2 soya, MON87751 soya or SYHT0H2 soya detection:

3 reactions for controls (1x negative control, 1x extraction control, 1x Positive Control)

For each sample: at least 1 reaction for each sample DNA

It is also recommended to prepare the master-mix with 10 % additional volume in order to compensate reagent loss. Allow the reagents to thaw, mix and centrifuge before opening and use.

Example for the calculation and preparation of 10 reactions:

Components of the master-mix	Amount per reaction	10 reactions (with 10% excess)
Reaction Mix	19.3 µl	212.3 µl
Taq Polymerase	0.7 µl	7.7 µl
Total volume	20 µl	220 µl

Mix each master-mix well and centrifuge shortly before use.

2.1.2 Preparation of the real-time PCR-mix

- Pipette 20 µl of the master-mix into appropriate tubes/wells.
- Close the negative control (the negative control is ready for PCR without any addition).
- Pipette 5 µl of sample DNA into the designated tubes/wells and close them.
- Pipette 5 µl of Positive Control into the designated tubes/wells and close them.
- Centrifuge all tubes/plates shortly at low speed.
- Place tubes/plates into the real-time PCR instrument and start the run according to the setup.

2.2 Interpretation of results

The evaluation has to be made according to the usual analysis program recommended by the real-time PCR instrument manufacturer.

The control reactions have to show the correct results.

MON87705 soya DNA is detected in the FAM-channel, DAS-81419-2 soya DNA is detected in the VIC/HEX - channel, MON87751 soya DNA is detected in the ROX-channel and SYHT0H2 soya DNA is detected in the Cy5-channel (see table).

A sample is stated **positive** for the respective parameter, if the sample DNA shows amplification in the respective channel.

A sample is stated **negative** for the respective parameter, if the sample DNA shows no amplification in the respective channel and if the external inhibition control of the sample is **positive** with a shift in Cp-value ≤ 2 compared to the negative control.

If the sample DNA in the external inhibition control shows **no amplification** or a shift in Cp-value > 2 compared to the negative control, it contains PCR inhibiting substances. A significant decrease in the fluorescence signal can also show the presence of PCR inhibiting substances. Under these circumstances DNA isolation and purification of the sample need to be improved. Alternatively the DNA can be diluted (recommendation 1:2 in PCR-water) and analysed again for inhibition. Please note that the dilution factor also affects the detection limit of the specific MON87705 soya, DAS-81419-2 soya, MON87751 soya or SYHT0H2 soya PCR assay.

Result in the respective channel				Result	Interpretation
FAM channel MON87705 soya	VIC/HEX channel DAS-81419-2 soya	ROX channel MON87751 soya	Cy5 channel SYHT0H2 soya	external inhibition- control	
positive	negative	negative	negative	positive	MON87705 soya DNA detected
negativ	positive	negative	negative	positive	DAS-81419-2 soya DNA detected
negativ	negative	positive	negative	positive	MON87751 soya DNA detected
negativ	negative	negative	positive	positive	SYHT0H2 soya DNA detected
negativ	negative	negative	negative	positive	Negative, target DNA is not detected
negativ	negative	negative	negative	negative	invalid

Note: The results displayed in the table above represent merely an example. Additional combinations are also possible.

3 Limitations of the method

- The presence of PCR inhibitors may cause invalid results.
- Extremely low levels of target below the limit of detection (LoD) may be detected, but results may not be reproducible.
- In highly processed samples, the limit of detection may be shifted. Factors such as high pressures, mechanical stresses, chemical treatments, extreme temperatures and/or extreme pH values during manufacturing process – such as in canning production – can damage or degrade nucleic acids. This means that the sensitivity of the test kit may be reduced and not all original components may be detected.

4 Further Information

4.1 Product Information

- Detailed information about setup of several real-time PCR devices
(Download: www.congen.de/en/company/downloads)
- Product-related documents (Download: www.congen.de/en/eifu/)
- Validation Report upon request

4.2 Technical Support

For further questions please contact your distributor or send an e-mail to sales@r-biopharm.de.

4.3 Distribution and Ordering

R-Biopharm AG
An der neuen Bergstrasse 17,
64297 Darmstadt, Germany
Phone: +49 (0) 61 51 - 81 02-0
Fax: +49 (0) 61 51 - 81 02-20
E-Mail: orders@r-biopharm.de
www.r-biopharm.com