CONGEN

SureFood® QUANTARD Allergen 40

Art. No. S3301
100 rxn

User Manual

October 2025

 Inhalt

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1 General Information

1.1 Description

Der SureFood® QUANTARD Allergen 40 is validated for quantitative determination of 1 mg to 400 mg allergenic substance / kg food sample.

The laboratory reference material SureFood® QUANTARD Allergen 40 is based on **corn flour** material. It contains all potentially allergenic food ingredients according to regulation (EU) No 1169/2011 (except for **lactose and sulphur dioxide**) in a concentration of **40 mg/kg**.

Content of allergenic substances:

Allergen	Material
Gluten containing cereals	Wheat flour type 405 (<i>Triticum</i> spp.)
Crustacean	Black Tiger shrimps (<i>Penaeus monodon</i>)
Molluscs	Sea scallop (<i>Placopecten magellanicus</i>)
Egg	Egg powder (Hen)
Fish	Humpback salmon (<i>Oncorhynchus gorbuscha</i>)
Peanut	Peanut (<i>Arachis hypogaea</i>)
Soy	Soybean (<i>Glycine max</i>)
Milk	Milkpowder (cow)
Edible Nut	Hazelnut (<i>Corylus avellana</i>), Almond (<i>Prunus dulcis</i>), Walnut (<i>Juglans regia</i>), Cashew (<i>Anacardium occidentale</i>), Pecan (<i>Carya illinoensis</i>), Brazil nut (<i>Bertholletia excelsa</i>), Pistachio (<i>Pistacia vera</i>), Macadamia (<i>Macadamia ternifolia</i>)
Celery	Celery seeds (<i>Apium graveolens</i>)
Mustard	Mustard seeds, yellow (<i>Sinapis alba</i>)
Sesame	Sesam seeds (<i>Sesamum indicum</i>)
Lupin	Lupin seeds (<i>Lupinus hartwegii</i>)
Buckwheat	Buckwheat (<i>Fagopyrum esculentum</i>)

The laboratory reference material SureFood® QUANTARD Allergen 40 is validated in combination with the DNA extraction method SureFood® PREP Advanced, protocol 1, and the SureFood® ALLERGEN real-time PCR detection methods.

Further information about the quantification of allergenic substances is given in the manual of the SureFood® ALLERGEN real-time PCR products.

In case of using alternative methods the sample weight and extraction method in combination with the detection system have to be validated.

1.2 DNA-preparation

SureFood® PREP Advanced (Art. No. S1053) is recommended for the DNA preparation.

As quantification is performed using an external reference, both the SureFood® QUANTARD Allergen 40 standard and the sample must undergo identical extraction procedures. The SureFood® QUANTARD Allergen 40 standard should be extracted in duplicate alongside the test samples to ensure accuracy and reproducibility.

Lysis, binding, and DNA purification steps must be conducted strictly in accordance with the established protocol.

1.3 Kit components and storage

Reagent	Amount
QUANTARD Allergen 40	1 x 2,0 g

Store all reagents dry at –28°C to – 16°C and protected from light.

Note: To avoid moisture absorption the unopened laboratory reference material should thaw at room temperature for at least one hour before use.

1.4 Additionally required equipment and materials

- DNA-Extraction kit (e.g. SureFood® PREP Advanced, Art. No.: S1053)
- Real-time PCR methods (SureFood® ALLERGEN series)

2 Further Information

2.1 Product Information

- Product-related documents (Download: www.congen.de/en/eifu/)

2.2 Technical Support

For further questions please contact your distributor or send an e-mail to sales@r-biopharm.de.

2.3 Distribution and Ordering

R-Biopharm AG
An der neuen Bergstrasse 17,
64297 Darmstadt, Germany
Phone: +49 (0) 61 51 - 81 02-0
Fax: +49 (0) 61 51 - 81 02-20
E-Mail: orders@r-biopharm.de
www.r-biopharm.com