CONGEN

SureFood® QUANTARD Allergen 40

Art. No. S3301

User Manual

February 2024

 **Inhalt**

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1 General Information

1.1 Description

The SureFood® QUANTARD Allergen 40 is validated for quantitative determination of 1 mg to 400 mg allergenic substance / kg food sample.

The laboratory reference material SureFood® QUANTARD Allergen 40 is based on **corn flour** material. It contains all potentially allergenic food ingredients according to regulation (EU) No 1169/2011 (except for **lactose and sulphur dioxide**) in a concentration of **40 mg/kg**.

Content of allergenic substances:

Allergen	Material
Gluten containing cereals	Wheat flour type 405 (<i>Triticum</i> spp.)
Crustacean	Black Tiger shrimps (<i>Penaeus monodon</i>), peeled, glazed, blanched, raw
Molluscs	Sea scallop (<i>Placopecten magellanicus</i>) *
Egg	Egg white powder , hen (<i>Gallus gallus</i>)
Fish	Humpback salmon filet (<i>Oncorhynchus gorbuscha</i>), untreated
Peanut	Peanut kernels (<i>Arachis hypogaea</i>), untreated, peeled
Soy	Soybean (<i>Glycine max</i>)
Milk	Milkpowder, cow (<i>Bos taurus</i>)
Edible Nut	Hazelnut kernels (<i>Corylus avellana</i>), without skin Almond (<i>Prunus dulcis</i>), whole, untreated, with skin Walnut kernels (<i>Juglans regia</i>), untreated, with skin Cashew kernels (<i>Anacardium occidentale</i>), untreated, without skin Pecan nuts (<i>Carya illinoensis</i>), untreated, with skin Brazil nut (<i>Bertholletia excelsa</i>), peeled, without skin Pistachio kernels (<i>Pistacia vera</i>), peeled, without skin Macadamia nut (<i>Macadamia ternifolia</i>), roasted in the shell, peeled
Celery	Celery seeds (<i>Apium graveolens</i>)
Mustard	Mustard seeds, yellow (<i>Sinapis alba</i>)
Sesame	Sesam seeds (<i>Sesamum indicum</i>), peeled
Lupin	Lupin (<i>Lupinus albus</i>), seed of white lupin, peeled
Buckwheat	Buckwheat (<i>Fagopyrum esculentum</i>)

***Note:** To be used from SureFood® ALLERGEN Molluscs Kit, Lot. 24493 (User Manual version January 2024).

The laboratory reference material SureFood® QUANTARD Allergen 40 is verified in combination with the DNA extraction method SureFood® PREP Advanced (S1053), protocol 1, and the SureFood® ALLERGEN real-time PCR detection methods.

Further information about the quantification of allergenic substances is given in the manual of the SureFood® ALLERGEN real-time PCR products.

In case of using alternative methods the sample weight and extraction method in combination with the detection system have to be verified.

1.2 DNA-preparation

For DNA preparation the use of SureFood® PREP Advanced Protocol 1 (Art. No. S1053) is recommended.

Because it is a quantitative determination with an external reference material the DNA preparation of sample and reference material have to be performed in the same way.

The reference material is supposed to be prepared parallel to the samples.

1.3 Kit components and storage

Reagent	Amount
QUANTARD Allergen 40	1 x 2.0 g

Store all reagents dry at –28°C to –16°C and protected from light.

Note: To avoid moisture absorption thaws the unopened laboratory reference material at room temperature for at least one hour before use.

1.4 Additionally required equipment and materials

- DNA-Extraction kit (e.g. SureFood® PREP Advanced, Art. No.: S1053)
- Real-time PCR methods (SureFood® ALLERGEN series)

2 Further Information

2.1 Technical Support

For further questions please contact your distributor or send an e-mail to sales@r-biopharm.de.

2.2 Distribution and Ordering

R-Biopharm AG
An der neuen Bergstrasse 17,
64297 Darmstadt, Germany
Phone: +49 (0) 61 51 - 81 02-0
Fax: +49 (0) 61 51 - 81 02-20
E-Mail: orders@r-biopharm.de
www.r-biopharm.com