CONGEN

SureFood® QUANTARD

Allergen 1 ppm

Art. No. S3305

User Manual

April 2026

 **Inhalt**

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1 General Information

1.1 Description

The SureFood® QUANTARD Allergen 1 ppm is a laboratory reference material for the determination of the limit of detection and the limit of quantification of the SureFood® ALLERGEN Kits.

The reference material SureFood® QUANTARD Allergen 1 ppm is based on **corn flour** material. It contains all potentially allergenic food ingredients according to regulation (EU) No 1169/2011 (except for **lactose and sulphur dioxide**) in a concentration of **1 mg/kg**.

Content of allergenic substances:

Allergen	Material
Buckwheat	Buckwheat (<i>Fagopyrum esculentum</i>)
Celery	Celery seeds (<i>Apium graveolens</i>)
Crustacean	Black Tiger shrimps (<i>Penaeus monodon</i>), peeled, glazed, blistered, raw, deveined
Edible Nut	Hazelnut kernels (<i>Corylus avellana</i>), without skin Almond (<i>Prunus dulcis</i>), whole, untreated, with skin Walnut kernels (<i>Juglans regia</i>), untreated, with skin Cashew kernels (<i>Anacardium occidentale</i>), untreated, without skin Pecan nuts (<i>Carya illinoensis</i>), untreated, with skin Brazil nut (<i>Bertholletia excelsa</i>), peeled, without skin Pistachio kernels (<i>Pistacia vera</i>), peeled, without skin Macadamia nut (<i>Macadamia ternifolia</i>), roasted in the shell, peeled
Egg	Egg white powder, hen (<i>Gallus gallus</i>)
Fish	Humpback salmon filet (<i>Oncorhynchus gorbuscha</i>), untreated
Gluten containing cereals	Wheat flour type 405 (<i>Triticum</i> spp.)
Lupin	Lupin (<i>Lupinus albus</i>), seed of white lupin, without skin
Milk	Milkpowder, cow (<i>Bos taurus</i>)
Mustard	Mustard seeds, yellow (<i>Sinapis alba</i>)
Molluscs	Sea scallop (<i>Placopecten magellanicus</i>), raw, glazed
Peanut	Peanut kernels (<i>Arachis hypogaea</i>), untreated, without skin
Sesame	Sesam seeds (<i>Sesamum indicum</i>), peeled
Soy	Soybean (<i>Glycine max</i>)

The reference material SureFood® QUANTARD Allergen 1 ppm is verified in combination with the DNA extraction method SureFood® PREP Advanced, protocol 1, and the SureFood® ALLERGEN real-time PCR detection methods.

In case of using alternative methods the sample weight and extraction method in combination with the detection system have to be validated.

1.2 DNA-preparation

For DNA preparation the use of SureFood® PREP Advanced Protocol 1 (Art. No. S1053) is recommended.

1.3 Kit components and storage

Reagent	Amount
QUANTARD Allergen 1 ppm	1 x 2.0 g

Store all reagents dry at -28 to -16°C and protected from light.

Note: To avoid moisture absorption thaws the unopened laboratory reference material at room temperature for at least one hour before use.

1.4 Additionally required equipment and materials

- DNA-Extraction kit (e.g. SureFood® PREP Advanced, Art. No. S1053)
- Real-time PCR methods (SureFood® ALLERGEN series)

2 Further Information

2.1 Product Information

- Product-related documents (Download: www.congen.de/en/eifu/)

2.2 Technical Support

For further questions please contact your distributor or send an e-mail to sales@r-biopharm.de.

2.3 Distribution and Ordering

R-Biopharm AG
An der neuen Bergstrasse 17,
64297 Darmstadt, Germany
Phone: +49 (0) 61 51 - 81 02-0
Fax: +49 (0) 61 51 - 81 02-20
E-Mail: orders@r-biopharm.de
www.r-biopharm.com