**CONGEN**

# **SureFood® ALLERGEN 4plex EU NUTS**

Art. No. S3404  
100 rxn

## **User Manual**

**December 2024**

 **Inhalt**

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## **Content**

1	General Information .....	10
1.1	Description .....	10
1.2	Limit of Detection .....	10
1.3	DNA preparation .....	11
1.4	Kit components and storage .....	11
1.5	Additionally required equipment and materials .....	11
1.6	Setup.....	11
1.7	Detection channel Set-up .....	12
2	Qualitative Analysis .....	13
2.1	Protocol .....	13
2.1.1	Preparation of the master-mix .....	13
2.1.2	Preparation of the real-time PCR-mix .....	13
2.2	Interpretation of results .....	14
3	Limitations of the method .....	15
4	Further Information .....	15
4.1	Product Information .....	15
4.2	Technical Support .....	15
4.3	Distribution and Ordering .....	15

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## **1 General Information**

### **1.1 Description**

The SureFood® ALLERGEN 4plex EU NUTS is a real-time PCR for the direct, qualitative detection and differentiation of specific almond (*Prunus dulcis*), cashew (*Anacardium occidentale*), pistachio (*Pistacia vera*), peanut (*Arachis hypogaea*), hazelnut (*Corylus avellana*), walnut (*Juglans regia*), pecan (*Carya illinoensis*), brazil nut (*Bertholletia excelsa*) and macadamia (*Macadamia tetraphylla*) DNA sequences according to directive (EC) 1169/2011 in food.

Each reaction contains an internal amplification control (IAC). If the DNA contains PCR inhibiting substances, the signal of the amplification control will be affected or the amplification will be suppressed. Examples for PCR inhibiting substances are alcohols (e.g. ethanol, isopropanol), surfactants (e.g. CTAB, SDS, Triton X100) and salts (e.g. sodium chloride). In addition spices, herbs, algae, cocoa and further sample matrices might have PCR inhibiting effects.

The real-time PCR assay can be performed with commonly used real-time PCR instruments, equipped for detection of four fluorescence emissions at the channels FAM, VIC/HEX, ROX and Cy5 at the same time. . The technical verification of instruments was performed on Qiagen Rotor-Gene Q, R-Biopharm RIDA®CYCLER, Bio-Rad CFX96Dx, Roche LightCycler® 480 II and Agilent AriaDx.

### **1.2 Limit of Detection**

The SureFood® ALLERGEN 4plex EU NUTS PCR has a limit of detection of  $\leq 1$  mg / kg using SureFood® PREP Advanced, protocol 1.

The limit of detection of the complete method (DNA extraction and real-time PCR) depends on sample matrix, processing grade, DNA preparation and DNA content.

The SureFood® PCR systems are very sensitive and therefore even a small amount of target DNA is sufficient for a successful analysis. The concentration of total DNA in the sample does not allow a conclusion on the quantity and quality of the target DNA.

**Note:** Inconsistent mixing ratios may cause a loss of sensitivity in the low concentration channel in mixed samples especially with high amplicon concentrations (Cp value < 20).

**1.3 DNA preparation**

For DNA preparation the use of SureFood® PREP Advanced (Art. No. S1053), protocol 1 is recommended.

**1.4 Kit components and storage**

Kit Code	Reagent	Amount	Lid Color
1	Reaction Mix	2 x 1050 µl	Yellow
2	Taq Polymerase	1 x 80 µl	Dark Red
3	Positive Control	1 x 200 µl	Light Blue

**Store all reagents at -28 to -16°C and protected from light. The Taq Polymerase can be stored at +2 to +8°C for multiple uses on the same day.**

**Note: The Taq Polymerase may be in a frozen or unfrozen state. This does not affect the quality of the Taq Polymerase or the performance of the real-time PCR.**

**1.5 Additionally required equipment and materials**

- DNA extraction kit (e.g. SureFood® PREP Advanced Art. No. S1053)
- real-time PCR instrument with four detection channels (510 nm, 580 nm, 610 nm and 660 nm)
- real-time PCR consumables (plates, tubes, foils, caps)
- pipettes with filter tips
- powder-free disposable gloves
- Vortex mixer
- micro centrifuge with a rotor for the reaction tubes

**1.6 Setup**

	<b>Blockcycler &amp; R-Biopharm RIDA®CYCLER</b>	<b>Rotorcycler</b>
Initial Denaturation (HOLD)	5 min, 95°C	1 min, 95°C
Cycles	35	35
Denaturation	15 sec, 95°C	10 sec, 95°C
Annealing/Extension (CYCLE)	30 sec, 60°C	15 sec, 60°C
Temperature Transition Rate/ Ramp Rate	Maximum	Maximum

**1.7 Detection channel Set-up**

<b>Real-time PCR device</b>	<b>Detection</b>	<b>Detection channel</b>	<b>Quencher</b>	<b>Note</b>
<b>Agilent AriaDx/Mx</b>	almond / cashew / pistachio	FAM	+	
	IAC	HEX	+	
	peanut / hazelnut / walnut	ROX	+	
	pecan / brazil nut / macadamia	Cy5	+	
<b>Bio-Rad CFX96/Dx/ Opus</b>	almond / cashew / pistachio	FAM	+	
	IAC	VIC/HEX	+	
	peanut / hazelnut / walnut	ROX	+	
	pecan / brazil nut / macadamia	Cy5	+	
<b>R-Biopharm RIDA®CYCLER</b>	almond / cashew / pistachio	green	+	<b>Ignore cycles before</b> , if there is a significant deviation in the baseline at the start of the run. Please see page 45 of the cyclers operating instructions, section 12.1.2 Cycling analysis parameter
	IAC	yellow	+	
	peanut / hazelnut / walnut	orange	+	
	pecan / brazil nut / macadamia	red	+	
<b>Qiagen Rotor- Gene Q</b>	almond / cashew / pistachio	green	+	<b>Note:</b> Please use only 0.1 ml reaction tubes. The gain settings must be set to 5 (factory default) for all channels.
	IAC	yellow	+	
	peanut / hazelnut / walnut	orange	+	
	pecan / brazil nut / macadamia	red	+	
<b>Roche LightCycler® 480 II</b>	almond / cashew / pistachio	465-510	+	The SureCC Color Compensation Kit I (Art. No. F4009) is required.
	IAC	533-580	+	
	peanut / hazelnut / walnut	533-610	+	
	pecan / brazil nut / macadamia	618-660	+	

## 2 Qualitative Analysis

### 2.1 Protocol

#### 2.1.1 Preparation of the master-mix

Calculate the total number of reactions needed (samples and control reactions) for the specific PCR assay. Recommended control reactions for the specific PCR assay: negative control, extraction control, Positive Control.

The reaction mix contains an internal amplification control (Inhibitioncontrol) per reaction.

#### Reactions needed for the qualitative EU Nuts detection:

3 reactions for controls (1x negative control, 1x extraction control, 1x Positive Control)

For each sample: at least 1 reaction for each sample DNA

It is also recommended to prepare the master-mix with 10% additional volume in order to compensate reagent loss. Allow the reagents to thaw, mix and centrifuge before opening and use.

#### Example for the calculation and preparation of 10 reactions:

Components of the master-mix	Amount per reaction	10 reactions (with 10% excess)
Reaction Mix	19.3 µl	212.3 µl
Taq Polymerase	0.7 µl	7.7 µl
<b>Total volume</b>	<b>20 µl</b>	<b>220 µl</b>

**Mix each master-mix well and centrifuge shortly before use.**

#### 2.1.2 Preparation of the real-time PCR-mix

- Pipette 20 µl of the master-mix into appropriate tubes/wells.
- Close the negative control (the negative control is ready for PCR without any addition).
- Pipette 5 µl of sample DNA into the designated tubes/wells and close them.
- Pipette 5 µl of Positive Control into the designated tubes/wells and close them.
- Centrifuge all tubes/plates shortly at low speed.
- Place tubes/plates into the real-time PCR instrument and start the run according to the setup.

**2.2 Interpretation of results**

The evaluation has to be made according to the usual analysis program recommended by the real-time PCR instrument manufacturer.

The control reactions have to show the correct results.

Almond, cashew and/or pistachio DNA is detected in the FAM-channel, peanut, hazelnut and/or walnut DNA is detected in the ROX-channel and pecan, brazil nut and/or macadamia DNA is detected in the Cy5-channel (see table). In the VIC/HEX-channel the amplification control (IAC) is detected.

A sample is stated **positive** for the respective parameter, if the sample DNA shows amplification in the respective channel. High amplicon concentrations can result in a weak or absent signal of the IAC.

A Cp value for the IAC is not needed to obtain a positive result of the positive control.

A sample is stated **negative** for the respective parameter, if the sample DNA shows no amplification in the respective channel and if the internal control (VIC/HEX-channel) of the sample is **positive** with a shift in Cp-value  $\leq 2$  compared to the negative control.

If the sample DNA in the VIC/HEX-Channel shows **no amplification** or a shift in Cp-value  $> 2$  compared to the negative control, it contains PCR inhibiting substances. A significant decrease in the fluorescence signal can also show the presence of PCR inhibiting substances. Under these circumstances DNA isolation and purification of the sample need to be improved. Alternatively the DNA can be diluted (recommendation 1:2 in PCR-water) and analysed again for inhibition. Please note that the dilution factor also affects the detection limit of the specific PCR assay.

Result in the respective channel				VIC/HEX channel IAC	Interpretation
FAM channel almond / cashew / pistachio	ROX channel peanut / hazelnut / walnut	Cy5 channel pecan / brazil nut / macadamia			
<b>positive</b>	negative	negative	<b>positive/negative</b>	almond-, cashew- and/or pistachio-DNA detected	
negative	<b>positive</b>	negative	<b>positive/negative</b>	peanut-, hazelnut- and/or walnut-DNA detected	
negative	negative	<b>positive</b>	<b>positive/negative</b>	pecan-, brazil- nut and/or macadamia-DNA detected	
negative	negative	negative	<b>positive</b>	negative, target- DNA is not detected	
negative	negative	negative	negative	invalid	

**Note: The results displayed in the table above represent merely an example. Additional combinations are also possible.**

### **3 Limitations of the method**

- The presence of PCR inhibitors may cause invalid results.
- Extremely low levels of target below the limit of detection (LoD) may be detected, but results may not be reproducible.
- Samples tested negative still could contain an allergen contamination below the limit of detection of the assay, or they might contain other allergen components like proteins or lipids for example.
- The pecan detection system shows 100% cross reactivity to shagbark hickory (*Carya ovata*).

### **4 Further Information**

#### **4.1 Product Information**

- Detailed information about setup of several real-time PCR devices (Download: [www.congen.de/en/downloads](http://www.congen.de/en/downloads))
- Product-related documents (Download: [www.congen.de/en/eifu/](http://www.congen.de/en/eifu/))
- Validation Report upon request

#### **4.2 Technical Support**

For further questions please contact your distributor or send an e-mail to [sales@r-biopharm.de](mailto:sales@r-biopharm.de).

#### **4.3 Distribution and Ordering**

R-Biopharm AG  
An der neuen Bergstrasse 17,  
64297 Darmstadt, Germany  
Phone: +49 (0) 61 51 - 81 02-0  
Fax: +49 (0) 61 51 - 81 02-20  
E-Mail: [orders@r-biopharm.de](mailto:orders@r-biopharm.de)  
[www.r-biopharm.com](http://www.r-biopharm.com)

r-biopharm®