**CONGEN**

# **SureFood® ALLERGEN Lupin**

Art. No. S3611  
100 rxn

## **User Manual**

**August 2025**

 **Inhalt**

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## Content

|       |                                                      |    |
|-------|------------------------------------------------------|----|
| 1     | General Information .....                            | 14 |
| 1.1   | Description .....                                    | 14 |
| 1.2   | Limit of Detection and Limit of Quantification ..... | 14 |
| 1.3   | DNA preparation .....                                | 15 |
| 1.4   | Kit components and storage .....                     | 15 |
| 1.5   | Additionally required equipment and materials .....  | 15 |
| 1.6   | Precautions for users .....                          | 15 |
| 1.7   | Setup.....                                           | 16 |
| 1.8   | Detection channel Set-up .....                       | 16 |
| 2     | Qualitative Analysis .....                           | 18 |
| 2.1   | Protocol .....                                       | 18 |
| 2.1.1 | Preparation of the master-mix .....                  | 18 |
| 2.1.2 | Preparation of the real-time PCR-mix .....           | 18 |
| 2.2   | Interpretation of results .....                      | 19 |
| 3     | Quantitative Analysis .....                          | 20 |
| 3.1   | Protocol .....                                       | 20 |
| 3.1.1 | Preparation of the master-mix .....                  | 20 |
| 3.1.2 | Preparation of the standard DNA dilutions .....      | 21 |
| 3.1.3 | Preparation of the real-time PCR-mix .....           | 21 |
| 3.2   | Interpretation of results .....                      | 22 |
| 4     | Limitations of the method .....                      | 24 |
| 5     | Further Information .....                            | 24 |
| 5.1   | Product Information .....                            | 24 |
| 5.2   | Technical Support .....                              | 24 |
| 5.3   | Distribution and Ordering .....                      | 24 |

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## 1 General Information

### 1.1 Description

The SureFood® ALLERGEN Lupin is a real-time PCR for the direct, qualitative and / or quantitative detection of specific lupin (*Lupinus* spp.) DNA sequences according to directive (EC) 1169/2011 in food and feed.

Each reaction contains an internal amplification control (IAC). If the DNA contains PCR inhibiting substances, the signal of the amplification control will be affected, or the amplification will be suppressed. Examples for PCR inhibiting substances are alcohols (e.g. ethanol, isopropanol), surfactants (e.g. CTAB, SDS, Triton X100) and salts (e.g. sodium chloride). In addition, spices, herbs, algae, cocoa and further sample matrices might have PCR inhibiting effects.

For the quantitative determination the use of the laboratory reference material SureFood® QUANTARD Allergen 40 (Art. No. S3301) containing 40 mg lupin / kg food sample and a detection system including a standard series is required.

The real-time PCR assay can be performed with commonly used real-time PCR instruments, equipped for detection of two fluorescence emissions at the channels FAM and VIC/HEX at the same time. The internal technical verification of instruments was performed on Agilent Mx3005P, Agilent AriaDx, Bio-Rad CFX96, Applied Biosystems 7500, Bio Molecular Systems MIC, Qiagen Rotor-Gene Q, Roche cobas® z 480 Analyzer and Roche LightCycler® 480 II.

In general this procedure is applicable for all homogenous food samples. The quantitative determination is inapplicable for swabs, sponges and inhomogeneous fluids/liquids.

The method is verified for quantitative determination of 1 mg to 400 mg allergenic substance / kg food sample using the reference material SureFood® QUANTARD Allergen 40 containing 40 mg allergenic substance / kg food sample.

### 1.2 Limit of Detection and Limit of Quantification

The SureFood® ALLERGEN Lupin PCR has a limit of detection of  $\leq 0.4$  mg / kg using SureFood® PREP Advanced, protocol 1.

The limit of quantification is 1 mg / kg using SureFood® PREP Advanced, protocol 1.

The limit of detection and quantification were determined using the SureFood® QUANTARD Allergen 40 (matrix: corn flour). The limit of detection and quantification of the complete method (DNA extraction and real-time PCR) depends on sample matrix, processing grade, DNA preparation and DNA content.

The SureFood® PCR systems are very sensitive and therefore even a small amount of target DNA is sufficient for a successful analysis. The concentration of total DNA in the sample does not allow a conclusion on the quantity and quality of the target DNA.

## 1.3 DNA preparation

For DNA preparation the use of SureFood® PREP Advanced (Art. No. S1053) or SureFast® Mag PREP Food (Art. No. F1060) is recommended.

## 1.4 Kit components and storage

| Kit Code | Reagent          | Amount      | Lid Color  |
|----------|------------------|-------------|------------|
| 1        | Reaction Mix     | 2 x 1050 µl | Yellow     |
| 2        | Taq Polymerase   | 1 x 80 µl   | Dark Red   |
| 3        | Positive Control | 1 x 200 µl  | Light Blue |
| 4        | Dilution Buffer  | 1 x 1600 µl | White      |
| 5        | Standard DNA     | 1 x 200 µl  | Dark Blue  |

**Store all reagents at -28 to -16°C and protected from light. The Taq Polymerase can be stored at +2 to +8°C for multiple uses on the same day.**

**Note: The Taq Polymerase may be in a frozen or unfrozen state. This does not affect the quality of the Taq Polymerase or the performance of the real-time PCR.**

## 1.5 Additionally required equipment and materials

- DNA extraction kit  
(e.g. SureFood® PREP Advanced Art. No. S1053 or SureFast® Mag PREP Food Art. No. F1060)
- reference material for quantification (SureFood® QUANTARD Allergen 40, Art. No. S3301)
- real-time PCR instrument with two detection channels (510 nm and 580 nm)
- real-time PCR consumables (plates, tubes, foils, caps)
- pipettes with filter tips
- powder-free disposable gloves
- Vortex mixer
- micro centrifuge with a rotor for the reaction tubes

## 1.6 Precautions for users

- Extraction, PCR preparation and the PCR run should be separated in different rooms to avoid cross-contaminations.
- This test must only be performed by laboratory personnel trained in molecular biology methods.
- Strictly follow the working instructions.
- When handling samples, wear disposable gloves. After finishing the test, wash your hands.
- Do not smoke, eat or drink in areas where samples or test reagents are being used.
- Do not use the kit after the expiration date.
- All reagents and materials used have to be disposed properly after use. Please refer to the relevant national regulation for disposal.

### 1.7 Setup

|                                        | <b>Blockcycler &amp;<br/>R-Biopharm RIDA®CYCLER</b> | <b>Rotorcycler</b> |
|----------------------------------------|-----------------------------------------------------|--------------------|
| Initial Denaturation (HOLD)            | 5 min, 95°C                                         | 1 min, 95°C        |
| Cycles                                 | 45                                                  | 45                 |
| Denaturation                           | 15 sec, 95°C                                        | 10 sec, 95°C       |
| Annealing/Extension (CYCLE)            | 30 sec, 60°C                                        | 15 sec, 60°C       |
| Temperature Transition Rate/ Ramp Rate | Maximum                                             | Maximum            |

### 1.8 Detection channel Set-up

| <b>Real-time PCR device</b>    | <b>Detection</b> | <b>Detection channel</b> | <b>Quencher</b> | <b>Note</b>                                                                                                                                                                                                   |
|--------------------------------|------------------|--------------------------|-----------------|---------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| <b>Agilent Mx3005P</b>         | Lupin            | FAM                      | +               |                                                                                                                                                                                                               |
|                                | IAC              | HEX                      | +               |                                                                                                                                                                                                               |
| <b>Agilent AriaMx / Dx</b>     | Lupin            | FAM                      | +               |                                                                                                                                                                                                               |
|                                | IAC              | HEX                      | +               |                                                                                                                                                                                                               |
| <b>Applied Biosystems 7500</b> | Lupin            | FAM                      | None            | Check the passive reference option ROX is none.                                                                                                                                                               |
|                                | IAC              | VIC                      | None            |                                                                                                                                                                                                               |
| <b>Bio-Rad CFX96/ Dx/ Opus</b> | Lupin            | FAM                      | +               | Baseline Settings:<br><ul style="list-style-type: none"> <li>• Baseline subtracted curve fit</li> <li>• Apply fluorescence drift correction</li> </ul>                                                        |
|                                | IAC              | VIC/HEX                  | +               |                                                                                                                                                                                                               |
| <b>Qiagen Rotor-Gene Q</b>     | Lupin            | green                    | +               | <b>Note:</b> Please use only 0.1 ml reaction tubes.<br>The gain settings must be set to 5 (factory default) for all channels.                                                                                 |
|                                | IAC              | yellow                   | +               |                                                                                                                                                                                                               |
| <b>R-Biopharm RIDA®CYCLER</b>  | Lupin            | green                    | +               | <b>Ignore cycles before,</b> if there is a significant deviation in the baseline at the start of the run. Please see page 45 of the cycler operating instructions, section 12.1.2 Cycling analysis parameter. |
|                                | IAC              | yellow                   | +               |                                                                                                                                                                                                               |

| Real-time PCR device        | Detection | Detection channel | Quencher | Note |
|-----------------------------|-----------|-------------------|----------|------|
| Roche LightCycler® 480 II   | Lupin     | 465-510           | +        |      |
|                             | IAC       | 533-580           | +        |      |
| Roche cobas® z 480 Analyzer | Lupin     | 465-510           | +        |      |
|                             | IAC       | 540-580           | +        |      |

## 2 Qualitative Analysis

### 2.1 Protocol

#### 2.1.1 Preparation of the master-mix

Calculate the total number of reactions needed (samples and control reactions) for the specific PCR assay.

Recommended control reactions for the specific PCR assay: negative control, extraction control and Positive Control.

The reaction mix contains an internal amplification control (inhibition control) per reaction.

#### Reactions needed for the qualitative detection of lupin:

3 reactions for controls\* (1x negative control, 1x extraction control, 1x Positive Control)

For each sample: at least 1 reaction for each sample DNA

It is also recommended to prepare the master-mix with 10% additional volume in order to compensate reagent loss. Allow the reagents to thaw, mix and centrifuge before opening and use.

#### Example for the calculation and preparation of 10 reactions:

| Components of the master-mix | Amount per reaction | 10 reactions (with 10% excess) |
|------------------------------|---------------------|--------------------------------|
| Reaction Mix                 | 19.3 µl             | 212.3 µl                       |
| Taq Polymerase               | 0.7 µl              | 7.7 µl                         |
| <b>Total volume</b>          | <b>20 µl</b>        | <b>220 µl</b>                  |

**Mix each master-mix well and centrifuge shortly before use.**

#### 2.1.2 Preparation of the real-time PCR-mix

- Pipette 20 µl of the master-mix into appropriate tubes/wells.
- Close the negative control.
- Pipette 5 µl of sample DNA into the designated tubes/wells and close them.
- Pipette 5 µl of Positive Control into the designated tubes/wells and close them.
- Centrifuge all tubes/plates shortly at low speed.
- Place tubes/plates into the real-time PCR instrument and start the run according to the setup.

#### \* Description of the controls

- Negative control: only master-mix
- Extraction control: the extraction is performed without the sample – components from used Prep Kit
- Positive Control: master-mix and within the kit's provided Positive Control

**2.2 Interpretation of results**

The evaluation must be made according to the usual analysis program recommended by the real-time PCR instrument manufacturer.

The control reactions must show the correct results.

DNA of lupin is detected in the FAM-channel. In the VIC/HEX-channel the internal amplification control (IAC) is detected.

A sample is stated **positive**, if the sample DNA shows amplification in the detection system with a Cp value of  $\leq 35$ . High amplicon concentrations can result in a weak or absent signal of the IAC.

A Cp value for the IAC is not needed to obtain a positive result of the Positive Control.

A sample is stated **negative**, if the sample DNA shows no amplification in the detection system or the obtained Cp value is  $> 35$ . The internal amplification control (inhibition control) of the sample must be **positive** (VIC/HEX) with a shift in Cp-Value  $\leq 2$  compared to the negative control.

If the sample DNA in the VIC/HEX-Channel shows **no amplification** or a shift in Cp-value  $> 2$  compared to the negative control, it contains PCR inhibiting substances. A significant decrease in the fluorescence signal can also show the presence of PCR inhibiting substances. Under these circumstances DNA isolation and purification of the sample need to be improved. Alternatively, the DNA can be diluted (recommendation 1:2 in PCR-water) and analysed again for inhibition. Please note that the dilution factor also affects the detection limit of the specific lupin PCR assay.

Alternatively, a self-selected reference sample at the detection limit can be used to assess the sample. This reference sample is used outside CONGEN's quality and evaluation guidelines.

**The following table shows the specification ranges of the kit controls**

|                                   | Specification range  | Internal amplification control (IAC) |
|-----------------------------------|----------------------|--------------------------------------|
| Negative control                  | negative             | $25 \leq Cp \leq 34$                 |
| Positive Control<br>(FAM – Lupin) | $25 \leq Cp \leq 33$ | not relevant                         |

### 3 Quantitative Analysis

#### 3.1 Protocol

##### 3.1.1 Preparation of the master-mix

Calculate the total number of reactions needed (samples and control reactions) for the specific PCR assay.

Recommended control reactions for the specific PCR assay: negative control, extraction control, standard curve as positive control and for quantification.

The reaction mix contains an internal amplification control (inhibition control) per reaction.

##### **Reactions needed for the quantitative detection of lupin:**

5 reactions for the standard curve\*

2 reactions for controls\* (1x negative control, 1x extraction control)

2 reactions for reference material (SureFood® QUANTARD Allergen 40 DNA)

For each sample: at least 1 reaction for each sample DNA

It is also recommended to prepare the master-mix with 10% additional volume in order to compensate reagent loss. Allow the reagents to thaw, mix and centrifuge before opening and use.

##### **Example for the calculation and preparation of 10 reactions:**

| Components of the master-mix | Amount per reaction | 10 reactions (with 10% excess) |
|------------------------------|---------------------|--------------------------------|
| Reaction Mix                 | 19.3 µl             | 212.3 µl                       |
| Taq Polymerase               | 0.7 µl              | 7.7 µl                         |
| <b>Total volume</b>          | <b>20 µl</b>        | <b>220 µl</b>                  |

**Mix each master-mix well and centrifuge shortly before use.**

##### **\*Description of the reactions**

- Negative control: only master-mix
- Extraction control: the extraction is performed without the sample – components from used Prep Kit
- Positive control: master-mix and within the kit's provided standards

**3.1.2 Preparation of the standard DNA dilutions**

Use the undiluted Standard DNA (**Code 5**) and dilute the Standard DNA (**Code 5**) in 1:10 steps in Dilution Buffer (**Code 4**) in order to prepare different concentrations. Prepare a dilution series of 4 steps. Prepare 4 reaction tubes (labelled S2 to S5) and add 45 µl Dilution Buffer each.

**The following procedure is recommended:**

| Standard | Dilution                               | Copy number per µl | Final copy number per reaction* |
|----------|----------------------------------------|--------------------|---------------------------------|
| S1       | Standard DNA undiluted                 | 100,000 copies     | 500,000 copies                  |
| S2       | 45 µl Dilution Buffer + 5 µl DNA of S1 | 10,000 copies      | 50,000 copies                   |
| S3       | 45 µl Dilution Buffer + 5 µl DNA of S2 | 1,000 copies       | 5,000 copies                    |
| S4       | 45 µl Dilution Buffer + 5 µl DNA of S3 | 100 copies         | 500 copies                      |
| S5       | 45 µl Dilution Buffer + 5 µl DNA of S4 | 10 copies          | 50 copies                       |

**\*Note:** 5 µl of standard DNA are used for each calibration point. The final copy number per reaction is to be entered in the analysis software of the real-time PCR detection system.

**3.1.3 Preparation of the real-time PCR-mix**

- Pipette 20 µl of the master-mix into appropriate tubes/wells.
- Close the negative control.
- Pipette 5 µl of sample DNA into the designated tubes/wells and close them.
- Pipette 5 µl of the SureFood® QUANTARD 40 Allergen DNA and the standard dilutions into the designated tubes/wells and close them.
- Centrifuge all tubes/plates shortly at low speed.
- Place tubes/plates into the real-time PCR instrument and start the run according to the setup.

**3.2 Interpretation of results**

The evaluation must be made according to the usual analysis program recommended by the real-time PCR instrument manufacturer.

The control reactions must show the correct results.

DNA of lupin is detected in the FAM-channel. In the VIC/HEX-channel the internal amplification control (IAC) is detected.

A sample is stated **positive** and can be **quantified**, if the sample DNA shows amplification in the detection system with a Cp value of  $\leq 35$ . High amplicon concentrations can result in a weak or absent signal of the IAC.

A Cp value for the internal amplification control is not needed to obtain a positive result of the Positive control / Standard DNA.

A sample is stated **negative**, if the sample DNA shows no amplification in the detection system or the obtained Cp value is  $> 35$ . The internal amplification control (inhibition control) of the sample has to be **positive** (VIC/HEX) with a shift in Cp-Value  $\leq 2$  compared to the negative control.

If the sample DNA in the VIC/HEX-Channel shows **no amplification** or a shift in Cp-value  $> 2$  compared to the negative control, it contains PCR inhibiting substances. A significant decrease in the fluorescence signal can also show the presence of PCR inhibiting substances. Under these circumstances DNA isolation and purification of the sample need to be improved. Alternatively, the DNA can be diluted (recommendation 1:2 in PCR-water) and analysed again for inhibition. Please note that the dilution factor also affects the detection limit of the specific lupin PCR assay.

For quantification mark the standards, controls and samples and make the evaluation according to the analysis program recommended by the real-time PCR instrument manufacturer. The value for the slope of the standard curve must be between -3.1 and -3.6 and the correlation coefficient  $R^2 > 0.98$ . In case of different values for the standard curve, it should not be used for calculation.

**The following table shows the specification ranges of the kit controls**

|                                       | Specification range  | Internal amplification control (IAC) |
|---------------------------------------|----------------------|--------------------------------------|
| Negative control                      | negative             | $25 \leq Cp \leq 34$                 |
| Standard DNA, undiluted (FAM – Lupin) | $16 \leq Cp \leq 24$ | not relevant                         |

For a final calculation of mg allergenic substance / kg food sample following formula is introduced:

$$X[mg/kg] = \frac{10^{\left(\frac{Cp_{sample}}{s}\right)}}{10^{\left(\frac{Cp_{Quantard}}{s}\right)}} * c_{Quantard}$$

s = slope | c = concentration | Cp<sub>Quantard</sub> = Ø of double assay

**Example:**

$$X[mg/kg] = \frac{10^{\left(\frac{28.5}{-3.251}\right)}}{10^{\left(\frac{29.2}{-3.251}\right)}} * 40 = 65.7 \text{ mg/kg}$$

Cp<sub>sample</sub> = 28.5 | Cp<sub>Quantard</sub> = Ø 29.2 | s = - 3.251 | c = 40 mg/kg

For this example an amount of **65.7 mg lupin / kg food sample** is calculated.

## 4 Limitations of the method

- The presence of PCR inhibitors may cause invalid results.
- Extremely low levels of target below the limit of detection (LoD) may be detected, but results may not be reproducible.
- Samples tested negative still could contain an allergen contamination below the limit of detection of the assay, or they might contain other allergen components like proteins or lipids for example.
- In highly processed samples, the detection limit may be shifted. Factors such as high pressures, mechanical stresses, chemical treatment, extreme temperatures and/or extreme pH values during manufacturing process – such as in canning production – can damage or degrade nucleic acids. This means that the sensitivity of the test kit may be reduced and not all original components may be detected.

## 5 Further Information

### 5.1 Product Information

- Microsoft Excel template of calculation and detailed information about setup of several real-time PCR devices (Download: [www.congen.de/en/downloads](http://www.congen.de/en/downloads))
- Product-related documents (Download: [www.congen.de/en/eifu/](http://www.congen.de/en/eifu/))
- Validation Report upon request

### 5.2 Technical Support

For further questions please contact your distributor or send an e-mail to [sales@r-biopharm.de](mailto:sales@r-biopharm.de).

### 5.3 Distribution and Ordering

R-Biopharm AG  
An der neuen Bergstrasse 17,  
64297 Darmstadt, Germany  
Phone: +49 (0) 61 51 - 81 02-0  
Fax: +49 (0) 61 51 - 81 02-20  
E-Mail: [orders@r-biopharm.de](mailto:orders@r-biopharm.de)  
[www.r-biopharm.com](http://www.r-biopharm.com)