CONGEN

SureFood® ANIMAL ID

Pork SENS PLUS

Art. No. S6017
100 rxn

User Manual

July 2025

 Inhalt

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1 General Information

1.1 Description

The SureFood® ANIMAL ID Pork SENS PLUS is a real-time PCR for the direct, qualitative detection of a specific pork (*Sus scrofa*) DNA sequence in food and feed.

Each reaction contains an internal amplification control (IAC). If the DNA contains PCR inhibiting substances, the signal of the amplification control will be affected or the amplification will be suppressed. Examples for PCR inhibiting substances are alcohols (e.g. ethanol, isopropanol), surfactants (e.g. CTAB, SDS, Triton X100) and salts (e.g. sodium chloride). In addition spices, herbs, algae, cocoa and further sample matrices might have PCR inhibiting effects.

The real-time PCR assay can be performed with commonly used real-time PCR instruments, equipped for detection of two fluorescence emissions at the channels FAM and VIC/HEX at the same time. The internal technical verification of instruments was performed on Roche LightCycler® 480 II, Roche cobas® z 480 Analyzer, Qiagen Rotor-Gene Q, Applied Biosystems 7500, Bio-Rad CFX96, R-Biopharm RIDA®CYCLER, Agilent AriaDx and Agilent Mx3005P.

1.2 Limit of Detection

The SureFood® ANIMAL ID Pork SENS PLUS real-time PCR is developed for the detection of pork DNA in muscle meat mixture at a relative amount of $\leq 0.0001\%$.

Due to the high sensitivity the method is suitable for the analysis of highly processed products like e.g. gelatin, sweets or meat bone meal.

The limit of detection of the complete method (DNA extraction and real-time PCR) depends on sample matrix, processing grade, DNA-preparation and DNA-content.

The SureFood® PCR systems are very sensitive and therefore even a small amount of target DNA is sufficient for a successful analysis. The concentration of total DNA in the sample does not allow a conclusion on the quantity and quality of the target DNA.

SureFood® ANIMAL ID Pork SENS PLUS (100 rxn)

Art. No. S6017

July 2025

1.3 DNA-preparation

For DNA-preparation of raw material the use of SureFood® PREP Basic (Art. No. S1052) and for highly processed food and feed the use of SureFood® PREP Advanced (Art. No. S1053) is recommended.

1.4 Kit components and storage

Kit Code	Reagent	Amount	Lid Color
1	Reaction Mix	2 x 1050 µl	Yellow
2	Taq Polymerase	1 x 11 µl	Red
3	Positive Control	1 x 200 µl	Light Blue

Store all reagents at -28 to -16°C and protected from light. The tube of the Taq Polymerase should not be thawed.

1.5 Additionally required equipment and materials

- DNA-Extraction kit
(e.g. SureFood® PREP Basic Art. No. S1052 / SureFood® PREP Advanced Art. No. S1053)
- real-time PCR instrument with two detection channels (510 nm and 580 nm)
- real-time PCR consumables (plates, tubes, foils, caps)
- pipettes with filter tips
- powder-free disposable gloves
- Vortex mixer
- micro centrifuge with a rotor for the reaction tubes

1.6 Precautions for users

- Extraction, PCR preparation and the PCR run should be separated in different rooms to avoid cross-contaminations.
- This test must only be performed by laboratory personnel trained in molecular biology methods.
- Strictly follow the working instructions.
- When handling samples, wear disposable gloves. After finishing the test, wash your hands.
- Do not smoke, eat or drink in areas where samples or test reagents are being used.
- Do not use the kit after the expiration date.
- All reagents and materials used have to be disposed properly after use. Please refer to the relevant national regulation for disposal.

1.7 Setup

	Blockcycler & R-Biopharm RIDA®CYCLER	Rotorcycler
Initial Denaturation (HOLD)	5 min, 95°C	1 min, 95°C
Cycles	45	45
Denaturation	15 sec, 95°C	10 sec, 95°C
Annealing/Extension (CYCLE)	30 sec, 60°C	15 sec, 60°C
Temperature Transition Rate/ Ramp Rate	Maximum	Maximum

1.8 Detection channel Set-up

Real-time PCR device	Detection	Detection channel	Quencher	Note
Agilent Mx3005P	pork	FAM	+	
	IAC	HEX	+	
Agilent AriaMx / Dx	pork	FAM	+	
	IAC	HEX	+	
Applied Biosystems 7500	pork	FAM	None	Check the passive reference option ROX is none.
	IAC	VIC	None	
Bio-Rad CFX96 / Dx / Opus	pork	FAM	+	Baseline Settings: • Baseline subtracted curve fit • Apply fluorescence drift correction
	IAC	HEX	+	
R-Biopharm RIDA®CYCLER	pork	green	+	Ignore cycles before , if there is a significant deviation in the baseline at the start of the run. Please see page 45 of the cyclor operating instructions, section 12.1.2 Cycling analysis parameter.
	IAC	yellow	+	
Qiagen Rotor-Gene Q	pork	green	+	Note: Please use only 0.1 ml reaction tubes. The gain settings must be set to 5 (factory default) for all channels.
	IAC	yellow	+	
Roche LightCycler® 480 II	pork	465-510	+	
	IAC	533-580	+	
Roche cobas® z 480 Analyzer	pork	465-510	+	
	IAC	540-580	+	

2 Qualitative Analysis

2.1 Protocol

2.1.1 Preparation of the master-mix

Calculate the total number of reactions needed (samples and control reactions) for the specific PCR assay.

Recommended control reactions for the specific PCR assay: negative control, extraction control and Positive Control.

The reaction mix contains an internal amplification control (inhibition control) per reaction.

Reactions needed for the qualitative detection of pork:

3 reactions for controls (1x negative control, 1x extraction control, 1 Positive Control)

For each sample: at least 1 reaction for each sample DNA

It is also recommended to prepare the master-mix with 10% additional volume in order to compensate reagent loss. Allow the reagents to thaw, mix and centrifuge before opening and use.

The tube of the Taq Polymerase should not be thawed and not be mixed by vortexing.

Example for the calculation and preparation of 10 reactions:

Components of the master-mix	Amount per reaction	10 reactions (with 10% excess)
Reaction Mix	19.9 µl	218.9 µl
Taq Polymerase	0.1 µl	1.1 µl
Total volume	20 µl	220 µl

Mix each master-mix well and centrifuge shortly before use.

2.1.2 Preparation of the real-time PCR-mix

- Pipette 20 µl of the master-mix into appropriate tubes/wells.
- Close the negative control.
- Pipette 5 µl of sample DNA into the designated tubes/wells and close them.
- Pipette 5 µl of Positive Control into the designated tubes/wells and close them.
- Centrifuge all tubes/plates shortly at low speed.
- Place tubes/plates into the real-time PCR instrument and start the run according to the setup.

*** Description of the controls**

- Negative control: only master-mix
- Extraction control: the extraction is performed without the sample – components from used Prep Kit
- Positive Control: master-mix and within the kit's provided Positive Control

2.2 Interpretation of results

The evaluation has to be made according to the usual analysis program recommended by the real-time PCR instrument manufacturer.

The control reactions have to show the correct results.

Pork DNA is detected in the FAM-channel. In the VIC/HEX-channel the amplification control is detected.

A sample is stated **positive** for pork, if the sample DNA shows amplification in the FAM-channel. High amplicon concentrations can result in a weak or absent signal of the internal amplification control (IAC).

A Cp value for the IAC is not needed to obtain a positive result of the positive control.

A sample is stated **negative** for pork, if the sample DNA shows no amplification in the FAM-channel and if the internal control (VIC/HEX-channel) of the sample is **positive** with a shift in Cp-Value ≤ 2 compared to the negative control.

If the sample DNA in the VIC/HEX-Channel shows **no amplification** or a shift in Cp-value > 2 compared to the negative control, it contains PCR inhibiting substances. A significant decrease in the fluorescence signal can also show the presence of PCR inhibiting substances. Under these circumstances DNA isolation and purification of the sample need to be improved. Alternatively the DNA can be diluted (recommendation 1:2 in PCR-water) and analysed again for inhibition. Please note that the dilution factor also affects the detection limit of the specific PCR assay.

3 Limitations of the method

- The presence of PCR inhibitors may cause invalid results.
- Extremely low levels of target below the limit of detection (LoD) may be detected, but results may not be reproducible.
- In highly processed samples, the limit of detection may be shifted. Factors such as high pressures, mechanical stresses, chemical treatment, extreme temperatures and/or extreme pH values during manufacturing process – such as in canning production – can damage or degrade nucleic acids. This means that the sensitivity of the test kit may be reduced and not all original components may be detected.

4 Further Information

4.1 Product Information

- Detailed information about setup of several real-time PCR devices
(Download: www.congen.de/en/downloads)
- Product-related documents (Download: www.congen.de/en/eifu/)
- Validation Report upon request

4.2 Technical Support

For further questions please contact your distributor or send an e-mail to sales@r-biopharm.de.

4.3 Distribution and Ordering

R-Biopharm AG
An der neuen Bergstrasse 17,
64297 Darmstadt, Germany
Phone: +49 (0) 61 51 - 81 02-0
Fax: +49 (0) 61 51 - 81 02-20
E-Mail: orders@r-biopharm.de
www.r-biopharm.com