CONGEN

SureFood® ALLERGEN 4plex Cereals

Art. No. S7006
100 rxn

User Manual

February 2026

 Inhalt

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Content

1	General Information	12
1.1	Description	12
1.2	Limit of Detection	12
1.3	DNA preparation	13
1.4	Kit components and storage	13
1.5	Additionally required equipment and materials	13
1.6	Precautions for users	13
1.7	Setup.....	14
1.8	Detection channel Set-up	14
2	Qualitative Analysis	16
2.1	Protocol	16
2.1.1	Preparation of the master-mix	16
2.1.2	Preparation of the real-time PCR-mix	16
2.2	Interpretation of results	17
3	Limitations of the method	18
4	Further Information	19
4.1	Product Information	19
4.2	Technical Support.....	19
4.3	Distribution and Ordering	19

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1 General Information

1.1 Description

The SureFood® ALLERGEN 4plex Cereals is a real-time PCR for the direct, qualitative detection and differentiation of specific wheat (*Triticum* spp.), barley (*Hordeum vulgare*) and rye (*Secale cereale*) DNA sequences according to directive (EC) 1169/2011 in food.

The test contains an internal amplification control (IAC). If the DNA contains PCR inhibiting substances, the signal of the amplification control will be affected or the amplification will be suppressed. Examples for PCR inhibiting substances are alcohols (e.g. ethanol, isopropanol), surfactants (e.g. CTAB, SDS, Triton X100) and salts (e.g. sodium chloride). In addition spices, herbs, algae, cocoa and further sample matrices might have PCR inhibiting effects.

The real-time PCR assay can be performed with commonly used real-time PCR instruments, equipped for detection of four fluorescence emissions at the channels FAM, VIC/HEX, ROX and Cy5 at the same time. The internal technical verification of instruments was performed on Agilent Mx3005P, Agilent AriaDx, Applied Biosystems 7500, Bio-Rad CFX96, Qiagen Rotor-Gene Q, R-Biopharm RIDA®CYCLER and Roche LightCycler® 480 II.

1.2 Limit of Detection

The SureFood® ALLERGEN 4plex Cereals real-time PCR has a limit of detection of ≤ 1 mg / kg using SureFood® PREP Advanced, protocol 1.

The limit of detection of the complete method (DNA extraction and real-time PCR) depends on sample matrix, processing grade, DNA preparation and DNA content.

The SureFood® PCR systems are very sensitive and therefore even a small amount of target DNA is sufficient for a successful analysis. The concentration of total DNA in the sample does not allow a conclusion on the quantity and quality of the target DNA.

Note: Inconsistent mixing ratios may cause a loss of sensitivity in the low concentration channel in mixed samples especially with high amplicon concentrations (Cp value < 20).

*e.g. 99.9 % wheat and 0.1 % rye.

1.3 DNA preparation

For DNA preparation the use of SureFood® PREP Basic (Art. No. S1052) or SureFood® PREP Advanced (Art. No. S1053) is recommended.

1.4 Kit components and storage

Kit Code	Reagent	Amount	Lid Color
1	Reaction Mix	2 x 1050 µl	Yellow
2	Taq Polymerase	1 x 80 µl	Dark Red
3	Positive Control	1 x 200 µl	Light Blue

Store all reagents at -28 to -16°C and protected from light. The Taq Polymerase can be stored at +2 to +8°C for multiple uses on the same day.

Note: The Taq Polymerase may be in a frozen or unfrozen state. This does not affect the quality of the Taq Polymerase or the performance of the real-time PCR.

1.5 Additionally required equipment and materials

- DNA extraction kit
(e.g. SureFood® PREP Basic Art. No. S1052 or SureFood® PREP Advanced Art. No. S1053)
- real-time PCR instrument with four detection channels (510 nm, 580 nm, 610 nm and 660 nm)
- real-time PCR consumables (plates, tubes, foils, caps)
- pipettes with filter tips
- powder-free disposable gloves
- Vortex mixer
- micro centrifuge with a rotor for the reaction tubes

1.6 Precautions for users

- Extraction, PCR preparation and the PCR run should be separated in different rooms to avoid cross-contaminations.
- This test must only be performed by laboratory personnel trained in molecular biology methods.
- Strictly follow the working instructions.
- When handling samples, wear disposable gloves. After finishing the test, wash your hands.
- Do not smoke, eat or drink in areas where samples or test reagents are being used.
- Do not use the kit after the expiration date.
- All reagents and materials used have to be disposed properly after use. Please refer to the relevant national regulation for disposal.

1.7 Setup

	Blockcycler & R-Biopharm RIDA®CYCLER	Rotorcycler
Initial Denaturation (HOLD)	5 min, 95°C	1 min, 95°C
Cycles	35	35
Denaturation	15 sec, 95°C	10 sec, 95°C
Annealing/Extension (CYCLE)	30 sec, 60°C	15 sec, 60°C
Temperature Transition Rate/ Ramp Rate	Maximum	Maximum

1.8 Detection channel Set-up

For additional details see cycler operating instructions for real-time PCR device.

Real-time PCR device	Detection	Detection channel	Quencher	Note
Agilent AriaDx/Mx	wheat	FAM	+	
	IAC	HEX	+	
	barley	ROX	+	
	rye	Cy5	+	
Agilent Mx3005P	wheat	FAM	+	
	IAC	HEX	+	
	barley	ROX	+	
	rye	Cy5	+	
Applied Biosystems 7500	wheat	FAM	+	Check the passive reference option ROX is none.
	IAC	HEX	+	
	barley	ROX	+	
	rye	Cy5	+	
Bio-Rad CFX96/Dx/Opus	wheat	FAM	+	Baseline Settings: • Baseline subtracted curve fit • Apply fluorescence drift correction
	IAC	VIC/HEX	+	
	barley	ROX	+	
	rye	Cy5	+	
R-Biopharm RIDA®CYCLER	wheat	green	+	Ignore cycles before , if there is a significant deviation in the baseline at the start of the run. Please see page 45 of the cycler operating instructions, section 12.1.2 Cycling analysis parameter.
	IAC	yellow	+	
	barley	orange	+	
	rye	red	+	

SureFood® ALLERGEN 4plex Cereals
(100 rxn)

Art. No. S7006

February 2026

Real-time PCR device	Detection	Detection channel	Quencher	Note
Qiagen Rotor-Gene Q	wheat	green	+	Note: Please use only 0.1 ml reaction tubes. The gain settings must be set to 5 (factory default) for all channels.
	IAC	yellow	+	
	barley	orange	+	
	rye	red	+	
Roche LightCycler® 480 II	wheat	465-510	+	The SureCC Color Compensation Kit I (Art. No. F4009) is required.
	IAC	533-580	+	
	barley	533-610	+	
	rye	618-660	+	

2 Qualitative Analysis

2.1 Protocol

2.1.1 Preparation of the master-mix

Calculate the total number of reactions needed (samples and control reactions) for the specific PCR assay.

Recommended control reactions for the specific PCR assay: negative control, extraction control and Positive Control.

The reaction mix contains an internal amplification control (inhibition control) per reaction.

Reactions needed for the qualitative wheat, barley and rye detection:

3 reactions for controls* (1x negative control, 1x extraction control, 1x Positive Control)

For each sample: at least 1 reaction for each sample DNA

It is also recommended to prepare the master-mix with 10 % additional volume in order to compensate reagent loss. Allow the reagents to thaw, mix and centrifuge before opening and use.

Example for the calculation and preparation of 10 reactions:

Components of the master-mix	Amount per reaction	10 reactions (with 10 % excess)
Reaction Mix	19.3 µl	212.3 µl
Taq Polymerase	0.7 µl	7.7 µl
Total volume	20 µl	220 µl

Mix each master-mix well and centrifuge shortly before use.

2.1.2 Preparation of the real-time PCR-mix

- Pipette 20 µl of the master-mix into appropriate tubes/wells.
- Close the negative control.
- Pipette 5 µl of sample DNA into the designated tubes/wells and close them.
- Pipette 5 µl of Positive Control into the designated tubes/wells and close them.
- Centrifuge all tubes/plates shortly at low speed.
- Place tubes/plates into the real-time PCR instrument and start the run according to the setup.

*** Description of the controls**

- Negative control: only master-mix
- Extraction control: the extraction is performed without the sample - components from used Prep Kit
- Positive Control: master-mix and within the kit's provided Positive Control

2.2 Interpretation of results

The evaluation has to be made according to the usual analysis program recommended by the real-time PCR instrument manufacturer.

The control reactions have to show the correct results.

Wheat DNA is detected in the FAM-channel, barley DNA is detected in the ROX-channel and rye DNA is detected in the Cy5-channel (see table). In the VIC/HEX-channel the internal amplification control (IAC) is detected.

A sample is stated **positive** for the respective parameter, if the sample DNA shows amplification in the respective channel. High amplicon concentrations can result in a weak or absent signal of the IAC.

A Cp value for the IAC is not needed to obtain a positive result of the Positive Control.

A sample is stated **negative** for the respective parameter, if the sample DNA shows no amplification in the respective channel and if the internal control (VIC/HEX-channel) of the sample is **positive** with a shift in Cp-value ≤ 2 compared to the negative control.

If the sample DNA in the VIC/HEX-Channel shows **no amplification** or a shift in Cp-value > 2 compared to the negative control, it contains PCR inhibiting substances. A significant decrease in the fluorescence signal can also show the presence of PCR inhibiting substances. Under these circumstances DNA isolation and purification of the sample need to be improved. Alternatively the DNA can be diluted (recommendation 1:2 in PCR-water) and analysed again for inhibition. Please note that the dilution factor also affects the detection limit of the specific PCR assay.

Result in the respective channel				Interpretation
FAM channel wheat	ROX channel barley	Cy5 channel rye	VIC/HEX channel IAC	
positive	negative	negative	positive/negative	wheat DNA detected
negative	positive	negative	positive/negative	barley DNA detected
negative	negative	positive	positive/negative	rye DNA detected
negative	negative	negative	positive	Negative, target DNA is not detected
negative	negative	negative	negative	invalid

Note: The results displayed in the table above represent merely an example. Additional combinations are also possible.

The following table shows the specification ranges of the kit controls

	Specification range	Internal amplification control (IAC)
negative control	negative	$25 \leq C_p \leq 34$
Positive Control (FAM – wheat)	$22 \leq C_p \leq 30$	not relevant
Positive Control (ROX – barley)	$22 \leq C_p \leq 30$	not relevant
Positive Control (Cy5 – rye)	$22 \leq C_p \leq 30$	not relevant

3 Limitations of the method

- The presence of PCR inhibitors may cause invalid results.
- Extremely low levels of target below the limit of detection (LoD) may be detected, but results may not be reproducible.
- Samples tested negative still could contain an allergen contamination below the limit of detection of the assay, or they might contain other allergen components like proteins or lipids for example.
- In highly processed samples, the detection limit may be shifted. Factors such as high pressures, mechanical stresses, chemical treatment, extreme temperatures and/or extreme pH values during manufacturing process – such as in canning production – can damage or degrade nucleic acids. This means that the sensitivity of the test kit may be reduced and not all original components may be detected.

4 Further Information

4.1 Product Information

- Detailed information about setup of several real-time PCR devices (Download: www.congen.de/en/downloads)
- Product-related documents (Download: www.congen.de/en/eifu/)
- Validation Report upon request

4.2 Technical Support

For further questions please contact your distributor or send an e-mail to sales@r-biopharm.de.

4.3 Distribution and Ordering

R-Biopharm AG
An der neuen Bergstrasse 17,
64297 Darmstadt, Germany
Phone: +49 (0) 61 51 - 81 02-0
Fax: +49 (0) 61 51 - 81 02-20
E-Mail: orders@r-biopharm.de
www.r-biopharm.com

r-biopharm®