## RIDASCREEN® Histamine (enzymatic)

**REF** **R1605**

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# RIDASCREEN® Histamine (enzymatic)

## Brief information

RIDASCREEN® Histamine (enzymatic) (Art. No. R1605) is an enzymatic test in microtiter plate format for the quantitative determination of histamine in fresh fish, canned fish, fish meal, wine (use RIDA® Sample Decolorant, Art. No. R1699, for sample preparation), and cheese.

All reagents required for the enzyme assay, including standards, are contained in the test kit. The test kit is sufficient for a maximum of 96 determinations (including standards). A microtiter plate spectrophotometer is required for quantification.

The RIDASCREEN® Histamine (enzymatic) test is:

- certified at AOAC-RI (PTM No. 031901)

Sample preparation:	Homogenization and extraction
Time requirement:	Sample preparation (10 fish samples) approx. 35 min Test implementation (incubation time) ..... 15 min
Standard:	Histamine
Limit of detection:	Fresh fish, canned fish ..... 0.75 mg/kg Fish meal ..... 3.75 mg/kg Wine..... 0.54 mg/kg Cheese.... 0.75 mg/kg
Limit of quantification:	Fresh fish, canned fish ..... 2 mg/kg Fresh fish incl. ascorbic acid ..... 3 mg/kg Fish meal ..... 10 mg/kg Wine..... 1.4 mg/kg Cheese... 2 mg/kg
Recovery:	Approx. 85 - 105 %
Specificity:	The enzyme is specific for histamine. There are no side reactions to L-histidine, L-tryptophan, serotonin, L-tyrosine, tyramine, L-lysine, trimethylamine, putrescine and cadaverine.

Further information is contained in the validation report.

The specificity of RIDASCREEN® Histamine (enzymatic) has been evaluated by the determination of the side activities of the respective substances in a buffer system (see validation report).

## **Related products for histamine determination**

RIDA® Sample Decolorant (Art. No. R1699) for wine analysis  
RIDASCREEN® Histamine (competitive ELISA, Art. No. R1601)

### **1. Intended use**

RIDASCREEN® Histamine (enzymatic) (Art. No. R1605) is used for the quantitative determination of histamine in fresh fish, canned fish, fish meal, wine and cheese.

### **2. General information**

Histamine is a product of the decomposition of histidine caused by the growth of certain bacteria in protein rich food like fish, cheese, and wine.

The amount of histamine formed depends on the bacterial species, the temperature, and the time of exposure and may exceed in fish 1000 mg/kg.

Fish of good quality contains less than 10 mg/kg histamine. The limit value for fish and products thereof is between 50 and 200 mg/kg depending on the fish species and the country. Similar high concentrations can be found in cheese, while histamine in wine usually does not exceed 15 mg/L.

### **3. Test principle**

The enzymatic determination is based on the histamine dehydrogenase which catalyzes the oxidative deamidation of histamine in the presence of an electron carrier that converts a dye to a color product. The color intensity is directly proportional to histamine concentration and is measured at 450 nm. The electron carrier and dye are coated on the microtiter plate.

## 4. Reagents provided

Each kit contains sufficient materials for a maximum of 96 measurements (including standard analyses). Each test kit contains:

Component	Cap color	Format	Histamine conc.	Volume
Microtiter plate	-	Ready to use		96 wells
Buffer	Transparent	Ready to use		15 mL
Standard 1	White	Ready to use	0 mg/L	1.3 mL
Standard 2	White	Ready to use	1 mg/L	1.3 mL
Standard 3	White	Ready to use	5 mg/L	1.3 mL
Standard 4	White	Ready to use	10 mg/L	1.3 mL
Standard 5	White	Ready to use	15 mg/L	1.3 mL
Standard 6	White	Ready to use	20 mg/L	1.3 mL
Enzyme solution	Red	Ready to use		1 mL
Spiking solution	Blue	Ready to use	500 mg/L	3 mL
Catalase Catalase to remove ascorbic acid	Black	Ready to use		1 mL

## 5. Reagents and equipment required but not provided

### 5.1. Equipment and reagents

- Distilled or deionized water
- Vortex
- Mixer, blender (e.g. Mr. Magic)
- Centrifuge, centrifuge vials
- Variable 100 - 200 µL micropipettes
- Multistep-pipette with 10 µL, 100 µL, 150 µL, 200 µL pipetting volume (e.g. Eppendorf combitips)
- Microtiter plate spectrophotometer (450 nm)
- Optional: RIDASOFT® Win.NET (Art. No. Z9996FF); use **version 1.104** or later for calculation

### Fresh fish analysis

- 50 mL polypropylene (PP) screw-cap vials (e.g. Greiner; Art. No. 227261)
- Homogenizer: laboratory mincer / grinder, Ultra-Turrax or mixer
- Water bath (100 °C / 212 °F)
- Paper filter (e.g. Whatman no.1 or similar quality)

### **Fresh fish analysis: As required for degradation of ascorbic acid**

- 1 N NaOH (e.g. Roth, Art. No. CN 58.1)
- 1 N HCl (e.g. Roth, Art. No. CN 63.1)
- H<sub>2</sub>O<sub>2</sub> solution, 30 % in water (e.g. Sigma, Art. No. 31642)
- Paper filter (e.g. Whatman no.1 or similar quality)

### **Wine analysis**

- 1.5 - 2 mL polypropylene vials (e.g. Eppendorf tube)
- RIDA® Sample Decolorant (Art. No. R1699) for sample preparation

### **Cheese analysis**

- 1.5 - 2 mL polypropylene vials (e.g. Eppendorf tube)
- 1 N perchloric acid (e.g. Bernd Kraft, Art. No. 20126.2700)
- 1 N KOH (e.g. Carl Roth, Art. No. K017.1)

## **6. Warnings and precautions for the users**

This test should only be carried out by trained laboratory personnel. The instruction for use must be strictly followed.

Use clean equipment and wear gloves before starting and during the assay in order to avoid cross-contamination.

Take precautions when using the boiling water bath to avoid being burnt or scalded.

This kit may contain hazardous substances. For hazard notes on the contained substances please refer to the appropriate material safety data sheets (SDS) for this product, available online at [www.r-biopharm.com](http://www.r-biopharm.com).

All reagents and materials must be recovered or disposed after use at customers own responsibility according to the protection of human health and the environment. Please observe the applicable national regulations concerning waste disposal (e.g. Waste Management Act, Regulations on Dangerous Chemicals, etc.).

## **7. Storage instructions**

Store the kit at 2 - 8 °C (35 - 46 °F). Do not freeze any test kit components.

Return any unused microwells to their original foil bag, reseal them together with the desiccant provided and further store at 2 - 8 °C (35 - 46 °F).

Do not use the test kit after the expiration date (see test kit label).

Do not interchange individual reagents between kits of different lot numbers.

## 8. Indication of instability or deterioration of reagents

- If the difference between absorption of  $A_2 - A_1 (= \Delta A)$  of standard 6 (e.g.  $\Delta A < 1.7$ ) is differing strongly from the certificate of analysis, then a deterioration of reagents has likely occurred
- The slope of the standard curve should lie in the range of 0.0847 - 0.1025 (mean 0.0936)

## 9. Preparation of Samples

Wear gloves before starting and during the assay to avoid cross-contamination.

The histamine concentration increases also in refrigerated food, therefore it is important to carry out the laboratory analysis directly after sampling.

Use only **plastic vials** (e.g. polypropylene, PP) for the sample preparation. Histamine adsorbs on glass surfaces which causes reduced recoveries.

### 9.1. Fresh fish and canned fish

- Weigh 5 g of homogenized sample into a 50 mL PP-screw cap vial.
- Pour 20 mL of dist. water over the sample.
- Close the vial and shake or vortex until the sample is **evenly suspended**.
- Incubate the sample for 20 min in a boiling water bath, take vials out of the water bath while wearing safety gloves after 10 min and shake immediately for 3 s so that samples are well suspended.
- Allow the sample to cool to room temperature in an ice bath for at least 2 min.
- Centrifuge at least 2,500 x g for 10 min.  
(Alternatively transfer 2 mL of the extract into centrifuge vials and centrifuge at high speed ( $> 10,000 \times g$ ) for 2 min in a micro-centrifuge. The extract should be as clear as possible. If particles are present, then filtrate the extract.)
- If a layer of fat has formed on top of the extract after centrifugation then hold the vial angular and take the extract from the side. Alternatively, carefully withdraw with a pipette the lower layer and deliver to a new vial, filter / centrifuge again.
- Use 100  $\mu\text{L}$  of the undiluted **clear extract** per well.
- The extracted sample is stable at room temperature (20 - 25 °C / 68 - 77 °F) for 2 h, or at 2 - 8 °C (36 - 46 °F) for 7 days.

### 9.1.1. Degradation of ascorbic acid (> 250 mg/kg) in fish

- If fresh fish has been treated with ascorbic acid (analysis using test kits Art. No. E1267 or 10409677035), an intense yellow color may occur (due to a creep reaction) during the subsequent enzymatic analysis within seconds ( $A_1$  measurement e.g. 3.0).

The ascorbic acid is degraded by  $H_2O_2$  / catalase as follows:

- Transfer 1 mL of the clear supernatant (see chapter 9.1) in a 50 mL PP screw cap vial.
- Add 100  $\mu$ L 1 N NaOH.
- Add 100  $\mu$ L 30 %  $H_2O_2$ .
- Mix softly and incubate 10 min at room temperature (20 - 25 °C / 68 - 77 °F).
- Add 100  $\mu$ L 1 N HCl.
- Mix gently.
- Add 10  $\mu$ L Catalase.
- Mix softly and incubate 10 min at room temperature (20 - 25 °C / 68 - 77 °F).
- Put the mixture for 5 min into a boiling water bath.
- Let cool down.
- Use 100  $\mu$ L of the **clear extract** per well.
- The extracted sample is stable at room temperature (20 - 25 °C / 68 - 77 °F) for 2 h, or at 2 - 8 °C (36 - 46 °F) for 7 days.

For validation spike the sample and determine the recovery.

### 9.2. Fish meal

- Weigh 1 g of homogenized fish meal into a 50 mL PP-screw cap vial.
- Pour 25 mL of dist. water over the sample.
- Close the vial and shake or vortex until the sample is **evenly suspended**.
- Incubate the sample for 20 min in a boiling water bath, take vials out of the water bath while wearing safety gloves after 10 min and shake immediately for 3 s so that samples are well suspended.
- Allow the sample to cool to room temperature in an ice bath for at least 2 min.
- Centrifuge at least 2,500 x g for 10 min.  
(Alternatively transfer 2 mL of the extract into centrifugal vials and centrifuge at high speed (> 10,000 x g) for 2 min in a micro-centrifuge. The extract should be as clear as possible. If particles are present, then filtrate the extract.

- If a layer of fat has formed on top of the extract after centrifugation then hold the vial angular and take the extract from the side. Alternatively, carefully withdraw with a pipette the lower layer and deliver to a new vial, filter / centrifuge again.
- Use 100  $\mu\text{L}$  of the undiluted **clear extract** per well.
- The sample extracts can be stored at room temperature (20 - 25  $^{\circ}\text{C}$  / 68 - 77  $^{\circ}\text{F}$ ) for 2 h, at 2 - 8  $^{\circ}\text{C}$  (36 - 46  $^{\circ}\text{F}$ ) for 3 days or at -20  $^{\circ}\text{C}$  (- 4  $^{\circ}\text{F}$ ) for 3 months.

### 9.3. Wine

Wine (red wine, white wine, rosé) contains anthocyanins and polyphenols which have to be removed prior to the enzymatic analysis.

- Add 200  $\mu\text{L}$  Reagent 1 of RIDA<sup>®</sup> Sample Decolorant to 200  $\mu\text{L}$  wine (rinse pipette tip).
- Add 200  $\mu\text{L}$  Reagent 2 of RIDA<sup>®</sup> Sample Decolorant, vortex, and incubate 5 min at room temperature (20 - 25  $^{\circ}\text{C}$  / 68 - 77  $^{\circ}\text{F}$ ).
- Centrifuge 2 min at a minimum of 14,000 x g at room temperature (20 - 25  $^{\circ}\text{C}$  / 68 - 77  $^{\circ}\text{F}$ ).
- Transfer 500  $\mu\text{L}$  of the supernatant in a new vial.
- Add 100  $\mu\text{L}$  Reagent 3 of RIDA<sup>®</sup> Sample Decolorant, vortex and incubate 5 min at room temperature (20 - 25  $^{\circ}\text{C}$  / 68 - 77  $^{\circ}\text{F}$ ).
- Centrifuge at 14,000 x g at room temperature (20 - 25  $^{\circ}\text{C}$  / 68 - 77  $^{\circ}\text{F}$ ) for 2 min.
- Use 100  $\mu\text{L}$  of the **clear extract** per well.
- The sample extract can be stored at room temperature (20 - 25  $^{\circ}\text{C}$  / 68 - 77  $^{\circ}\text{F}$ ) for 2 h or at 2 - 8  $^{\circ}\text{C}$  (36 - 46  $^{\circ}\text{F}$ ) for 1 day.

### 9.4. Cheese

If the cheese rind is edible then it should also be part of the representative sample.

- Weigh 60 g of cheese into the plastic cup for blending.
- Add 140 mL of water and mix for 15 s to suspend all particles. Repeat mixing two more times to suspend all particles.
- Remove the fat layer with a spatula.
- Transfer 50 mL of the unclear solution to a plastic vial.
- Centrifuge at 3,500 x g for 10 min (at 4  $^{\circ}\text{C}$  if possible; otherwise cool sample down in an ice bath).
- Remove the fat layer with a spatula.

- Transfer 0.8 mL of the clear solution to a plastic vial.
- Add 200 µL 1 N perchloric acid.
- Mix well.
- Centrifuge at 20,000 x g for 2 min at room temperature (20 - 25 °C / 68 - 77 °F).
- Transfer 500 µL of the clear supernatant to a new centrifugal plastic vial.
- Add 100 µL 1 N KOH and mix.
- Centrifuge 2 min at of least 20,000 x g at room temperature (20 - 25 °C / 68 - 77 °F).
- Use 100 µL of **clear extract** per well.
- The sample extract can be stored at room temperature (20 - 25 °C / 68 - 77 °F) for 2 h or at 2 - 8 °C (36 - 46 °F) for 18 h.

## 10. Test implementation

### 10.1. Test preparation

Bring all reagents to room temperature (20 - 25 °C / 68 - 77 °F) and mix carefully before use.

### 10.2. Test implementation

Insert a sufficient number of wells into the microwell holder for all standard and samples to be run in duplicate. Record standard and sample positions.

1. Using a multi-stepper, add 150 µL of buffer to the wells and carefully shake the plate manually for 3 sec.
2. Add 100 µL of standards, controls or samples (prepared according to chapter 9) to separate wells in duplicate. Flush the pipette tips. Thereafter, carefully shake the plate manually for 3 sec.
3. After 3 min\* (for wine samples after 10 min), the absorption  $A_1$  is measured at 450 nm **without** a reference filter (e.g. 620 nm) (avoid any bubbles in the wells).
4. Add 10 µL of the blue dyed enzyme solution with a multi-stepper to each well. Thereafter, carefully shake the plate manually for 3 sec.
5. After 10 min, the absorption  $A_2$  is measured at 450 nm **without** a reference filter (e.g. 620 nm). For evaluation use the RIDASOFT® Win.NET, select  $A_1$  and  $A_2$  and calculate.

\* For **fish** an **increase of yellow color** may occur (creep reaction).

- If  $A_1$  is e.g. 0.3 then the incubation time is increased to 10 min.

- If  $A_1$  is e.g. 3.0 then the fish sample should be prepared with catalase (see 9.1.1).

## 11. Evaluation

At R-Biopharm AG a special software, the **RIDASOFT® Win.NET (Art. No. Z9996FF)**, is available for evaluation of the RIDASCREEN® assays. The calculation is be done by using a linear regression, the **dilution factor** of the samples has to be **taken into account**. (From software version 1.105 onwards, the default factor is 5 which needs to be modified if required).

Sample	Dilution factor	Measuring range
Fresh / canned fish	5	2 - 100 mg/kg
Fish incl. degradation of ascorbic acid	6.55	2.62 - 131 mg/kg
Fish meal	25	10 - 500 mg/kg
Wine	3.6	1.4 - 72 mg/kg
Cheese	5	2 - 100 mg/kg

For the evaluation it should be clarified, that the quality criteria are fulfilled for the current test run. The course of the standard curve can be taken from the attached analysis certificate.

The slope of the standard curve should lie in the range of 0.0847 - 0.1025 (mean 0.0936). The residuals for  $\Delta A$  should lie between - 0.05 und 0.05 (refer to the residual plot, last page of the RIDASOFT® Win.NET report). Larger deviation of single wells are indicating (almost) always an insufficient precision of pipetting, therefore the pipetting technique needs to be improved.

The coefficient of variation (CV) should not exceed 10 % for samples and standards (except for CVs calculated from standard 1 and samples close to standard 1). The recovery of check samples is usually between 75 % and 125 %.

Higher values of the absorbance ( $A_{450\text{ nm}}$ ) for the calibration curve, especially for the zero standard, may be a result of histamine contamination. If absorbance values ( $A_{450\text{ nm}}$ ) > standard 6, dilute the final test extract with water and perform another assay.

## 12. Recommendation

In order to ensure a high analytical performance we recommend:

- Each final test extract should be assayed in duplicate.
- Ensure accurate pipetting.
- Carry along test controls for quality control and to ensure an accurate and correct test procedure. Histamine-free and histamine-containing (naturally contaminated or spiked) samples (e.g. FAPAS® reference samples) should be used.
- For spiking, the spike solution contained in the kit should be used.

### 10 mg/L spike control:

- 1:50 dilution of the spiking solution Histamine (500 mg/L) e.g. 100 µL + 4.9 mL water.
- Use 100 µL in the assay.

### 50 mg/kg histamine spiked fish sample:

- Pipette 500 µL of the spiking solution Histamine (500 mg/L) to 5 g homogenized fish.
- Extract the sample with 19.5 mL of water (see chapter 9.2.).

### 250 mg/kg histamine spike fish meal sample:

- Pipette 500 µL of the spiking solution Histamine (500 mg/L) to 1 g homogenized fish meal sample.
- Extract the sample with 24.5 mL of water (see chapter 9.3.).

### 10 mg/L histamine spiked wine sample:

- Pipette 200 µL of the spiking solution Histamine (500 mg/L) to 9.8 mL wine.
- Extract the sample with RIDA® Sample Decolorant (see chapter 9.4.).

## 13. Further application notes

- Wine – More sensitive analysis with RIDASCREEN® Histamine (enzymatic).
- Reduced sample preparation time for fresh fish and canned fish.

For additional applications (e.g. milk, yeast extract, pet food) and product information (e.g. automation, training videos) please contact your local distributor or R-Biopharm AG at this address: [sales@r-biopharm.de](mailto:sales@r-biopharm.de).

## Version overview

Version number	Chapter and title
2019-02-26	Release version
2022-06-13	Current version Changes made: <ul style="list-style-type: none"><li>– General linguistic revision</li><li>– Addition of AOAC-RI certificate (certified since 2019)</li></ul>

## Explanation of symbols

General symbols:

Follow the instructions for use

Batch number

Expiry date (YYYY-MM)

Storage temperature

Article number

Number of test determinations

Manufacturing date (YYYY-MM)

Manufacturer + address

## Disclaimer

The user assumes all risk in using R-Biopharm AG's products and services.

R-Biopharm AG will warrant that its products and services meet all quality control standards set by R-Biopharm AG, and R-Biopharm AG will, at its option, replace or repair any components, product or repeat services which prove to be defective in workmanship or material within product specific warranty periods or expiration dates and which our examination shall disclose to our satisfaction to be defective as such.

This warranty is expressly in lieu of all other warranties, expressed or implied, as to quality, description, fitness for any particular purpose, merchantability, productiveness, or any other matter. R-Biopharm AG shall be in no way responsible for the proper use of its products and hereby disclaims all other remedies, warranties, guarantees or liabilities, expressed or implied, arising by law or otherwise, and it shall have no liability for any lost profits or damage, direct, indirect or otherwise, to person or property, in connection with the use of any of its products or services.

This warranty shall not be extended, altered or varied except by a written instrument signed by an authorized representative of R-Biopharm AG.

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are registered trademarks of R-Biopharm AG.