Chicken Pancreatin (γ - Glutamylhydrolase)

Art. No.: P2002

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Chicken Pancreatin (γ - Glutamylhydrolase)

Brief information

Chicken pancreatin (γ - glutamylhydrolase) for the enzymatic treatment of food and animal feed to determine the natural folic acid content. Chicken pancreatin is used for the enzymatic treatment of food prior to the determination of its natural folic acid content. The natural folic acid is hydrolysed with chicken pancreatin into foylmonoglutamate or foyldiglutamate. Chicken pancreatin is ideal for microbiological folic acid determination in microtiter plate format (VitaFast[®]). It is also suitable for the determination of the natural folic acid content using other methods. One bottle contains 500 mg chicken pancreatin and is sufficient for 50 sample treatments.

1. Reagents provided

Chicken pancreatin (γ - glutamylhydrolase) 500 mg (solid), shelf life 1 year at - 20 to - 30 °C (- 4 to - 22 °F).

2. Warning and precautions for the user

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3. Preparation

- open the chicken pancreatin bottle, place the rubber stopper with the open side up
- add 5 ml redist. or deionized water
- close the bottle with the rubber stopper and dissolve by shaking
- prepare 100 μ l aliquots from the chicken pancreatin solution in 0.5 ml (or 1.0 ml) reaction vials and close the reaction vials (sufficient for 50 sample treatments)
- store the reaction vials at - 20 to - 30 °C (- 4 to - 22 °F) (shelf life 6 months)
- one 100 μ l aliquot is sufficient for 1 g sample extraction with phosphate buffer (0.05 mol / l; 0.1 % ascorbate; pH 7.2)