## SureFood® VIRUS Hepatitis A PLUS (100 Reakt.)

Art.-Nr. S7125

Version 1.0

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3. Preparation of the PCR-mix
  - Pipette 20 µl of the master-mix into appropriate tubes or wells.
  - For the negative control pipette 5 µl of PCR Water into the designated tubes and close them.
  - Pipette 5 µl of sample RNA into the designated tubes and close them.
  - Pipette 5 µl of the Positive Control into the designated tubes and close them.
  - Centrifuge all tubes/plates shortly at low speed.
  - Place tubes/plates into the real-time PCR instrument and start the run according to the setup.

#### Interpretation of results

The evaluation has to be made according to the usual analysis program recommended by the real-time PCR instrument manufacturer. The control reactions need to give the correct results.

A sample is stated **positive** for hepatitis A, if the sample RNA shows an amplification in the FAM-channel. A sample is stated **negative** for hepatitis A, if the sample RNA shows no amplification in the FAM-channel and the internal amplification control (VIC/HEX-channel) of the sample is **positive**.

If the sample RNA and the internal amplification control are **negative** the sample contains PCR-inhibiting substances or the sample preparation was not successful. Under these circumstances an evaluation of the samples is not possible. RNA isolation and purification for the sample need to be improved.

#### Product Information

- Validation Report
- Material Safety Data Sheet
- Product Information

#### Technical Support

For further questions please contact your distributor or send an e-mail to sales@r-biopharm.de.

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**Description**

The test is a real-time RT-PCR for the direct and qualitative detection of hepatitis A virus. The test contains an Internal Control RNA (ICR) as an internal control of sample preparation procedure and to determine possible PCR-inhibition. The real-time PCR assay can be used with established real-time PCR instruments, equipped for detection of two fluorescence emissions at 522 nm and 553 nm (FAM and VIC) at the same time (Rotor-Gene Q, Stratagene MxSeries, Roche LightCycler® 480/2.0\*/1.5\* etc.).

**Limit of Detection**

The SureFood® VIRUS Hepatitis A PLUS real-time RT-PCR has a limit of detection of ≤ 50 RNA copies. The assay limit of detection depends on RNA preparation and RNA content.

**RNA-preparation**

For RNA-preparation the use of SureFood® PREP DNA/RNA Virus is recommended.

The test assay contains an Internal Control RNA (ICR), which can either be used as PCR inhibition control or as extraction control for the sample preparation procedure and as a PCR inhibition control.

If the ICR is used only as a PCR inhibition control, 1 µl of the ICR should be added to the Master-Mix.

If the ICR is used as an extraction control for the sample preparation procedure and as PCR inhibition control, 20 µl of the ICR should be added during extraction procedure. The ICR should always be added to the specimen-lysis buffer mixture and must not be added directly to the specimen.

**Kit components and storage**

2x	Reaction Mix (700 µl)	<b>(Code 1)</b>
1x	PP-Mix (770 µl)	<b>(Code 2)</b>
1x	Enzyme Mix (80 µl)	<b>(Code 3)</b>
2x	Internal Control RNA (1800 µl)	<b>(Code R)</b>
1x	PCR Water (500 µl)	<b>(Code N)</b>
1x	Positive Control (100 µl)	<b>(Code P)</b>

Store all reagents at -20°C and protected from light.

**Additionally required equipment and materials**

- Real-time PCR instrument, equipped with two detection channels at 522 nm and 553 nm
- Real-time PCR consumables (plates, tubes, foils, caps)
- Pipettes with filter tips

**Protocol**

1. Preparation of the master-mix

Calculate the total number of reactions needed (samples and control reactions). Recommended control reactions: negative control, positive control and extraction control. The test assay contains an Internal Control RNA (ICR), which can either be used as PCR inhibition control or as extraction control.

\*note: For users of the Roche LightCycler® 2.0 and 1.5 a Color Compensation is necessary. The use of the SureCC Color Compensation Kit II (Art.-No. S4010) is recommended for the color compensation of such devices.

It is also recommended to prepare the master-mix with 10 % additional volume in order to compensate reagent loss. Allow the reagents to thaw, mix by vortexing and centrifuge before opening and use.

Example for the calculation and preparation of 10 reactions for ICR as extraction and PCR inhibition control:

Components for master-mix	Amount per reaction	10 reactions (with 10% excess)
Reaction Mix	12.5 µl	137.5 µl
PP-Mix	6.9 µl	75.9 µl
Enzyme Mix	0.7 µl	7.7 µl
<b>Total volume</b>	<b>20.1 µl</b>	<b>221.1 µl</b>

Mix each master-mix well and centrifuge shortly before use.

Example for the calculation and preparation of 10 reactions for ICR only as PCR inhibition control:

Components for master-mix	Amount per reaction	10 reactions (with 10% excess)
Reaction Mix	12.5 µl	137.5 µl
PP-Mix	6.9 µl	75.9 µl
Internal Control RNA	1.0 µl	11.0 µl
Enzyme Mix	0.7 µl	7.7 µl
<b>Total volume</b>	<b>21.1 µl</b>	<b>232.1 µl</b>

Mix each master-mix well and centrifuge shortly before use.

2. Setup

	<b>Blockcycler/ Rotorcycler</b>
RT-Reaktion	10 min, 58°C
Denaturierung	1 min, 95°C
Cycles	45
Denaturation	15 sec, 95°C
Annealing/Extension	30 sec, 55°C
Temperature Transition Rate/ Ramp Rate	Maximum
Fluorescence Detection Setup	Detection: End of extension phase Detection System: Various devices <b>FAM</b> -channel, Quencher: BHQ Rotor-Gene Q Green LightCycler 480 II 465 nm - 510 nm  Internal Amplification Control: Various devices <b>VIC/HEX</b> -channel, Quencher: BHQ Rotor-Gene Q Yellow LightCycler 480 II 533 nm - 580 nm
Detailed information on the setup of several real-time PCR devices is available at the Congen homepage: <a href="http://www.congen.de/en/information/device-settings-for-surefood-real-time-pcr-kits.html">http://www.congen.de/en/information/device-settings-for-surefood-real-time-pcr-kits.html</a>	