English

Deutsch

Français

Compact Dry ETC is a ready to use, selective and chromogene plate for the detection and enumeration of Enterococcus

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Specimen pretreatment

Viable count in water or liquid foodstuff

Drop 1 ml of specimen (dilute if necessary) on the middle of the Compact Dry plate.

Viable count in solid foodstuff

Add buffer solution to the specimen and homogenize by stomacher®. Drop 1 ml of specimen (dilute if necessary) on the middle of the dry sheet of the Compact Dry plate.

Viable count in swab test specimen

Use a steril wet swab to wipe the surface which should be analyzed and put it back into the device with wiping solution. Drop 1 ml of wiping solution (dilute if necessary) on the middle of the Compact Dry plate. Use the Compact Dry Swab (Art. No. ZCS1002953, content 40 pcs) as single application (1 ml) or Promedia ST-25 (Art. No. Z0302, content 10 pcs) as multiple application (10 ml) for this procedure.

Test instructions

1. Open the cap and drop 1 ml of specimen on the middle of the Compact Dry plate.
2. Specimen diffuses automatically and evenly into the sheet and transforms the dried sheet into a gel within seconds.
3. Put the cap again on the plate and write the information needed on the memorandum section.
4. Turn over the capped plate and put in the incubator.
5. After incubation count the number of blue/blue green colored colonies underneath the plate. White paper placed under the plate helps to count the colonies.

Incubation time 20 - 24 hours
Incubation temperature 37 ± 1 °C

Please use the incubation time / temperature according to the national food analysis recommended for *Enterococcus*

Interpretation of the results

Colonies grown are almost all blue / blue-green. Bacteria other than *Enterococcus* are inhibited to grow and they do not form any colonies.

Storage and shelf life

Keep at room temperature (+ 1 to + 30 °C). Total shelf life 18 months after manufacturing.

Notes

- Some colonies might not be clearly blue/blue-green colored.
- High concentrations on plates (> 300 cfu) will cause the entire growth area to become blue / blue-green. In this case dilute the specimen.
- After use please follow the current disposal regulations.
- The growth area is 20 cm². The back of the plate has a grid carved of 1 cm x 1 cm to make the colony counting easier. In case of any difficulties to count colonies due to large number of colonies grown, total viable count can be obtained by multiplying 20 by an average number of colonies per grid counted from several grids.
- Compact Dry plates are produced at an ISO 9001 certified site.

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