SureFood® GMO ID MON863 Corn (100 React.)

Art.-No. S2037

Version 1.0

2. Setup

	Blockcycler	Rotorcycler
Initial Denaturation (HOLD) Cycles	5 min, 95°C	1 min, 95°C
Denaturation	45	45
Annealing/Extension (CYCLE)	15 sec, 95°C	10 sec, 95°C
	30 sec, 60°C	15 sec, 60°C
Temperature Transition Rate/Ramp Rate	Maximum	Maximum
Fluorescence Detection Setup	Detection: End of extension phase Reporter Dye: FAM Quencher Dye: BHQ Passive Reference: none	LightCycler® Channel: 530 or F1 Acquisition mode: Single in extension phase Rotor-Gene Q Reporter Dye: FAM (Green)

Detailed information on the setup of several real-time PCR devices is available at the Congen homepage: <http://www.congen.de/en/information/device-settings-for-surefood-real-time-pcr-kits.html>

3. Preparation of the PCR-mix

- Pipette 20 µl of the master-mix into appropriate tubes/wells or capillaries.
- Close the negative control (the negative control is ready for PCR without any addition).
- Pipette 5 µl of sample DNA into the designated tubes/wells or capillaries and close them.
- Pipette 5 µl of the Positive Control into the designated tubes/wells or capillaries and close them.
- Centrifuge all tubes/wells or capillaries shortly at low speed.
- Place tubes/wells or capillaries into the PCR instrument and start the run according to the setup.

Interpretation of results

The evaluation has to be made according to the usual analysis program recommended by the real-time PCR instrument manufacturer. The control reactions need to give the correct results.

A sample is stated **positive**, if the sample DNA shows an amplification signal in the MON863 Corn system.

A sample is stated **negative**, if the sample DNA shows no amplification in the MON863 Corn system and a performed amplification control (inhibition control) with S2026 or S2126 of the sample is **positive**. If the sample DNA and a performed amplification control are **negative** the sample contains PCR-inhibiting substances. Under these circumstances an evaluation of the samples is not possible. DNA isolation and purification for the sample need to be improved.

Product Information

- Validation Report
- Product Information
- Material Safety Data Sheet

Technical Support

For further questions please contact your distributor or send an e-mail to sales@r-biopharm.de.

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2x	Reaction Mix (1,1 ml)	(Code 1)
1x	Taq Polymerase (11 µl)	(Code 2)
1x	Positive Control (200 µl)	(Code 3)

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	Blockcycler	Rotorcycler
Initial Denaturation (HOLD) Cycles	5 min, 95°C 45	1 min, 95°C 45
Denaturation Annealing/Extension (CYCLE)	15 sec, 95°C 30 sec, 60°C	10 sec, 95°C 15 sec, 60°C
Temperature Transition Rate/ Ramp Rate	Maximum	Maximum
Fluorescence Detection Setup	Detection: End of extension phase Reporter Dye: FAM Quencher Dye: BHQ Passive Reference: none	LightCycler® Channel: 530 oder F1 Acquisition mode: Single in extension phase Rotor-Gene Q Reporter Dye: FAM (Green)

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Description

The test detects MON863 Corn DNA. Therefore the kit contains an event specific real-time PCR system for the MON863 Corn (OECD unique identifier MON-ØØ863-5). The real-time PCR assay can be used with established real-time PCR instruments (Roche LightCycler®, Rotor-Gene Q, ABI PRISM, Eppendorf realplex, BioRad CFX96, Stratagene MxSeries etc.).

Limit of Detection

The SureFood® GMO ID MON863 Corn assay has a limit of detection of ≤ 5 DNA copies. The assay limit of detection depends on sample matrix, processing grade, DNA preparation and DNA content.

DNA-preparation

For DNA-preparation of raw material the use of SureFood® PREP Plant and for highly processed food and feed the use of SureFood® PREP Plant X is recommended.

Kit components and storage

2x	Reaction Mix (1.1 ml)	(Code 1)
1x	Taq Polymerase (11 µl)	(Code 2)
1x	Positive Control (200 µl)	(Code 3)

Store all reagents at –20°C and protected from light.

Additionally required equipment and materials

- Real-time PCR instrument
- Real-time PCR consumables (plates, tubes, foils, capillaries)
- Pipettes with filter tips

Protocol

1. Preparation of the master-mix

Calculate the total number of reactions needed (samples and control reactions). Recommended control reactions: negative control, positive control, extraction control and an inhibition control. For the preparation of the inhibition control the use of the SureFood® GMO 4plex 35S/NOS/FMV+IAC kit (Art.-No. S2126) or the SureFood® GMO 35S/NOS/FMV kit (Art.-No. S2026), respectively, is recommended.

It is also recommended to prepare the master-mix with 10 % additional volume in order to compensate reagent loss. Allow the reagents to thaw, mix by vortexing and centrifuge before opening and use.

Example for the calculation and preparation of 10 reactions:

Components for master-mix	Amount per reaction	10 reactions (with 10% excess)
Reaction Mix	19.9 µl	218.9 µl
Taq Polymerase	0.1 µl	1.1 µl
Total volume	20 µl	220 µl

Mix each master-mix well and centrifuge shortly before use.

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R-Biopharm AG