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Set of 3 processed Soya Assay Controls

1. Intended Use

The processed assay controls are used to control the extraction, the test implementation and the handling of the RIDASCREEN[®]FAST Soya (R7102).

2. Content

Each assay control contains 2 g homogenized cookie.

The Set of 3 processed assay controls is composed of homogenized cookies with a defined content of soya. Assay control A consists of a soya-free cookie. To produce the assay controls B and C different concentrations of soya-free and a soya-containing cookie were mixed.

3. Storage

Store sample at 2 - 8 °C (35.6 - 46.4 °F) in the dark.

4. Preparation of samples and test procedure

To avoid cross contamination, it is recommended to extract the negative control sample first. Extract the samples and carry out the ELISA procedure as explained in the product information of RIDASCREEN[®]FAST Soya (R7102).

Remark:

The supernatant obtained after centrifugation should be transferred into a separate vial. The supernatant can be stored up to eight weeks in a tightly closed vial in the dark at - 20 °C / - 4 °F.

5. Results

The ELISA RIDASCREEN[®]FAST Soya quantifies soy proteins in raw products as well as in **processed** food. If the results do not correlate with the values in the table, a trouble shooting of the ELISA procedure has to be carried out.

Assay control	Dilution factor (final)	MV (ppm soya protein)	CV (%)	Range (ppm soya protein)
A (n=10; r=30)	100	< LOQ		
B (n=10; r=30)	100	4.6	9.4	3.2 - 5.9
C (n=10; r=30)	100	13.5	6.3	9.3 - 17.3

n = number of extractions; r = number of repeated measurements
 MV : mean value
 CV : coefficient of variation
 LOQ: Limit of Quantification

Remark

The soya assay controls are **not** a reference material. For validation of the RIDASCREEN®FAST Soya Kit spiked samples should be used.

The data corresponds to our present state of technology and provides information on our products and their uses. R-Biopharm makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. Defective products will be replaced. There is no warranty of merchantability of this product, or of the fitness of the product for any purpose. R-Biopharm shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.

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are registered trademarks of R-Biopharm AG