CONGEN

SureFast[®]
Histomonas meleagridis ONE

Art. No. F5213
100 rxn

User Manual

July 2019

 **Inhalt**

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Content

1	General Information	7
1.1	Description	7
1.2	Limit of Detection	7
1.3	Kit components and storage	7
1.4	Additionally required equipment and materials	7
1.5	Setup	8
1.6	Detection channel Set-up	8
2	Qualitative Analysis	9
2.1	Protocol	9
2.1.1	DNA-preparation	9
2.1.2	Preparation of the master-mix	9
2.1.3	Preparation of the real-time PCR-mix	10
2.2	Interpretation of results	10
3	Further Information	10
3.1	Product Information	10
3.2	Technical Support	10

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1 General Information

1.1 Description

The SureFast® *Histomonas meleagridis* ONE can be applied for the fast and simple isolation and detection of *Histomonas meleagridis* DNA in tissue samples.

This method comprises the following three steps: DNA-preparation, specific real-time PCR detection and interpretation of results.

Each reaction contains an internal amplification control (IAC). If the DNA contains PCR inhibiting substances, the signal of the amplification control will be affected or the amplification will be suppressed. Examples for PCR inhibiting substances are alcohols (e.g. ethanol, isopropanol), surfactants (e.g. CTAB, SDS, Triton X100) and salts (e.g. sodium chloride).

The real-time PCR assay can be performed with commonly used real-time PCR instruments, equipped for detection of two fluorescence emissions at the channels FAM and VIC/HEX at the same time. The technical validation of instruments was performed on Agilent Mx3005P, Bio-Rad CFX96, Roche LightCycler® 480 II, Roche LightCycler® 2.0, Applied Biosystems 7500, Qiagen Rotor-Gene Q and R-Biopharm RIDA®CYCLER.

1.2 Limit of Detection

The SureFast® *Histomonas meleagridis* ONE real-time PCR has a limit of detection of ≤ 5 DNA copies. The assay limit of detection depends on sample matrix, processing grade, DNA preparation and DNA content.

1.3 Kit components and storage

Kit Code	Reagent	Amount	Lid Color
1	Reaction Mix	2 x 1100 μ l	Yellow
2	Taq Polymerase	1 x 80 μ l	Dark Red
3	Positive Control	1 x 200 μ l	Light Blue
L	Lysis Buffer	2 x 25 ml	Clear

Store all reagents at -20°C and protected from light. The Taq Polymerase can be stored at $+2$ to $+8^{\circ}\text{C}$ for multiple uses on the same day.

The Lysis Buffer should be stored at -20 to $+8^{\circ}\text{C}$. Reagents should be used until the expiry date indicated on the label.

Note: The Taq Polymerase may be in a frozen or unfrozen state. This does not affect the quality of the Taq Polymerase or the performance of the real-time PCR.

1.4 Additionally required equipment and materials

- real-time PCR instrument with two detection channels (510 nm and 580 nm)
- real-time PCR consumable (plates, tubes, capillaries, foils, caps)
- pipettes with filter tips
- 2 ml tubes
- powder-free disposable gloves
- Vortex mixer
- centrifuge with a rotor for the reaction tubes
- Thermomixer/heating block (up to 95°C)

1.5 Setup

	Blockcycler & R-Biopharm RIDA®CYCLER	Rotorcycler & LightCycler® 480 II & LTF MyGo Pro
Initial Denaturation (HOLD)	1 min, 95°C	1 min, 95°C
Cycles	45	45
Denaturation	15 sec, 95°C	10 sec, 95°C
Annealing/Extension (CYCLE)	30 sec, 60°C	15 sec, 60°C
Temperature Transition Rate/ Ramp Rate	Maximum	Maximum

1.6 Detection channel Set-up

Real-time PCR device	Detection	Detection channel	Quencher	Note
Agilent Mx3005P	<i>Histomonas meleagridis</i>	FAM	+	
	IAC	HEX	+	
Agilent AriaDx	<i>Histomonas meleagridis</i>	FAM	+	
	IAC	HEX	+	
Applied Biosystems 7500	<i>Histomonas meleagridis</i>	FAM	None	Check the passive reference option ROX is none.
	IAC	VIC	None	
Bio-Rad CFX96	<i>Histomonas meleagridis</i>	FAM	+	
	IAC	VIC	+	
R-Biopharm RIDA®CYCLER	<i>Histomonas meleagridis</i>	green	+	
	IAC	yellow	+	
LTF MyGo Pro	<i>Histomonas meleagridis</i>	FAM	+	
	IAC	VIC	+	
Qiagen Rotor-Gene Q	<i>Histomonas meleagridis</i>	green	+	
	IAC	yellow	+	
Roche LightCycler® 2.0	<i>Histomonas meleagridis</i>	530	None	The SureCC Color Compensation Kit II (Art. No. F4010) is required. Check the Seek Temperature is 58°C.
	IAC	560	None	
Roche LightCycler® 480 II	<i>Histomonas meleagridis</i>	465-510	+	
	IAC	533-580	+	
Roche cobas® z 480 Analyzer	<i>Histomonas meleagridis</i>	465-510	+	
	IAC	540-580	+	

2 Qualitative Analysis

2.1 Protocol

2.1.1 DNA-preparation

Transfer 50 mg tissue sample into a 2.0 ml reaction tube (not provided with the kit) and add 500 µl of Lysis Buffer (**Code L**) to this sample.

Vortex the tube briefly and put the tube in the heating block.

Incubate for 10 minutes at 95 °C without shaking.

Remove the tube from the heating block and leave the tube at room temperature for 1 minute.

Note: It is necessary to ensure that the sediment is not stirred up and no particles are pipetted in the PCR reaction.

The lysate is ready-to-use for PCR. If the lysate is not used immediately in the PCR or is intended for a longer storage, transfer 100 µl of the lysate in a new tube (not provided with the kit) and store at -20 °C.

2.1.2 Preparation of the master-mix

Calculate the total number of reactions needed (samples and control reactions) for the specific PCR assay as well as for the inhibition control.

Recommended control reactions for the specific PCR assay: negative control, extraction control, positive control. For the analysis of enrichments additional controls are recommended: zero control (sample before enrichment) and medium control. The reaction mix contains an internal amplification control (IAC) per reaction.

Reactions needed for the qualitative *Histomonas meleagridis* detection:

3 reactions for controls (1x no-template control, 1x extraction control, 1x positive control)

For each sample: at least 1 reaction for each sample DNA

It is also recommended to prepare the master-mix with 10 % additional volume in order to compensate reagent loss. Allow the reagents to thaw, mix and centrifuge before opening and use.

Example for the calculation and preparation of 10 reactions:

Components of the master-mix	Amount per reaction	10 reactions (with 10% excess)
Reaction Mix	19.3 µl	212.3 µl
Taq Polymerase	0.7 µl	7.7 µl
Total volume	20 µl	220 µl

Mix each master-mix well and centrifuge shortly before use.

2.1.3 Preparation of the real-time PCR-mix

- Pipette 20 µl of the master-mix into appropriate tubes/wells.
- Close the negative control (the negative control is ready for PCR without any addition).
- Pipette 5 µl of sample DNA into the designated tubes/wells and close them.
- Pipette 5 µl of Positive Control into the designated tubes/wells and close them.
- Centrifuge all tubes/plates or capillaries shortly at low speed.
- Place tubes/plates into the real-time PCR instrument and start the run according to the setup.

2.2 Interpretation of results

The evaluation has to be made according to the usual analysis program recommended by the real-time PCR instrument manufacturer.

The control reactions have to show the correct results.

Histomonas meleagridis DNA is detected in the FAM-channel. In the VIC/HEX-channel the amplification control is detected.

A sample is stated **positive** for *Histomonas meleagridis*, if the sample DNA shows amplification in the FAM-channel.

A sample is stated **negative** for *Histomonas meleagridis*, if the sample DNA shows no amplification in the FAM-channel and if the internal control (VIC/HEX-channel) of the sample is **positive** with a shift in Cp-Value ≤ 2 compared to the negative control.

If the sample DNA in the VIC/HEX-Channel shows **no amplification** or a shift in Cp-value > 2 compared to the negative control, it contains PCR inhibiting substances. A significant decrease in the fluorescence signal can also show the presence of PCR inhibiting substances. Under these circumstances DNA isolation and purification of the sample need to be improved. Alternatively the DNA can be diluted (recommendation 1:2 in PCR-water) and analysed again for inhibition. Please note that the dilution factor also affects the detection limit of the specific *Histomonas meleagridis* PCR assay.

3 Further Information

3.1 Product Information

- Detailed information about setup of several real-time PCR devices
(Download: www.congen.de/en/company/downloads)
- Validation Report upon request

3.2 Technical Support

For further questions please send an e-mail to info@congen.de.