CONGEN

SureFood[®] ANIMAL ID 3plex Rat/Mouse+IAAC

Art. No. S6127
100 rxn

User Manual

June 2020

 **Inhalt**

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Content

1	General Information	8
1.1	Description	8
1.2	Limit of Detection	8
1.3	Cross reactivity	8
1.4	DNA-preparation	8
1.5	Kit components and storage	8
1.6	Additionally required equipment and materials	8
1.7	Setup.....	9
1.8	Detection channel Set-up	9
2	Qualitative Analysis	10
2.1	Protocol	10
2.1.1	Preparation of the master-mix	10
2.1.2	Preparation of the real-time PCR-mix	10
2.2	Interpretation of results	11
3	Further Information	12
3.1	Product Information	12
3.2	Technical Support	12
3.3	Distribution and Ordering	12

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1 General Information

1.1 Description

The SureFood® ANIMAL ID 3plex Rat/Mouse+IAAC is a real-time PCR for the direct, qualitative detection and differentiation of specific Brown rat (*Rattus norvegicus*)/Black rat (*Rattus rattus*) and House mouse (*Mus musculus*) DNA sequences.

Each reaction contains an internal amplification control and an internal detection assay for vertebrates DNA (IAAC). If the DNA contains PCR inhibiting substances, the signal of the amplification control will be affected or the amplification will be suppressed. Examples for PCR inhibiting substances are alcohols (e.g. ethanol, isopropanol), surfactants (e.g. CTAB, SDS, Triton X100) and salts (e.g. sodium chloride). In addition spices, herbs, algae, cocoa and further sample matrices might have PCR inhibiting effects.

The real-time PCR assay can be performed with commonly used real-time PCR instruments, equipped for detection of three fluorescence emissions at the channels FAM, VIC/HEX and Cy5 at the same time. The technical validation of instruments was performed on Agilent Mx3005P, Agilent AriaDx, Bio-Rad CFX96, BioRad CFX 96 Dx, R-Biopharm RIDA®CYCLER, Roche LightCycler® 480 II and Qiagen Rotor-Gene Q.

1.2 Limit of Detection

The SureFood® ANIMAL ID 3plex Rat/Mouse+IAAC real-time PCR is developed for the detection of rat and mouse DNA in muscle meat mixture at a relative amount of ≤ 500 DNA copies.

The limit of detection of the complete method (DNA extraction and real-time PCR) depends on sample matrix, processing grade, DNA-preparation and DNA-content.

The SureFood® PCR systems are very sensitive and therefore even a small amount of target DNA is sufficient for a successful analysis. The concentration of total DNA in the sample does not allow a conclusion on the quantity and quality of the target DNA.

Note: In mixed samples, inconsistent mixing ratios (e.g. 99.9 % rat and 0.1 % mouse) may cause a loss of sensitivity in the low concentration channel.

1.3 Cross reactivity

Cross reactivity was observed with DNA extracts from Turkestan rat (*Rattus pyctoris*), Asian house rat (*Rattus tanezumi*) and Algerian mouse (*Mus spretus*).

1.4 DNA-preparation

For DNA-preparation of raw material the use of SureFood® PREP Basic (Art. No. S1052) and for highly processed food and feed the use of SureFood® PREP Advanced (Art. No. S1053) is recommended.

1.5 Kit components and storage

Kit Code	Reagent	Amount	Lid Color
1	Reaction Mix	2 x 1100 µl	Yellow
2	Taq Polymerase	1 x 11 µl	Dark Red
3	Positive Control	1 x 200 µl	Light Blue

Store all reagents at –20°C and protected from light.

1.6 Additionally required equipment and materials

- DNA-Extraction kit
(e.g. SureFood® PREP Basic Art. No. S1052 / SureFood® PREP Advanced Art. No. S1053)
- real-time PCR instrument with three detection channels (510 nm, 580 nm and 660 nm)
- real-time PCR consumables (plates, tubes, foils, caps)

SureFood[®] ANIMAL ID 3plex Rat/Mouse+IAAC (100 rxn)

Art. No. S6127

June 2020

- pipettes with filter tips
- powder-free disposable gloves
- Vortex mixer
- micro centrifuge with a rotor for the reaction tubes

1.7 Setup

	Blockcycler & R-Biopharm RIDA[®]CYCLER	Rotorcycler
Initial Denaturation (HOLD)	5 min, 95°C	1 min, 95°C
Cycles	35	35
Denaturation	15 sec, 95°C	10 sec, 95°C
Annealing/Extension (CYCLE)	30 sec, 60°C	15 sec, 60°C
Temperature Transition Rate/ Ramp Rate	Maximum	Maximum

1.8 Detection channel Set-up

Real-time PCR device	Detection	Detection channel	Quencher	Note
Agilent Mx3005P	Rat	FAM	+	
	IAAC	HEX	+	
	Mouse	Cy5	+	
Agilent AriaDx	Rat	FAM	+	
	IAAC	HEX	+	
	Mouse	Cy5	+	
Bio-Rad CFX96	Rat	FAM	+	
	IAAC	VIC/HEX	+	
	Mouse	Cy5	+	
Bio-Rad CFX96 Dx	Rat	FAM	+	
	IAAC	VIC/HEX	+	
	Mouse	Cy5	+	
R-Biopharm RIDA[®]CYCLER	Rat	green	+	
	IAAC	yellow	+	
	Mouse	red	+	
Qiagen Rotor-Gene Q	Rat	green	+	Note: Please use only 0.1 ml reaction tube. The gain settings must be set to 5 (factory default) for all channels.
	IAAC	yellow	+	
	Mouse	red	+	
Roche LightCycler[®] 480 II	Rat	465-510	+	
	IAAC	533-580	+	
	Mouse	618-660	+	
Roche cobas[®] z 480 Analyzer	Rat	465-510	+	
	IAAC	540-580	+	
	Mouse	610-670	+	

2 Qualitative Analysis

2.1 Protocol

2.1.1 Preparation of the master-mix

Calculate the total number of reactions needed (samples and control reactions) for the specific PCR assay as well as for the inhibition control.

Recommended control reactions for the specific PCR assay: negative control, extraction control, positive control. The reaction mix contains an internal amplification control (IAC) per reaction.

Reactions needed for the qualitative detection of rat and mouse:

3 reactions for controls (1x no-template control, 1x extraction control, 1 positive control)

For each sample: at least 1 reaction for each sample DNA

It is also recommended to prepare the master-mix with 10 % additional volume in order to compensate reagent loss. Allow the reagents to thaw, mix and centrifuge before opening and use.

The tube of the Taq Polymerase should be kept at -20°C and not be mixed by vortexing.

Example for the calculation and preparation of 10 reactions:

Components of the master-mix	Amount per reaction	10 reactions (with 10% excess)
Reaction Mix	19.9 µl	218.9 µl
Taq Polymerase	0.1 µl	1.1 µl
Total volume	20 µl	220 µl

Mix each master-mix well and centrifuge shortly before use.

2.1.2 Preparation of the real-time PCR-mix

- Pipette 20 µl of the master-mix into appropriate tubes/wells.
- Close the negative control (the negative control is ready for PCR without any addition).
- Pipette 5 µl of sample DNA into the designated tubes/wells and close them.
- Pipette 5 µl of Positive Control into the designated tubes/wells and close them.
- Centrifuge all tubes/plates shortly at low speed.
- Place tubes/plates into the real-time PCR instrument and start the run according to the setup.

2.2 Interpretation of results

The evaluation has to be made according to the usual analysis program recommended by the real-time PCR instrument manufacturer.

The control reactions have to show the correct results.

Rat DNA is detected in the FAM-channel and mouse DNA is detected in the Cy5-channel. In the VIC/HEX-channel it is possible to detect animal DNA in the sample as well as the amplification control (IAC) in a sample with no animal DNA inside. (see also table below)

A sample is stated **positive** for the respective parameter, if the sample DNA shows amplification in the respective channel.

A sample is stated **negative** for the respective parameter, if the sample DNA shows no amplification in the respective channel. The internal amplification control (inhibition control) of the sample has to be **positive** with a shift in Cp-Value ≤ 2 compared to the negative control in the VIC/HEX-channel.

If the internal control signal in the VIC/HEX-channel of the sample DNA is detected significantly before the signal of the negative control the sample is stated **positive** for animal DNA.

Is the Cp-value of the internal control in the VIC/HEX-channel in the range of the negative control the sample contains no PCR-inhibiting substances but only a low amount or no animal DNA.

If the sample DNA in the VIC/HEX-channel shows **no amplification** or a shift in Cp-value > 2 compared to the negative control, it contains PCR inhibiting substances. A significant decrease in the fluorescence signal can also show the presence of PCR inhibiting substances. Under these circumstances DNA isolation and purification of the sample need to be improved. Alternatively the DNA can be diluted (recommendation 1:2 in PCR-water) and analysed again for inhibition. Please note that the dilution factor also affects the detection limit of the specific rat or mouse PCR assay.

result in the respective channel			Interpretation
FAM channel Rat	Cy5 channel Mouse	VIC/HEX channel animal + IAC	
positive	negative	positive	Rat DNA detected
negative	positive	positive	Mouse DNA detected
negative	negative	positive	vertebrates DNA detected
negative	negative	negative	invalid

Note: If the sample contains more than one animal species the DNA mixture can have a competitive influence on the absolute fluorescence. The lower the concentration of the determinant DNA is in a mixture of animal DNAs the lower is the fluorescence level of the amplification curve.

3 Further Information

3.1 Product Information

- Microsoft Excel template of calculation and detailed information about setup of several real-time PCR devices (Download: www.congen.de/en/company/downloads)
- Validation Report upon request

3.2 Technical Support

For further questions please contact your distributor or send an e-mail to sales@r-biopharm.de.

3.3 Distribution and Ordering

R-Biopharm AG
An der neuen Bergstrasse 17,
64297 Darmstadt, Germany
Phone: +49 (0) 61 51 - 81 02-0
Fax: +49 (0) 61 51 - 81 02-20
E-Mail: orders@r-biopharm.de
www.r-biopharm.com